

VELP Scientifica Solutions: Seal for SER 158

TYPE OF SOLVENTS	Type of Seal	Heating Level of Glass Beaker	Boiling Stone	Boiling Point (°C)	Ignition Temperature (°C)
Acetone	Grey Butyl Seal A00000298 	4	Required	56,5	538
Ethyl Acetate	Grey Butyl Seal A00000298 	8	Required	77	427
Ethanol	Grey Butyl Seal A00000298 	9	Required	78,5	423
Methanol	Grey Butyl Seal A00000298 	9	Required	64,7	470
Chloroform	Green Viton Seal A00000298 	4	Required	61-62	Not Combustible
Petroleum ether	Green Viton Seal A00000298 	3	-	30-80	Not Specified
n-Hexane	Green Viton Seal A00000298 	4	-	69	234
Diethyl ether	White Vaflon Seal A00000298 	1	-	34,6	180-190
Toluene	White Vaflon Seal A00000298 	12	Required	110,6	536
Acetone n-Hexane 1:1	White Vaflon Seal A00000298 	4	Required	-	-
Acetone n-Hexane 1:3	White Vaflon Seal A00000298 	4	Required	-	-

