Instruction Manual
Manuale di istruzioni
Manuel d'instructions
Manual de instrucciones
Bedienungsanleitung

FOC 120E Connect Cooled Incubator

F10300540

FOC 120I Connect Cooled Incubator

F10400541

FOC 200E Connect Cooled Incubator

F10300542

FOC 200I Connect Cooled Incubator

F10400543

FOC 200IL Connect Cooled Incubator

F10400544

General Information / Informazioni Generali / Informations Générales / Información General / Allgemeine Hinweise

Before using the unit, please read the following instruction manual carefully.

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There is a small quantity of isobutane gas (R600a) in the refrigerating system.

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Do not dispose of this equipment as urban waste, in accordance with EEC directive 2012/19/UE.

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This unit must be used for laboratory applications only.

The manufacturer declines all responsibility for any use of the unit that does not comply with these instructions.

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It is responsibility of the user appropriately decontaminate the instrument in case of dangerous substances fall on or in it accordingly to the safety datasheet of substances used and to the current laboratories safety standards. It is not possible to decontaminate the product under steam.

It is also responsibility of the user to use substances for cleaning or decontaminating which do not react with internal parts of the instrument or with the material contained in it. In case of doubts on the compatibility of a cleaning solution, contact the manufacturer or local distributor.

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Beware that liquids don't get in touch with power supply or with internal electronic components.

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Use only power cable with grounding.

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For the electrical connection, carefully follow the instructions provided in the following manual. It is advisable to wait at least four hours before connecting the appliance to allow the liquid refrigerant gas inside the refrigerant circuit to flow into the compressor.

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Adequate air circulation should be around the appliance, failure leads to abnormal overheating which can cause performance loss.

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If possible, the back of the product must be against a wall to avoid touching hot parts (compressor and condenser).

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The instrument must not be placed near radiators or heat sources

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Never cover the condenser fins on the back of the instrument.

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The unit offers excellent thermal stability and homogeneity at all set temperatures thanks to the dedicated “Auto Tuning” thermoregulation system.

The incubator constantly monitors also the ambient temperature and it adjusts the power of the heating and cooling system accordingly.

The internal temperature is measured by the internal probe and it is shown on the display.

The unit is equipped with two internal sockets, controlled by a switch on the frontal panel, to power instruments placed inside the incubator (e.g. VELP BOD and respirometers systems).

The FOC 120I, FOC 200I and FOC 200IL incubators have an additional internal transparent door that allows the user to visually check the samples without causing big internal temperature variations.

The FOC 200IL is also equipped with two lighted shelves. Each shelf has six led bars horizontally mounted on the lower part which illuminate the below shelves. Furthermore, a dedicated timer for the internal lightening allows the user to create several operating cycles with automatic switch off and on.

The integrated Wi-Fi module transmits and receives information in order to monitor and control the unit through the VELP Ermes cloud platform.

In accordance with current regulations, all VELP incubators are CFC-free (free from chloro fluoro carbons).

2. Installation

- Unpacking
 - Check the integrity of the unit after unpacking

- The box includes
 - FOC Cooled Incubator
 - Instruction manual
 - Power supply cable

- Installation
 - Make sure that the technical specification of the unit and the power supply line are the same
 - Connect the instrument to an easily accessible socket (compliant with the current safety norms), using only the provided power supply cable

NOTE: *if the instrument has been transported or stored horizontally, leave it in the upright position for at least 24 hours before switching it on. This will allow the cooling circuit to settle. If this procedure is not carried out, the refrigerating system could be damaged. This damage is not covered by warranty.*

NOTE: *do not install the equipment near sources of heat (ovens, etc.). The unit should be installed in a well-ventilated area, avoid niches or small rooms.*

3. Operation

Turn the unit on using the main switch on the frontal panel.

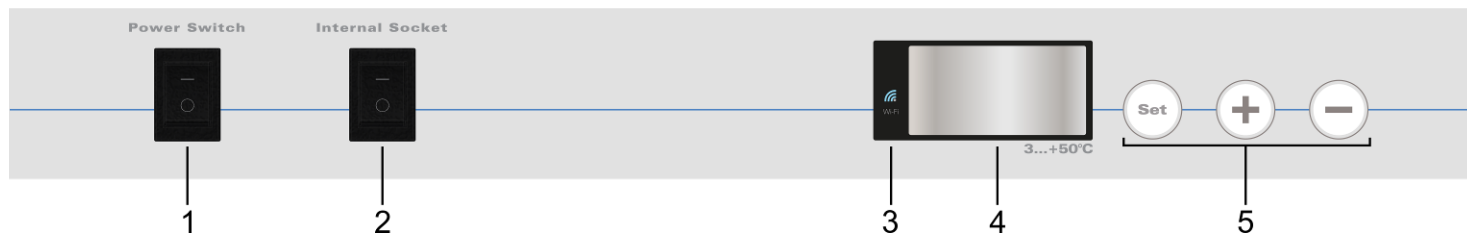
The display shows the software version and the last setpoint value. Automatically, the unit starts the thermoregulation to achieve and maintain the set temperature.

The internal temperature is measured by the internal probe and it is shown on the display.

To adjust the working temperature, press the “Set” button. Then use the “+” and “-” buttons to increase or decrease the setpoint value, respectively (by keeping the buttons pressed, the setpoint can be adjusted faster).

The new setpoint is confirmed by pressing again the “Set” button or automatically if no buttons are pressed.

Use the dedicated switch on the frontal panel to control the internal sockets.

- 1 Main switch
- 2 Internal sockets switch
- 3 Wi-Fi led
- 4 Display
- 5 Control buttons

Placing products / samples in the incubator

Do not place hot products in the incubator; otherwise, the average internal temperature will raise and the other products inside can be damaged.

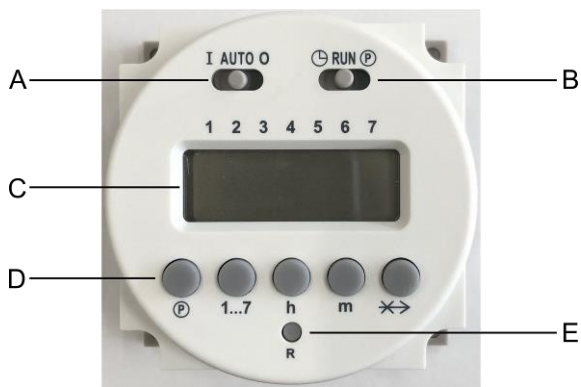
Do not often open the door to avoid big internal temperature variations and to reduce energy consumption.

To ensure good air circulation within the unit, avoid placing large quantity of products and do not place objects against the inside back panel.

Do not place any product too close to the temperature probe.

Connect only 230 V (current consumption 3 A max.) instruments to the internal sockets.

Programmable timer for internal lightening (only on FOC 200 IL)

- A Sliding switch for operating mode selection
- B Sliding switch for time and programs setting
- C Timer display
- D Control buttons
- E Reset button

- Current day and time setting:
 - Move the sliding switch “B” to the left position (🕒)
 - Use the “h” and “m” buttons to set the current hours and minutes, respectively
 - Use the “1...7” button to set the current week day
1 = Monday, 2 = Tuesday, ..., 7 = Sunday
An arrow on the display indicates the selection
 - Move the sliding switch “B” to the middle position (RUN).
- Operating mode selection:
 - Move the sliding switch “A”
AUTO: the lights will be turned on and off according to the set operating cycle
I: the lights are always on
O: the lights are always off
- Operating cycle setting:

It is possible to set up to 16 commutations (8 switching ON and 8 switching OFF).
The internal lights will be turned on at switching ON commutation times. Then, they will be turned off at switching OFF commutation times.
The switching ON commutation are identified with an odd number (1, 3, 5, ..., 15) and the lamp bulb symbol (💡).
The switching OFF commutation are identified with an even number (2, 4, 6, ..., 16).

 - Move the sliding switch “B” to the right position (🕒)
 - Press the “1...7” button to select the week day or a group of days:
1 = Monday, 2 = Tuesday, ..., 7 = Sunday
1 – 5 = From Monday to Friday 1 – 6 = From Monday to Saturday
6 – 7 = Saturday and Sunday 1 – 7 = The whole week
Arrows on the display indicate the selection.
 - Use the “P” button to select the desired commutation. Then use the “h” and “m” buttons to set the commutation time
Repeat these steps to set the needed commutations.

NOTE: the commutation identification number is showed in the right part of the display
NOTE: to delete a commutation press the “P” and the “↔” buttons for few seconds. The display will show “0:00”

 - Move the sliding switch “B” to the middle position (RUN).

NOTE: to avoid error, any switching ON commutation should be followed by a switching OFF commutation.

- Commutation swap function:

When an operating cycle is in progress, by pressing the “↔” button it is possible to swap the lights status.
If they are on, they will be switched off and vice versa.
By pressing again the same button, the light status will be restored as set in the operating cycle.

NOTE: the swap works only until the next commutation. Then, it will be cancelled and the operating cycle restored.

4. Setting mode

Press the “Set” button for 3 seconds to access the setting mode.

Use the “+” and “-” buttons to browse through the available menus, listed below.

Press the “Set” button to access the selected menu. Then, use the “+” and “-” buttons to modify the value of the parameter (when it is possible).

Press the “Set” button to confirm.

To exit the setting mode, do not press any buttons for 5 seconds.

Text shown on the display	Default value	Range	Menu's name and description
<i>CAL</i>	- 0.5 °C	- 9,9 ÷ 9,9 °C 0,1 °C steps	<u>Probe calibration</u> It allows to align the temperature probe reading to a reference thermometer
<i>ULI</i>	50 °C	<i>LLI</i> ÷ 50 °C 0,1 °C steps	<u>Setpoint upper limit</u> It limits the maximum value of temperature setpoint
<i>LLI</i>	3 °C	3 ÷ <i>ULI</i> °C 0,1 °C steps	<u>Setpoint lower limit</u> It limits the minimum value of temperature setpoint
<i>CCO</i>	---		<u>Compressor, heating element and fan operating time</u> It shows compressor, heating element and fan operating times. Operating times are showed in hours up to 999 hours (around 42 days). Then they are showed in days using different combinations of the digits' decimal point.
<i>CHE</i>			
<i>CFA</i>			
<i>UFI</i>	NO	YES – NO	<u>Wi-Fi connection</u> It allows to enable / disable the Wi-Fi connection
<i>OtC</i>	---		<u>One time code</u> It shows the One Time Code during the registration of the device on the VELP Ermes cloud platform
<i>AP</i>	NO	YES – NO	<u>Access point</u> It allows to enable / disable the Access Point mode of the Wi-Fi connection
<i>rSt</i>	NO	YES – NO	<u>Reset</u> It allows to reset all the system settings to the default values

5. Error and warning messages

When the display shows an error message, the instrument's functions are stopped automatically.

AL1	Internal temperature too low	Warning messages only on Ermes platform
AL2	Internal temperature too high	
AL3	Internal probe broken	
AL4	Time to reach setpoint too high	
AL5	Big internal temperature variation	

To remove the error message, restart the instrument.

If the alarm persists on the display, please contact VELP Scientifica's technical service department.

6. VELP Hermes cloud platform configuration

VELP Hermes is a revolutionary cloud platform that transforms and improves your laboratory experience.

VELP Hermes creates a connected ecosystem of devices, people and data that transmit information between each other cutting down distances and expanding your scientific potential.

VELP Hermes collects and stores your data with the maximum level of encryption following the highest cyber-security standards.

- Activate VELP Hermes account:
 - Go to www.velp.com through your internet browser
 - Create your VELP account or login if you already have it
 - Access to your profile page and select “Configure Your VELP Hermes Account”. Activate your VELP Hermes account after having accepted Terms and Conditions.
 - You will receive an email containing your ID, Password and a link to proceed to VELP Hermes authentication at the email address that you entered during the registration
 - At the first login on VELP Hermes platform, a new password will be requested

- Connect the incubator to your Wi-Fi network:
 - Turn on the FOC Connect
 - Enable the Wi-Fi connection and the access point mode as described above
 - Using a PC, tablet or smartphone select the *VELP_FOC_(Serial Number)* network available on the Wi-Fi list
 - Open a web browser on your PC, tablet or smartphone; insert 192.168.4.1 into the address bar and press “Enter” Enter “admin” as User ID and “velp_ermes” as Password
 - Configure the laboratory Wi-Fi network parameters. Then press “Push” to confirm.
NOTE: if necessary, contact your IT specialist to get all your laboratory Wi-Fi connection details
 - Restart the FOC Connect and wait until the Wi-Fi led stops blinking (it has to be constantly ON)
NOTE: if the led is still blinking after 2 minutes, please try to repeat the procedure and check that the network parameters are correct

- Register the incubator on VELP Hermes platform:
 - On the incubator, access the One Time Code (OTC) setting menu as described above
 - On the PC, tablet or smartphone, login on VELP Hermes with your account and select “Device Registration”
 - Enter the FOC Connect product code, serial number and purchase date.
Press “Next”
 - Enter the One Time Code showed on the display of the incubator (the first numbers are preceded by the letter “H” and the last ones by the letter “L”).
Press “Next” and wait for the connection
NOTE: if the OTC is not showed on the display of the incubator, please try to repeat the procedure and check that the network parameters are correct
 - Give a name to the FOC Connect and enter the address where the instrument is located
Press “Next”
 - Assign the FOC Connect to an existing virtual laboratory or create a new one through “Laboratory Management”

At the end of the procedure, you will be able to monitor and control the connected incubator through “Instruments” menu.

NOTE: the incubator can be connected only to a 2.4 GHz Wi-Fi network.

NOTE: for further information, please see VELP website.

7. MODBUS serial communication

The incubators are equipped with two RS485 female connectors which can be used for MODBUS serial communication. The table below contains the registers assignment:

Register address	Register name	Write	Byte number	Function
1	Instrument	NO	2	Instrument model
2	Serial number	NO	6	Instrument serial number
5	Product code	NO	12	Instrument product code
11	Mainboard SW	NO	8	Mainboard software version
24	Status	NO	2	Instrument status (heating/cooling)
25	Alarm	NO	2	Error message ID
27	Temperature	NO	2	Current internal temperature
33	Temperature setpoint	NO	2	Current temperature setpoint
50	Setpoint lower limit	NO	2	Current setpoint lower limit
51	Setpoint upper limit	NO	2	Current setpoint upper limit
101	Modbus address	YES	1	Modbus address
103	Temperature	YES	2	Working temperature setting

Virtual serial port:

Baudrate [bps]	9600
Data bits	8
Stop bit	1
Parity	None
Maximum number of registers for single request	24
Sampling time	≥ 1 second

Serial port pinout:

2	Data + (A)
3	GND
4	Data – (B)

Examples (16 bit CRC for standard MODBUS)

Reading single register: Temperature (°C)

Request		Reply		Description
Field	(Hex)	Field	(Hex)	
Address	0x64	Address	0x64	
Control command	0x03	Control command	0x03	
High start address	0x00	Number of bytes	0x02	
Low start address	0x1B	High register value	0x00	25 °C (250 / 10)
Number of High registers	0x00	Low register value	0xFA	
Number of Low registers	0x01	High CRC	0x74	
High CRC	0xFD	Low CRC	0x0F	
Low CRC	0xF8			

Reading multiple registers: Serial Number

Request		Reply		Description
Field	(Hex)	Field	(Hex)	
Address	0x64	Address	0x64	
Control command	0x03	Control command	0x03	
High start address	0x00	Number of bytes	0x06	
Low start address	0x02	High register value	0x31	12345 0x00 0x31 0x32 0x33 0x34 0x35
Number of High registers	0x00	Low register value	0x00	
Number of Low registers	0x03	High register value	0x33	
High CRC	0xAD	Low register value	0x32	
Low CRC	0xFE	High register value	0x35	
		Low register value	0x34	
		High CRC	0x0A	
		Low CRC	0x0A	

Writing single register: Working Temperature (°C)

Request		
Field	(Hex)	Description
Address	0x64	
Control command	0x06	
High start address	0x00	
Low start address	0x67	
Number of High registers	0x00	5 °C (50 / 10)
Number of Low registers	0x32	
High CRC	0xB0	
Low CRC	0x35	

8. Maintenance

No routine or extraordinary maintenance is necessary except periodically cleaning.

Cleaning

Disconnect the unit from the power supply and use a cloth dampened with a non-flammable non-aggressive detergent.

Repair

Repairs must be carried out by authorized Velp personnel only.

The instrument must be transported using the original shockproof packaging when freight forwarders, couriers or other are taking care of the unit transport. Follow the indications on the original packaging (e.g. place the instrument on a pallet).

It is the responsibility of the user, to properly decontaminate the unit in case of hazardous substances remaining on the surface or interior of the device. If in doubt about the compatibility of a cleaning or decontamination product, contact the manufacturer or distributor.

9. Technical data

Model	FOC 200E and FOC 200I	FOC 200IL	FOC 120E and FOC 120I
Power supply	230 V - 50/60 Hz		
Usable internal volume	169 l		109 l
Power	400 W		120 W
Dimensions (WxHxD)	540x1263x550 mm (21.3x49.7x21.7 in)	540x1300x550 mm (21.3x51.2x21.7 in)	540x912x550 mm (21.3x35.9x21.7 in)
Weight	46.3 Kg (102.1 lbs)	50 Kg (110.2 lbs)	36 Kg (79.4 lbs)
Internal temperature visualization	3-digits display		
Temperature setting	Digital		
Temperature range setting	3.0 – 50.0 °C		
Temperature probe	Pt100		
Display resolution	0.1 °C		
Sample temperature stability	± 0.5°C		
Sample homogeneity temperature	± 0.5°C		
Heating regulator type	P.I.D.		
Cooling regulator type	Electronic ON-OFF		
Thermostatting control type	Auto Tuning		
Interfaces	Wi-Fi, 2 x RS485		
Max altitude of installation	4000 m (13123 ft) above sea level		
Max humidity admitted	80%		
Storage temperature range	-10 - +60 °C (14 – 140 °F)		
Environmental temperature range	5 - 35 °C (41 – 104 °F)		
Luminous flux	-	30000 Lux/shelf	-
No. max switching on and off commutations for the lighted shelves	-	16	-
Max load (Kg) for each shelf	10 kg (Homogeneously)		

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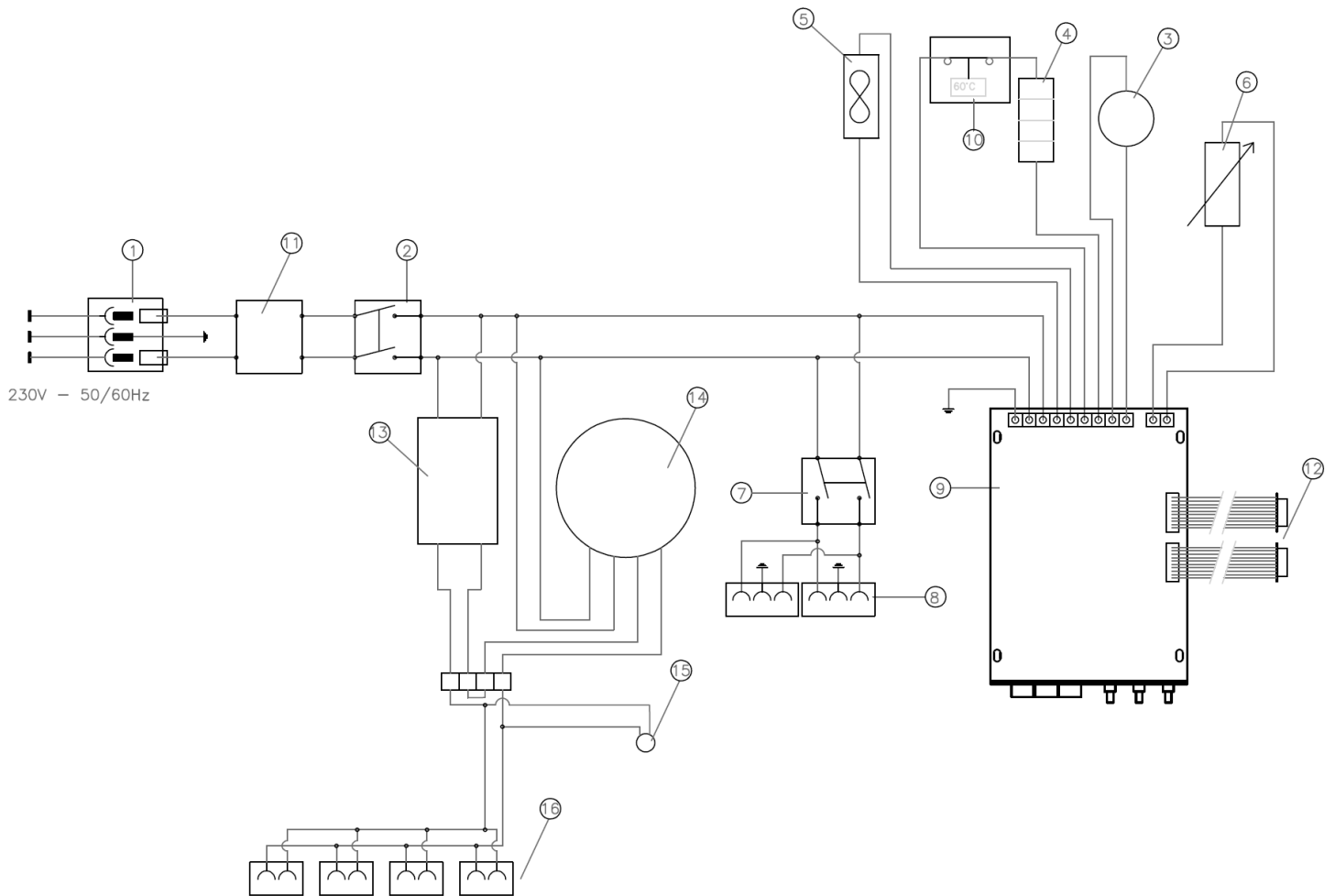
We, the manufacturer VELP Scientifica, under our responsibility declare that the product is manufactured in conformity with the following standards:

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EN 61010-1	Safety requirements for electrical equipment for measurement, control, and laboratory use Part 1: General requirements
EN 61010-2-010	Particular requirements for laboratory equipment for the heating of material
ETSI EN 301 489-1 V2.2.3	Electromagnetic compatibility (EMC) standard for radio equipment and services; Part 1: Common technical requirements
ETSI EN 301 489-17 V3.2.4	ElectroMagnetic Compatibility (EMC) standard for radio equipment and services; Part 17: Specific conditions for Broadband Data Transmission Systems
EN 61326-1	Electrical equipment for measurement, control and laboratory use - EMC requirements - Part 1: General requirements
ETSI EN 300 328 V1.9.1	Wideband transmission systems; Data transmission equipment operating in the 2,4 GHz band

and satisfies the essential requirements of the following directives:

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2006/42/EC	Machinery directive
2014/53/EU	Radio Equipment directive (RED)
2015/863/EU (RoHS III)	Restriction of the use of certain hazardous substances
2012/19/EU (WEEE)	Waste of electric and electronic equipment

plus modifications / più modifiche / plus modifications / más sucesivas modificaciones / in der jeweils gültigen Fassung.

13. Declaration of conformity ^{UK} _{CA}

We, the manufacturer VELD Scientifica, under our responsibility declare that the product is manufactured in conformity with the following regulations:

S.I. 2017/1206 Radio Equipment Regulations 2017

according to the relevant designated standards:

EN 61010-1 Safety requirements for electrical equipment for measurement, control, and laboratory use
Part 1: General requirements

EN 61010-2-010 Particular requirements for laboratory equipment for the heating of material

ETSI EN 301 489-1 V2.2.3 Electromagnetic compatibility (EMC) standard for radio equipment and services;
Part 1: Common technical requirements

ETSI EN 301 489-17 V3.2.4 ElectroMagnetic Compatibility (EMC) standard for radio equipment and services;
Part 17: Specific conditions for Broadband Data Transmission Systems

EN 61326-1 Electrical equipment for measurement, control and laboratory use - EMC requirements -
Part 1: General requirements

ETSI EN 300 328 V1.9.1 Wideband transmission systems; Data transmission equipment operating in the 2,4 GHz
band

and satisfies the essential requirements of regulations:

S.I. 2008/1597 Supply of Machinery (Safety) Regulations 2008

S.I. 2012/3032 Restriction of the Use of Certain Hazardous Substances in Electrical and Electronic
Equipment Regulations 2012

S.I. 2013/3113 Waste Electrical and Electronic Equipment Regulations 2013

plus modifications.

Thank you for having chosen VELP!

Established in 1983, VELP is today one of the world's leading manufacturer of analytical instruments and laboratory equipment that has made an impact on the world-wide market with Italian products renowned for innovation, design and premium connectivity. VELP works according to **ISO 9001**, **ISO 14001** and **ISO 45001** Quality System Certification.

Our instruments are manufactured in Italy according to the IEC 1010-1 and CE regulation.

Our product lines:

Analytical instruments

Elemental Analyzers
Digestion Units
Distillation Units
Solvent Extractors
Fiber Analyzers
Dietary Fiber Analyzers
Oxidation Stability Reactor
Consumables

Laboratory Equipment

Magnetic Stirrers
Heating Magnetic Stirrers
Heating Plates
Overhead stirrers
Vortex mixers
Dispersers
COD Thermoreactors
BOD and Respirometers
Cooled Incubators
Flocculators
Overhead Shakers
Turbidimeter
Open Circulating Baths
Pumps

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VELP Scientifica Srl