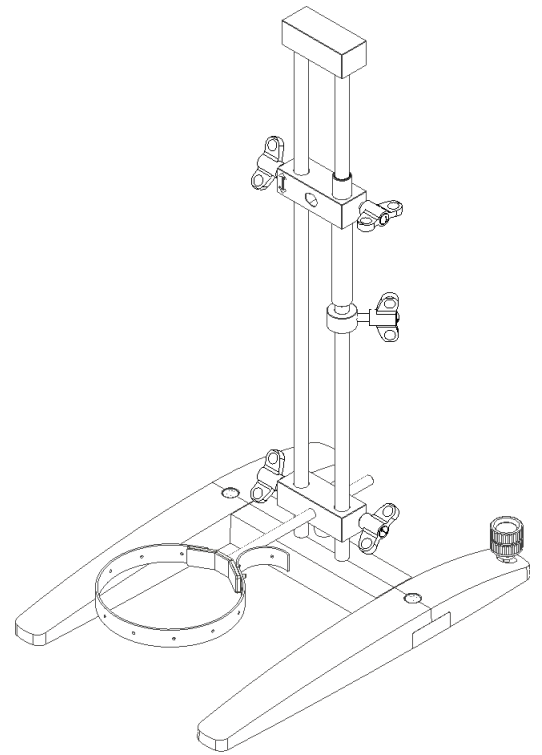
Instruction Manual
Manuale di istruzioni
Manuel d'instructions
Manual de instrucciones
Bedienungsanleitung
指导手册

Double Rod Mini H-Stand Set

A00000484

General Information / Informazioni Generali / Informations Générales / Información General / Allgemeine Hinweise / 一般信息

Before using the unit, please read the following instruction manual carefully.

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This unit must be used for indoor laboratory applications only. The manufacturer declines all responsibility for any use of the unit that does not comply with these instructions. If the product is used in a way not specified by the manufacturer or with non-specified accessories, product's safety may be compromised.

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This unit has been designed and manufactured in compliance with the following standards:

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The vibrations generated by the instrument may loosen the fixing knobs.

Check that the instrument and the sample container are securely mounted in position before use and at regular periods.

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In relation to the many possible applications and configurations, it is not possible to guarantee the safety of the user on the basis of the construction requirements established for the product. Therefore, additional user safety measures may be necessary depending on the specific application.

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The Double Rod Mini H-Stand Set is the ideal solution to firmly fix dispersers and overhead stirrers that require absolute stability during operation.

By adopting a double vertical rod configuration and a new sliding system, the Mini H-Stand is particularly stable and allows for precise and quick positioning of the instrument in the processing container.

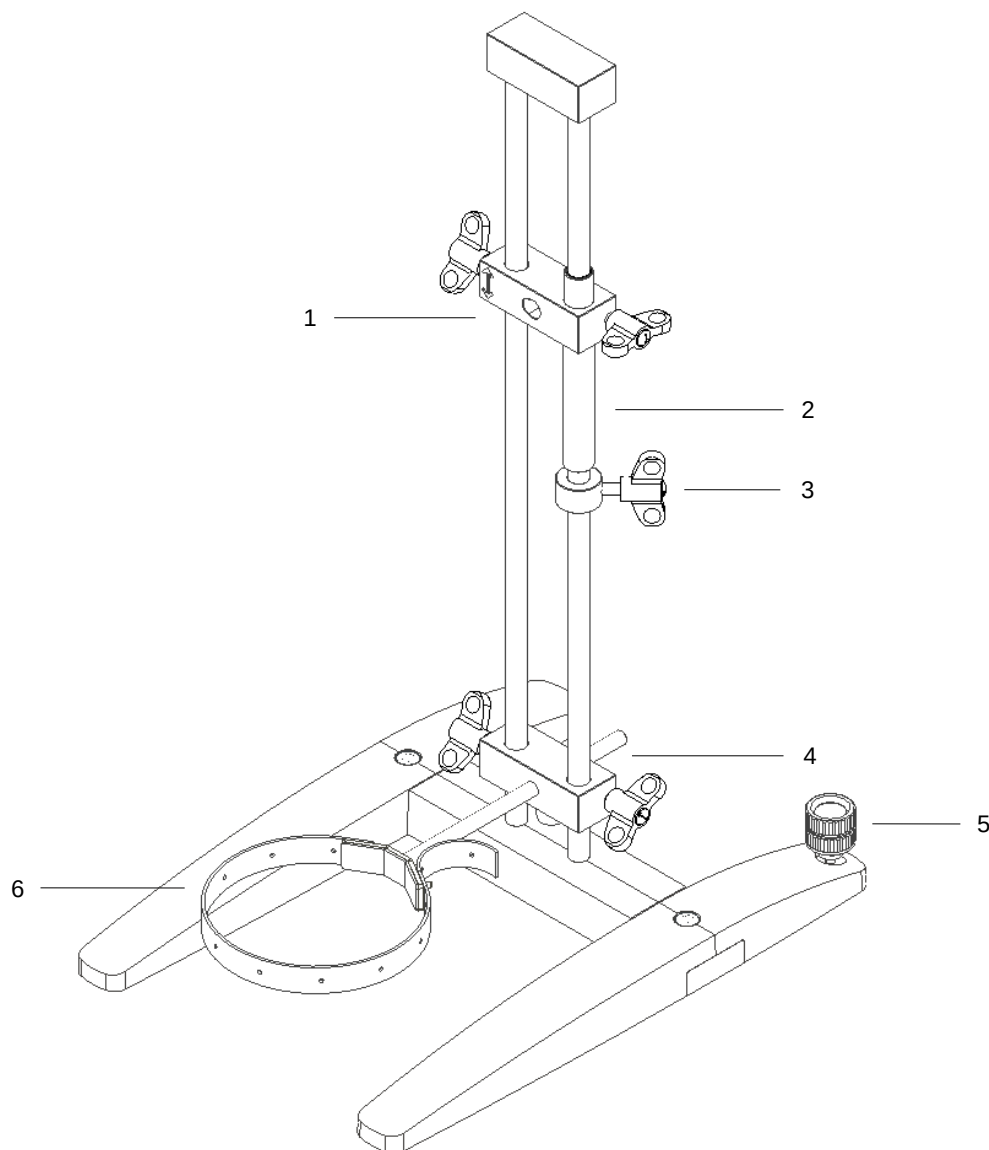
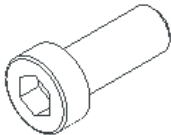
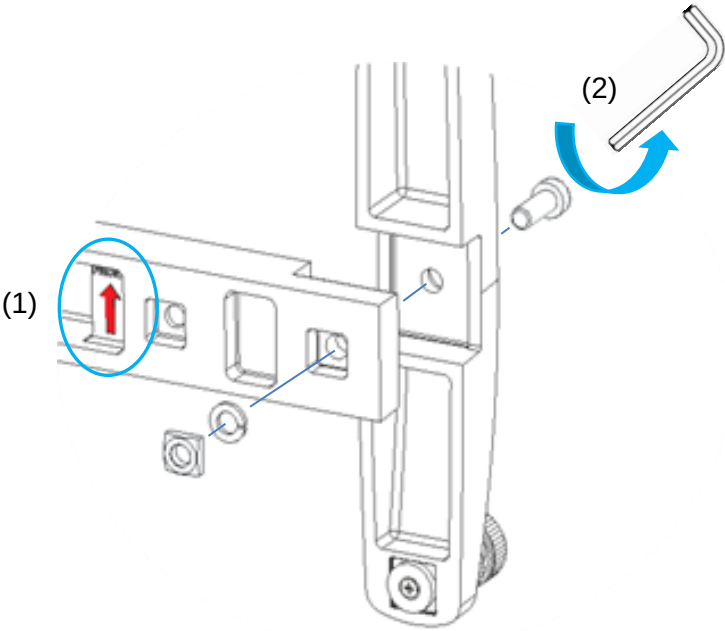
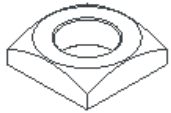
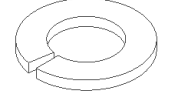
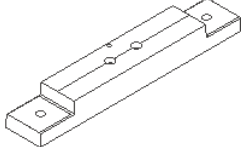
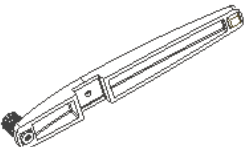
Figure 1

	Description	Function
1	Upper support block	Instrument support
2	Sliding tube	Instrument sliding device
3	Endpoint ring	Instrument vertical movement endpoint
4	Lower support block	Vessel holder with safety belt support
5	Levelling foot	Compensate for unevenness in the setup surface
6	Vessel holder with safety belt	Secure the sample container

2. Assembly

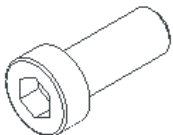
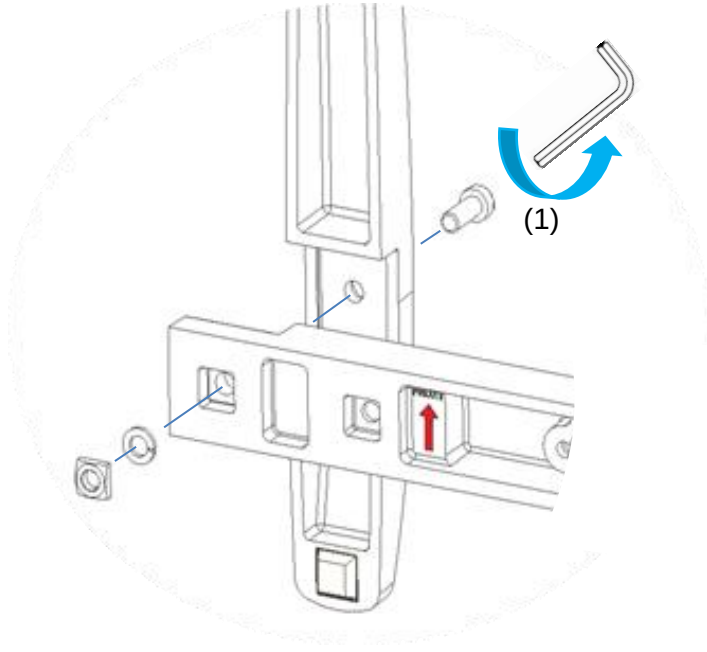
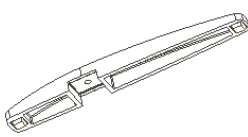
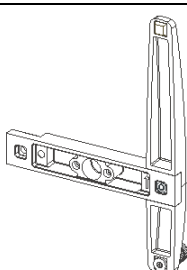
- Unpacking
 - Check the integrity of the unit after unpacking
- The box includes:
 - Double Rod Mini H-Stand Set:
 - Assembled double rod structure
 - Lateral support
 - Lateral support with levelling foot
 - Main support
 - N°2 - Screw M10
 - N°2 - Nut M10
 - N°2 - Spring washer M10
 - N°2 - Screw cover
 - N°2 - Screw M6
 - N°2 - Washer M6
 - N°2 - Spring washer M6
 - Vessel holder with safety belt (Code 40003259)
 - Instruction manual
 - Exagonal key 8 mm
 - Exagonal key 5 mm

1. STEP

N° pieces	Description	Part image	Assembly
1	Screw M10		
1	Nut M10		
1	Spring washer M10		
1	Main support		
1	Lateral support with levelling foot		
1	Exagonal key 8mm		

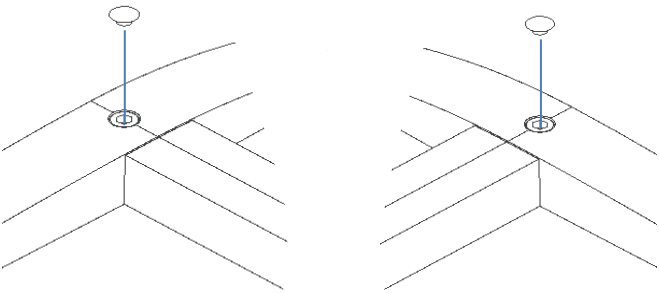
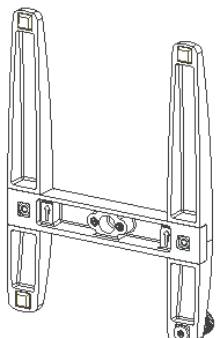
- (1) Position the main support with the arrow pointing forward
 (2) Tighten firmly

2. STEP

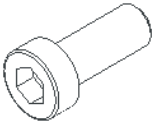
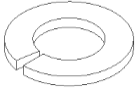
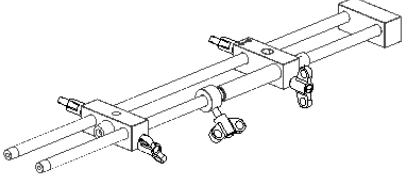
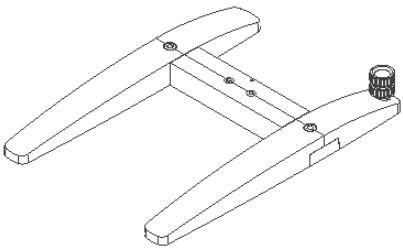
N° pieces	Description	Part image	Assembly
1	Screw M10		
1	Nut M10		
1	Spring washer M10		
1	Lateral support		
1	Pre-assembled base (Step 1)		
1	Exagonal key 8mm		

(1) Tighten firmly

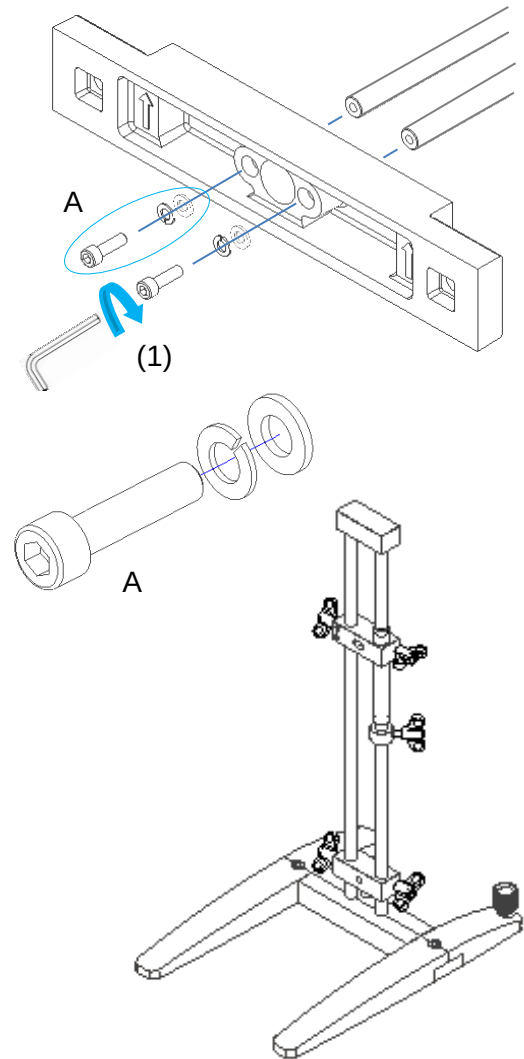
3. STEP

N° pieces	Description	Part image	Assembly
2	Screw cover		
1	Pre-assembled base (Step 2)		

4. STEP

N° pieces	Description	Part image
2	Screw M6	
2	Spring washer M6	
2	Washer M6	
1	Assembled double rod structure	
1	Pre-assembled base (Step 3)	
1	Exagonal key 5mm	

Assembly

(1) Tighten gradually by alternating between the two screws, then tighten firmly

3. Installation

Follow the Safety Regulations described in this Instruction Manual.

- Install the device on a flat, stable, clean, non-slip, dry and fire-resistant surface at 30 cm (at least) from the walls.
 - Adjust the levelling foot (Pos. 5 Fig. 1) to compensate for unevenness on the setup surface
 - Lock the levelling foot in place by tightening the lock nut
- Prior of each use, always check the good condition of the device.

4. Operation

- Fixing the instrument:
 - Move the upper support block to the desired height and tighten the left knob to lock it in place
 - Slide the instrument support shaft into the upper support block
 - Tighten the right knob to lock the instrument at the chosen depth

Note:

Respect the max permissible load as described in Chapter 6 "Technical data".

- Fixing the vessel holder with safety belt (Code 40003259):
 - Move the lower support block to the desired height and tighten the left knob to lock it in place
 - Slide the vessel holder into the lower support block
 - Tight the right knob to lock the vessel holder with safety belt at the chosen depth
- Securing the sample container:
 - Attach the belt to one side of the clamp through one of the first two holes in the belt ("A" or "B" - Fig. 2). The two holes allow for different levels of safety belt tension when securing the container.
 - The vessel holder with safety belt is required to hold the container in place while the instrument is working, thus ensuring safe operation
 - Stretch the tape slightly around the container and secure it to the clamp through the hole that fits best
 - When securing the receptacle, if the belt is too loose or too tight, change the first attachment hole between "A" and "B" (see Fig. 2)

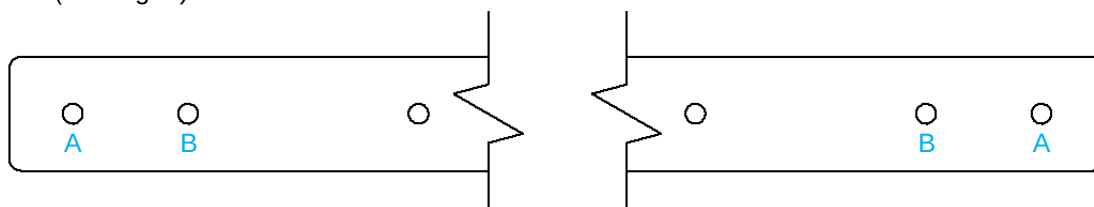
Figure 2 - (Belt tension adjusting – holes "A" and "B")

- Adjusting the Endpoint ring (Pos. 3 Fig. 1):

Note:

The Endpoint ring allows the operator to set an endpoint to the vertical movement of the instrument. In this way, the instrument cannot accidentally hit the bottom of the receptacle containing the sample to be processed, thus ensuring operator safety.

By setting a lower end point for the instrument, the ring also allows to simplify the operation when many samples need to be processed in series.

- Loosen the endpoint ring fixing knob
 - Fix the disperser at the minimum height needed for operation
 - Slide the safety ring upwards until it stops against the sliding tube
 - Firmly tighten the endpoint ring fixing knob
- Adjusting the height of the disperser:
 - Hold the instrument by hand and loosen the left knob of the upper support block
 - Slide the instrument to the desired height and tighten the left knob to lock it in place

Note:

1. *The vibrations generated by the instrument may loosen the fixing knobs. Check that the instrument and the sample container are securely mounted in position before use and at regular periods.*
2. *Before operating on the Mini H-Stand, always turn off the main switch of the instrument*

5. Maintenance and cleaning

No routine or extraordinary maintenance is necessary for the unit.

Cleaning

Disconnect the unit from the power supply (if mounted on the stand) and use a cloth dampened with a non-flammable non-aggressive detergent.

Repair

Repairs must be carried out by authorized Velp personnel only.

The transport of the device by freight forwarders, couriers or others must be carried out using the original shockproof packaging. Follow any instructions on the original packaging (e.g., palletizing).

It is the responsibility of the user to properly decontaminate the unit in case of hazardous substances remaining on the surface or interior of the device. If in doubt about the compatibility of a cleaning or decontamination product, contact the manufacturer or distributor.

6. Technical data

	Model	A00000484
General features	Dimensions (WXHxD):	300x650x430 mm (11,8x25,6x16,9 in)
	Weight	4,9 kg (10,8 lb)
	Construction materials	Aluminum / Stainless steel
	Environmental temperature admitted	+5...+40 °C
	Storage temperature admitted	-10...+60 °C
	Maximum load admitted	5 kg
Vessel holder with safety belt	Vessel diameter	40 – 200 mm

7. Spare parts

40003259	Vessel holder with safety belt
10007983	Adhesive foot

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We, the manufacturer VELP Scientifica, under our responsibility, declare that the product is manufactured in conformity with the following standards:

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EN ISO 12100 Safety of machinery

and satisfies the essential requirements of the following directives:

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2006/42/EC Machinery directive

9. Declaration of conformity

We, the manufacturer VELP Scientifica, under our responsibility declare that the product is manufactured in conformity with the following regulations:

EN ISO 12100 Safety of machinery

and satisfies the essential requirements of regulations:

S.I. 2008/1597 Supply of Machinery (Safety) Regulations 2008

Thank you for having chosen VELP!

Established in 1983, VELP is today one of the world's leading manufacturers of analytical instruments and laboratory equipment that has made an impact on the world-wide market with Italian products renowned for innovation, design and premium connectivity. VELP works according to **ISO 9001, ISO 14001 and ISO 45001** Quality System Certification. Our instruments are manufactured in Italy according to the IEC 1010-1 and CE regulation.

Our product lines:

Analytical instruments

Elemental Analyzers
Digestion Units
Distillation Units
Solvent Extractors
Fiber Analyzers
Dietary Fiber Analyzers
Oxidation Stability Reactor
Consumables

Laboratory Equipment

Magnetic Stirrers
Heating Magnetic Stirrers
Heating Plates
Overhead stirrers
Vortexers and Shakers
Dispersers
COD Thermoreactors
BOD and Respirometers
Cooled Incubators
Flocculators
Turbidimeter
Open Circulating Baths
Pumps

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