EN - Instruction Manual
IT - Manuale di istruzioni
FR - Manuel d'instructions
ES - Manual de instrucciones
DE - Bedienungsanleitung
CN - 指导手册

CONTROLLER Digital

F201A0505

General Information / Informazioni Generali / Informations Générales / Información General / Allgemeine Hinweise / 一般信息

Before using the unit, please read the following instruction manual carefully.
Prima dell'utilizzo dello strumento si raccomanda di leggere attentamente il seguente manuale operativo.
Avant d'utiliser l'instrument, il est recommandé de lire attentivement le présent manuel d'instructions.
Antes de utilizar el instrumento, le recomendamos que lea con atención el siguiente manual de funcionamiento.
Bitte lesen Sie vor Inbetriebnahme des Geräts diese Bedienungsanleitung sorgfältig durch
在使用本装置之前，请仔细阅读以下说明手册

Do not dispose of this equipment as urban waste, in accordance with EEC directive 2002/96/CE.
Non smaltire l'apparecchiatura come rifiuto urbano, secondo quanto previsto dalla Direttiva 2002/96/CE.
Ne pas recycler l'appareil comme déchet solide urbain, conformément à la Directive 2002/96/CE.
No tirar el aparato en los desechos urbanos, como exige la Directiva 2002/96/CE.
Dieses Gerät unterliegt der Richtlinie 2002/96/EG und darf nicht mit dem normalen Hausmüll entsorgt werden.
根据 EEC 指令 2002/96/CE，请不要将本设备作为城市垃圾处理。

This unit must be used for laboratory applications indoor only. The manufacturer declines all responsibility for any use of the unit that does not comply with these instructions. If the product is used in a not specified way by the manufacturer or with not specified accessories, product's safety may be compromised.

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This unit has been designed and manufactured in compliance with the following standards:

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Electrical equipment for laboratory use

UL 61010-1

General requirement - Canadian electrical code

CAN/CSA-C22.2 No.61010-1

VELP reserves the right to modify the characteristics of its products with the aim to constantly improving their quality.

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The values indicated on the rating plate of the instrument must correspond to those of the power supply.

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Fasten the unit to the support rod using the fixing block. Position the instrument with a distance from the wall of 30 cm (at least).

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The unit is fitted with two fuses (2xT5 A L 250 V), found in the socket on the back. To replace one or more, disconnect the mains cable and, using a screwdriver, lift up the small cover on the fuse box.

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Do not use with explosive or dangerous materials for which the equipment is not designed. The stirrer must not be used in explosive atmospheres, in bain-marie or to stir harmful liquids prior using protective measures according to the safety standards of the processed products and/or in force in the laboratories including personal protective equipment and the presence of an extraction hood which ensures at least 10-fold air change in accordance with the standards EN 14175 and DIN 12924.

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It is dangerous to run the motor free in air. Always fix the motor before turning the unit on.

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The solution may release toxic, dangerous or poisonous gases. Adequate safety measures must be taken, in accordance with the safety regulations in force, including the presence of hood and personal protective equipment (masks, gloves, goggles, etc.).

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CONTROLLER Digital is an innovative solution for control and data logging in process reactors, able to monitor the temperature in the reactor and the shaft rotation speed of brushless gear-motors for liquid volumes till 150L, even with viscosity changes. The motor is reversible.

Figure 1. CONTROLLER Digital

- A 7 segment display LED
- B Torque indicator bar
- C Clockwise rotation LED
- D Reverse key
- E Anticlockwise rotation LED
- F Speed control knob
- G Timer LED

- H Timer/Lock key
- I Lock LED
- J Main switch
- K USB door
- L Appliance inlet with fuse holder
- M Motor circular connector 6P *
- N Motor circular connector 4P **

*Circular connectors Model MIL-C-505 MS3102A-14S-6

**Circular connectors Model MIL-C-505 MS3102A -14S-2

2 EN - Installation

- Unpacking
 - Check the integrity of the unit after unpacking.
- The box includes
 - Controller Digital
 - Power supply cable
 - Instruction manual
- First installation
 - Place the unit on non-flammable surface
 - Fix the instrument to the support through the handle
 - Connect the instrument to one of the following brushless motors 4 poles, 24 V:
1/11HP (max 4.4 A) or 1/4HP (max 8,3 A) or 3/8HP (max 15 A)

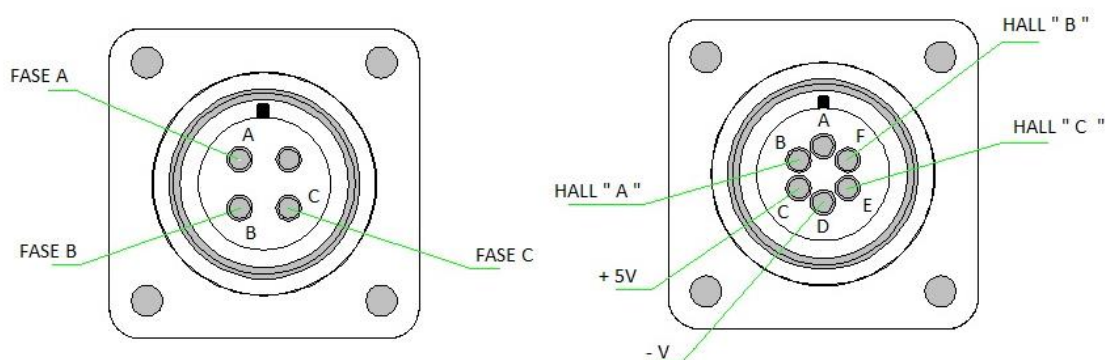
Figure 2. Connector side view

- Make sure that the rating value of the instrument corresponds to the one of the power supplies
- Ensure that the socket provided with grounding, is compliant with the current safety norms and easy to reach. Use only the cable provided with the instrument.
- Insert the power cable into the socket. and switch the instrument on.
- Select the motor type (see chapt.5).

NOTE: The cable can be substituted only by main cables with same features (T=70°C, connector C14).

3 EN - Working

Commissioning	➤ Switch on the instrument through the main switch (J) ➤ Display (A) shows software version, the last set point value and OFF (if "Mode" is set on Stop, see chapt.5)
Stirring	➤ Adjust speed set point by turning the speed control knob (F) ➤ Click the knob to start stirring ➤ The speed increases until set point achievement ➤ A microprocessor ensures constant speed even when the viscosity changes (counter-reaction) ➤ Switch off the stirring by clicking the knob
Timer	➤ Click Timer key (H). HH:MM appears on display ➤ Rotate speed control knob (F) to set the desired value. Click to confirm ➤ Once timer is set, timer LED (G) lights up ➤ The LED light is continuously until the instrument starts to stir. When the instrument stirs, the LED flashes. ➤ It's possible to visualize (or modify) countdown in each moment during the analysis by clicking Timer key (H) ➤ When countdown finishes, "End" is visualized on the display.

Reverse	➤ Set the stirring rotation direction by clicking Reverse key (D). ➤ The rotation direction set has its relative LED lighted on: clockwise (C) and anticlockwise (E)
Torque	➤ The torque indicator bar (B), made by four LEDs, shows the intensity of the torque applied ➤ CONTROLLER Digital model doesn't show a precise torque value as the Advance model
Lock	➤ Holding the Lock key (H) for 3 seconds it's possible to lock the control panel ➤ Unlock the control panel by holding key H for 3 seconds ➤ When the instrument is locked LED (I) lights up

4 EN - External Connections

USB	➤ for software upgrading, using a dedicated program to requested by to service@velp.it and to be installed in the PC
------------	--

5 EN - Menu

Hold the control knob (F) for 3 seconds to enter into the setting mode when stirring is off.

Turning the control knob (F) it is possible to pass from one of the following parameters to the next one. Click the knob to modify:

Parameter shown on display	Default value	Range	Description
mot	mot 1	mot 1 mot 2 mot 3 mot 4	Type motor: ➤ Mot 1: A00000415 ➤ Mot 2: A00000416 ➤ Mot 3: A00000417 ➤ Mot 4: A0000.....
mode	STOP	STOP - run	Set starting mode: ➤ Stop: when the instrument is switched on through the main switch, OFF is on display ➤ Run: when the instrument is switched on, it restarts to work with the last set points set
L INt	500	100 ÷ 1200 SPEED SELEcted Motor	Set full scale speed: ➤ It limits the maximum value of speed set point (from 100rpm, step 100rpm)
COun			Running time. Days (24H): ➤ Motor working time is displayed in hours until 9999 hours (around 416 days). After that, it's shown in days starting from 417 and a dot turns on as following, to differentiate days from hours ".417"
UPGr			Software upgrading: ➤ It allows to install a new software version
rSEt	n0	YES - n0	Reset: ➤ It allows to restore default setting parameters

Click the stirring control knob (F) to display the set value. Turn the knob to set the new value and click to confirm.

To leave settings mode wait 10 seconds without touching any knob or click the lock key.

6 EN - Error messages

When the display shows an error message, the stirring function stops automatically.

Error code	Cause
AL1	Motor doesn't start stirring
AL2	High internal motor temperature
AL3	Motor overload
AL4	High driver temperature
AL5	Safety relay fault

To remove the error message, switch the instrument off and then switch on it.

If alarm persists on the display, please contact VELP Scientifica technical service department. service@velp.it

7 EN - Maintenance

Maintenance	➤ No routine or extraordinary maintenance is necessary. ➤ Repairs must be carried out by authorized Velp personnel only. ➤ Instrument must be transported into its original packaging any indications present on the original packaging must be followed (e.g., palletized).
Cleaning	➤ Disconnect the unit from the power supply and use a cloth with a non-flammable non-aggressive detergent.

8 EN - Technical data

General features	Model	F201A0505
	Power supply	115-230V (+/-10%) – 50-60 Hz
	Dimensions (WxHxD)	90x180x170 mm (3.54x7.10x6.70 in)
	Weight	1,5 kg (3,30 lb)
	Power input	500 W
	Max power output	280 W (15A)
	Construction material (structure)	Stainless steel
	Working in continuous	Admitted
	Settable restart modality	Stop or work
	Noisiness	<< 60 dBa
	Environmental temperature admitted	+5...+40 °C
	Storage temperature admitted	-10...+60 °C
	Max humidity	80%
	Overvoltage category	II
	Pollution degree CEI EN61010-1	2
	Max altitude	2000 m
Stirring	Programmable speed range	35-500rpm (max speed depends to the motor selected)
	Speed selection	1 rpm step
	Stirring alarm	Motor fault
Torque	Max torque admitted	621 Ncm
Counters	Motor counter	Working hours

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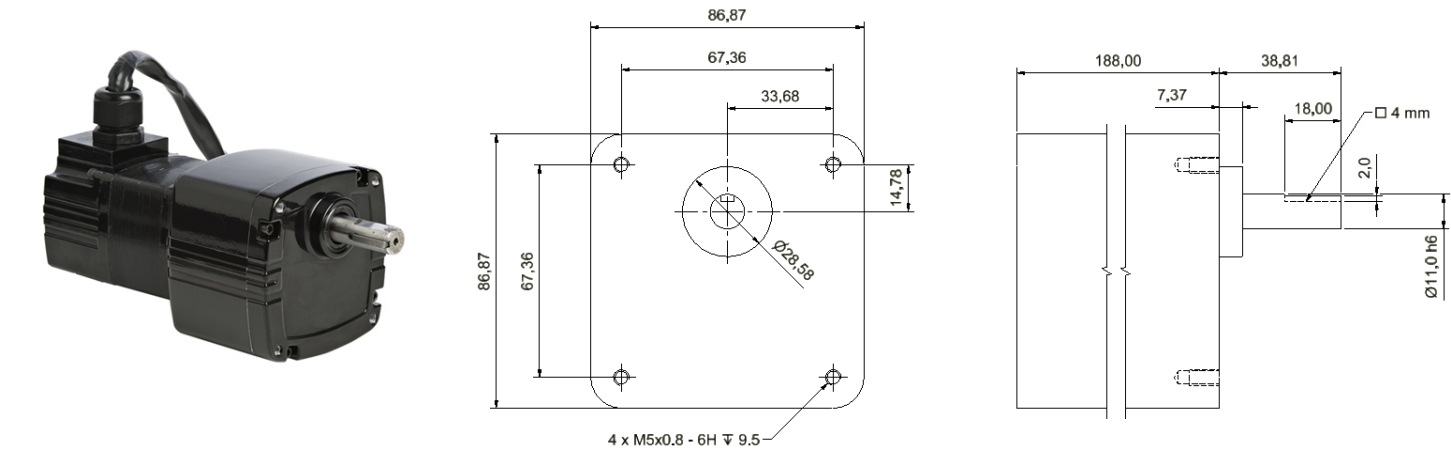
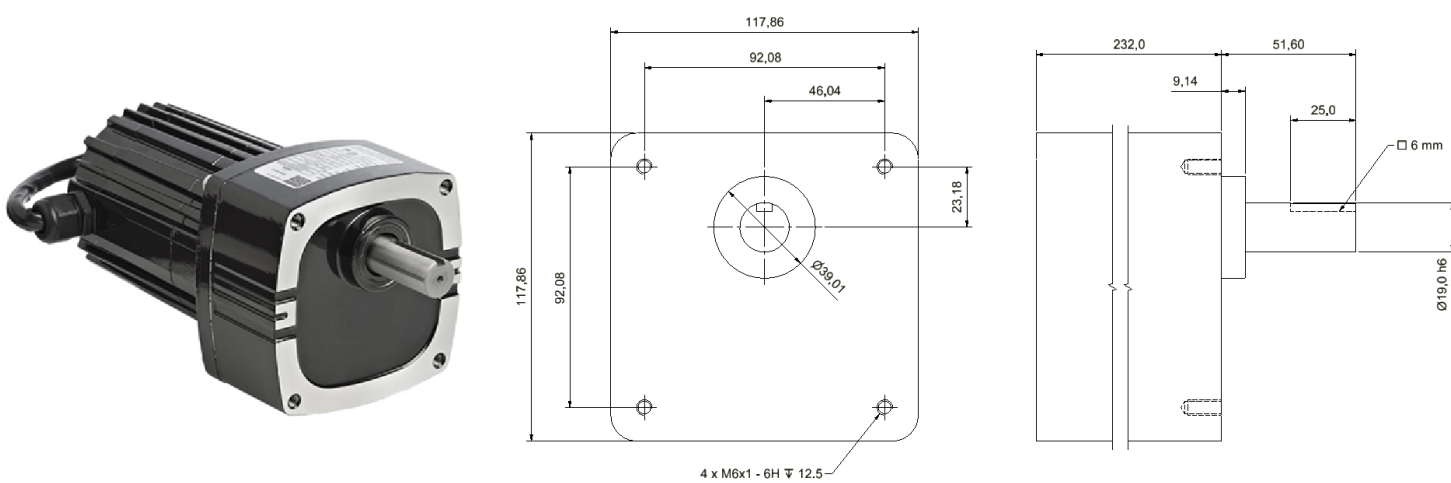
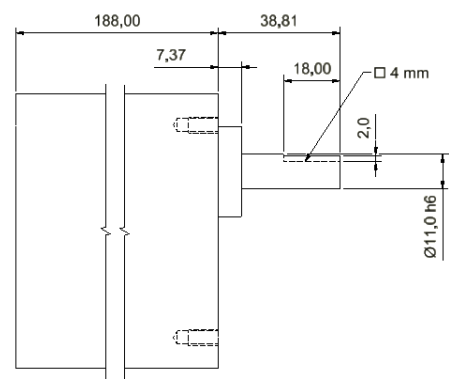
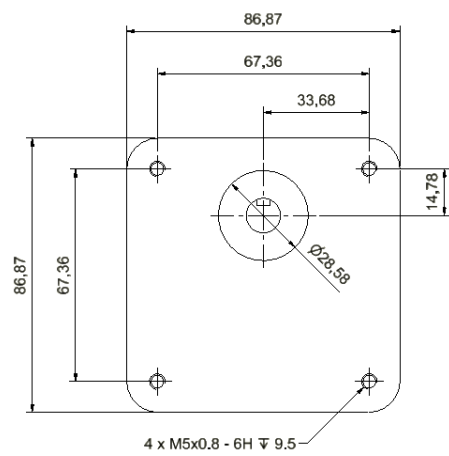
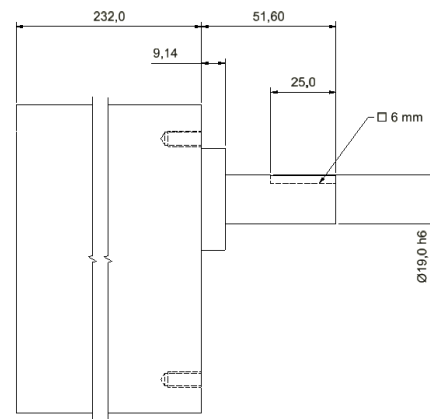
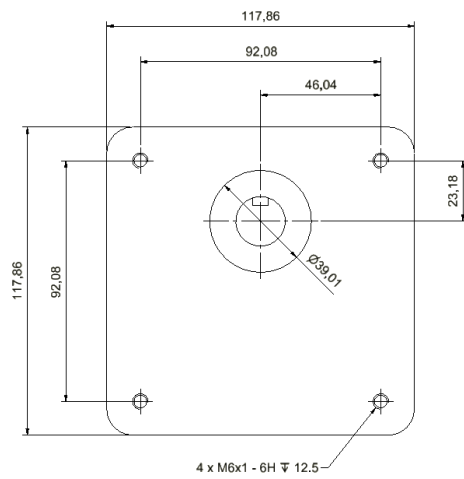
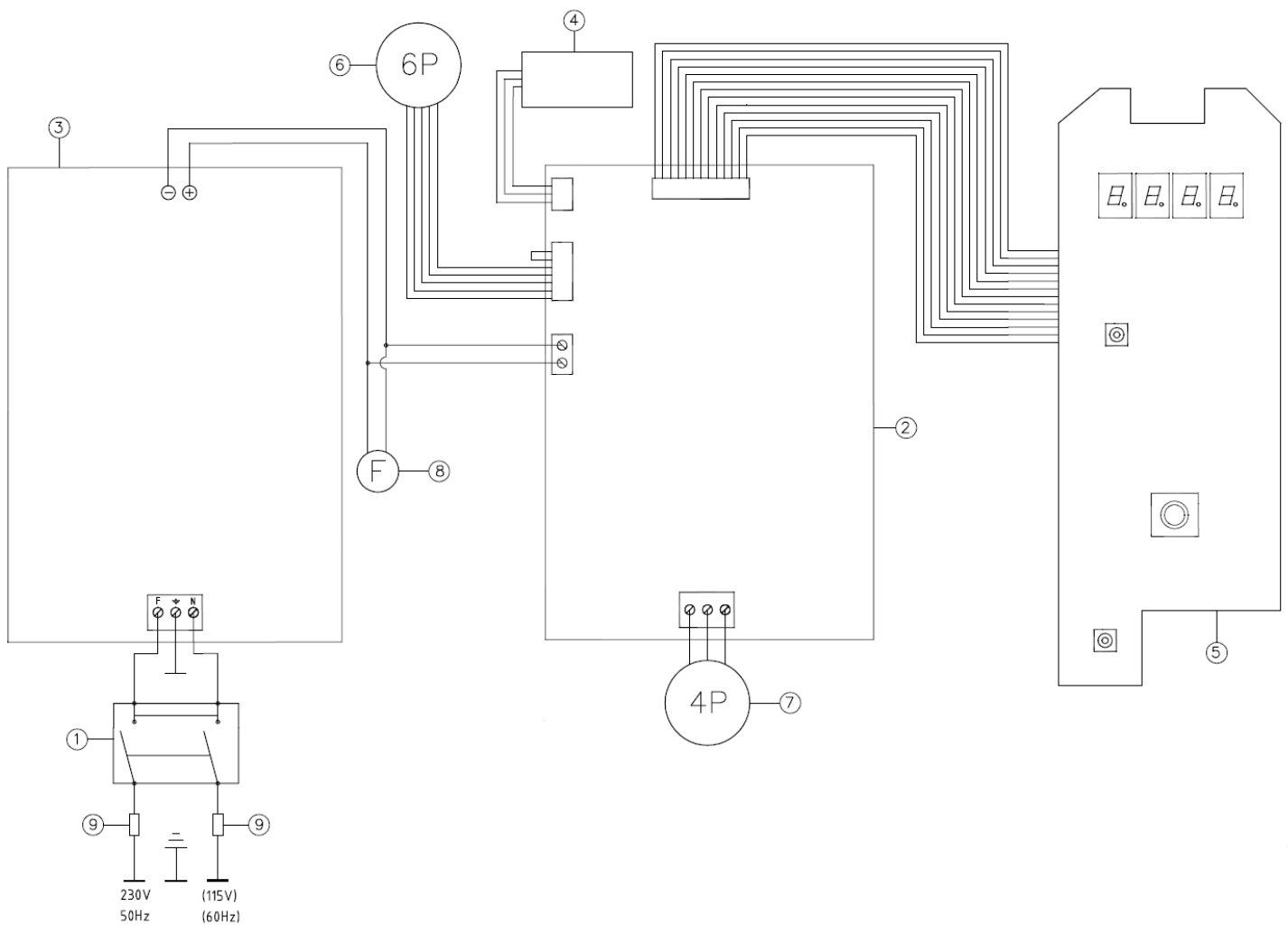
We, the manufacturer VELP Scientifica, under our responsibility declare that the product is manufactured in conformity with the following standards:

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EN 61010-1	Safety Requirements for Electrical Equipment for Measurement, Control, and Laboratory Use - Part 1: General Requirements.
EN 61010-2-051	Safety requirements for electrical equipment for measurement, control and laboratory use - Part 2-051: Particular requirements for laboratory equipment for mixing and stirring
EN 61326-1	Electrical equipment for measurement, control and laboratory use. EMC requirements. General requirements
2015/863/EU (RoHS III)	Restriction of the use of certain hazardous substances
2012/19/EU (WEEE)	On waste electrical and electronic equipment

and satisfies the essential requirements of the following directives:

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13 Declaration of conformity ^{UK}CA

We, the manufacturer VELP Scientifica, under our responsibility declare that the product is manufactured in conformity with the following regulations:

S.I. 2016/1101 Electrical Equipment (Safety) Regulations 2016

S.I. 2016/1091 Electromagnetic Compatibility Regulations 2016

according to the relevant designated standards:

EN 61010-1 Safety requirements for electrical equipment for measurement, control, and laboratory use
Part 1: General requirements

EN 61010-2-051 Particular requirements for laboratory equipment for mixing and stirring

EN 61326-1 Electrical equipment for measurement, control and laboratory use - EMC requirements -
Part 1: General requirements

and satisfies the essential requirements of regulations:

S.I. 2008/1597 Supply of Machinery (Safety) Regulations 2008

S.I. 2012/3032 Restriction of the Use of Certain Hazardous Substances in Electrical and Electronic
Equipment Regulations 2012

S.I. 2013/3113 Waste Electrical and Electronic Equipment Regulations 2013

plus modifications.

Thank you for having chosen VELP!

Established in 1983, VELP is today one of the world's leading manufacturer of analytical instruments and laboratory equipment that has made an impact on the world-wide market with Italian products renowned for innovation, design and premium connectivity. VELP works according to **ISO 9001**, **ISO14001** and **ISO 45001** Quality System Certification.

Our instruments are manufactured in Italy according to the IEC 1010-1 and CE regulation.

Our product lines:

Analytical instruments

Elemental Analyzers
Digestion Units
Distillation Units
Solvent Extractors
Fiber Analyzers
Dietary Fiber Analyzers
Oxidation Stability Reactor
Consumables

Laboratory Equipment

Magnetic Stirrers
Heating Magnetic Stirrers
Heating Plates
Overhead stirrers
Vortex mixers
Dispersers
COD Thermoreactors
BOD and Respirometers
Cooled Incubators
Flocculators
Overhead Shakers
Turbidimeter
Open Circulating Baths
Pumps

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