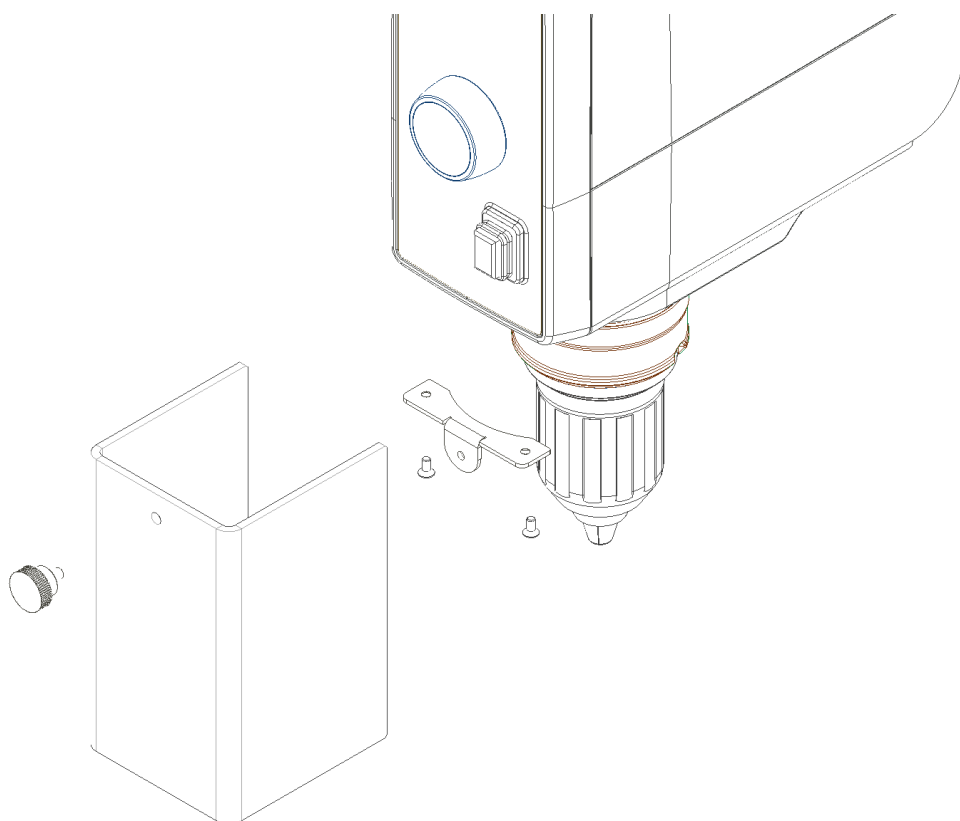
Instruction Manual
Manuale di istruzioni
Manuel d'instructions
Manual de instrucciones
Bedienungsanleitung

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Read the operating instructions carefully before use and follow the safety instructions.

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Keep the operating instructions in a place where it can be accessed by everyone.

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Safe operation of VELP Scientifica laboratory device with Stands and Double clamp is only guaranteed for combinations listed in the related device operating manual.

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The knob may come loose, resulting in risk to the user and trial set-up.

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Do not substitute knob or tighten it with a tool.

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Check that the Chuck guard is properly fixed before each use.

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Remove the Chuck guard only when the instrument is disconnected from the power supply.

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Use the chuck guard only with VELP Scientifica Laboratory devices.

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Do not use the chuck guard if damaged.

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Observe the technical specifications of this manual.

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Assembly / Assemblaggio / Assemblage / Montaje / Montage

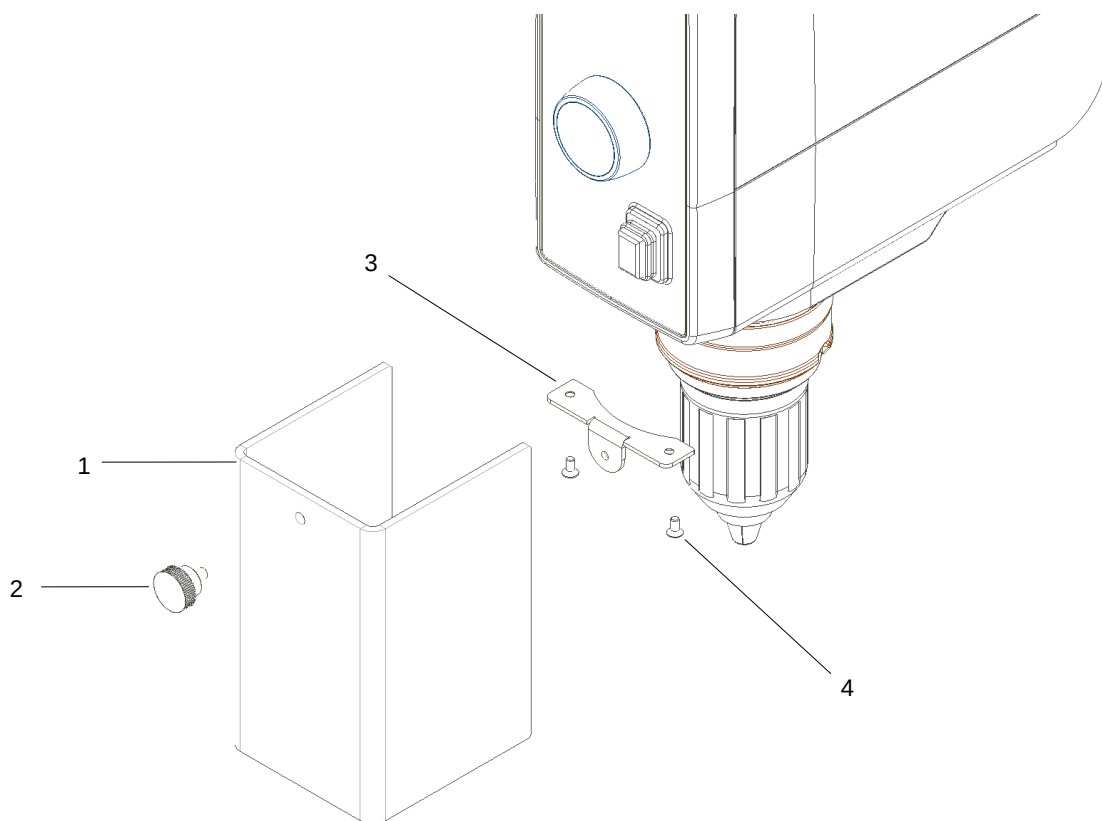
The assembly of the A00001312 stirring Chuck guard may be accomplished only with switched off equipment and with pulled power supply plug.

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Fix the support plate (Pos.3) using the two screws (Pos.4). Fix the chuck guard (Pos.1) using the knob (Pos.2).

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Thank you for having chosen VELP!

Established in 1983, VELP is today one of the world's leading manufacturer of analytical instruments and laboratory equipment that has made an impact on the world-wide market with Italian products renowned for innovation, design and premium connectivity. VELP works according to **ISO 9001**, **ISO 14001** and **OHSAS 18001** Quality System Certification.

Our instruments are manufactured in Italy according to the IEC 1010-1 and CE regulation.

Our product lines:

Analytical instruments

Elemental Analyzers
Digestion Units
Distillation Units
Solvent Extractors
Fiber Analyzers
Dietary Fiber Analyzers
Oxidation Stability Reactor
Consumables

Laboratory Equipment

Magnetic Stirrers
Heating Magnetic Stirrers
Heating Plates
Overhead stirrers
Vortex mixers
Homogenizers
COD Thermoreactors
BOD and Respirometers
Cooled Incubators
Flocculators
Overhead Shakers
Turbidimeter
Radiation Detector
Open Circulating Baths
Pumps

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