Instruction Manual
Manuale di istruzioni
Manuel d'instructions
Manual de instrucciones
Bedienungsanleitung
指导手册

ESP Ultraflat Magnetic Stirrer

F206A0179

**General Information / Informazioni Generali / Informations Générales / Información General /
Allgemeine Hinweise / 一般信息**

Before using the unit, please read the following instruction manual carefully.

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Do not dispose of this equipment as urban waste, in accordance with EEC directive 2002/96/CE.

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This unit must be used for laboratory applications indoor only. The manufacturer declines all responsibility for any use of the unit that does not comply with these instructions. If the product is used in a not specified way by the manufacturer or with non specified accessories, product's safety may be compromised.

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This unit has been designed and manufactured in compliance with the following standards:

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**IEC/EN 61010-1
IEC/EN 61010-2-051**

Electrical equipment for laboratory use

UL 61010-1

General requirement - Canadian electrical code

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Do not use with explosive and dangerous materials for which the equipment is not designed. The stirrer must not be used in explosive atmospheres.

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It is responsibility of the user appropriately decontaminate the instrument in case of dangerous substances fall on or in it.

It is also responsibility of the user to use safety substances for cleaning or decontaminating, which do not react with internal parts of the instrument or with the material contained in it. In case of doubts on the compatibility of a cleaning solution, contact the manufacturer or local distributor.

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Do not use with explosive and dangerous materials for which the equipment is not designed. The stirrer must not be used in explosive atmospheres, in bain-marie and to stir combustible liquids that have a low combustion temperature. The minimum fire point of flammable solution is 200 °C. Only small amounts (< 50 ml) of flammable liquid can be used with the device.

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In case of black-out, once the power is back the device will restart automatically.

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The magnetic stirrer without motor ESP is used for chemical, biotechnological, pharmaceutical, microbiological, medical applications, for example for growing microorganisms, dissolving nutrients and solids, avoiding settling of suspended matter or during titrations. The power of magnetic field allows to stir liquids in vessels with different shapes up to 5 liters.

Ultra-thin magnetic stirrer without moving mechanical components, not requiring maintenance. The stirring system is composed by coils inducing a rotating magnetic field which is controlled by a microprocessor; speed range 50 ÷ 1.100 rpm. By the rotation of the knob completely clockwise the maximum stirring speed (1.100 rpm) will be obtained. By a counter clockwise rotation a lower of speed will be obtained up to a complete stop. Minimum speed is 50 rpm.

A slow start allows an optimal progression of stirring speed. The modern and ergonomic structure is fabricated from materials with an high resistance.

2. Assembly and installation

Upon receipt and after having removed the packaging, please check the integrity of the instrument. The box includes:

- Magnetic stirrer ESP
- Plug for power supply
- Switching VELP 100-240V/12V without plug
- Instruction manual

2.1 Electrical connection

After having unpacked the instrument, place the unit on the laboratory bench.

Before connecting the instrument to the power supply, make sure that the values on the rating plate correspond to those of the power supply. Connect the unit to the power supply using the transformer supplied.

Ensure that the socket and the relative cut-off device conform to current safety norms and easy to reach.

2.2 Start-up

Rotate the speed knob completely to the left. Place the flasks containing the sample and a suitable magnetic stirring bar on the stirring plate. Then, set the speed by turning the speed knob.

3. Operating controls

It is possible to select speeds between 50 and 1100 rpm by rotating the knob.

4. Maintenance

No routine or extraordinary maintenance is necessary apart from periodically cleaning the unit as described in this manual. In compliance with the product guarantee law, repairs to our units must be carried out in our factory, unless previously agreed otherwise with local distributors. The instrument must be transported in its original packaging and any indications present on the original packaging must be followed (e.g. palletized).

It is the responsibility of the user, to properly decontaminate the unit in case of hazardous substances remaining on the surface or interior of the device. If in doubt about the compatibility of a cleaning or decontamination product, contact the manufacturer or distributor.

4.1 Cleaning

Disconnect the unit from the power supply and use a cloth dampened with an non-flammable non-aggressive detergent.

5. Technical data

Admitted power supply	0.5 A / 100-240 V / 50-60 Hz (+/-10%)
Power	12 Vdc 6 W
Dimensions (WxHxD)	160x35x230 mm
Weight (with plug)	1 kg
Speed range	50 ÷ 1100 rpm
Stirring capacity	5 L
Operating mode	Continuous
Construction material	Technopolymer
Environmental ambient temperature	5 - 40 °C (41 - 104°F)
Storage temperature range	-10... + 60 °C (14 - 140°F)
Pollution degree	2
Overvoltage category	II
Sound level	35 dBa
Max humidity	80%
Altitude	Up to 2000 m.

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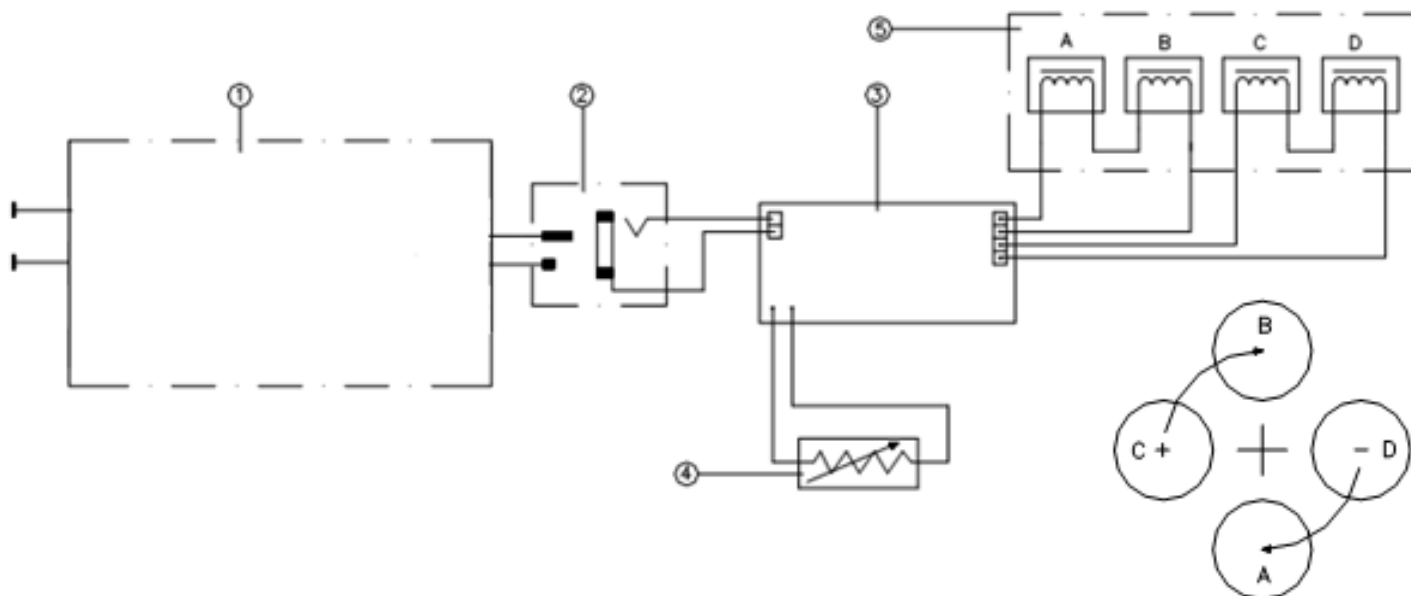
We, the manufacturer VELP Scientifica, under our responsibility declare that the product is manufactured in conformity with the following standards:

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2015/863/EU (RoHS III)

2012/19/EU (WEEE)

and satisfies the essential requirements of the following directives:

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9. Declaration of conformity ^{UK}CA

We, the manufacturer VELP Scientifica, under our responsibility declare that the product is manufactured in conformity with the following regulations:

S.I. 2016/1101 Electrical Equipment (Safety) Regulations 2016

S.I. 2016/1091 Electromagnetic Compatibility Regulations 2016

according to the relevant designated standards:

EN 61010-1 Safety requirements for electrical equipment for measurement, control, and laboratory use
Part 1: General requirements

EN 61010-2-051 Particular requirements for laboratory equipment for mixing and stirring

EN 61326-1 Electrical equipment for measurement, control and laboratory use - EMC requirements -
Part 1: General requirements

and satisfies the essential requirements of regulations:

S.I. 2008/1597 Supply of Machinery (Safety) Regulations 2008

S.I. 2012/3032 Restriction of the Use of Certain Hazardous Substances in Electrical and Electronic
Equipment Regulations 2012

S.I. 2013/3113 Waste Electrical and Electronic Equipment Regulations 2013

plus modifications.

Thank you for having chosen VELP!

Established in 1983, VELP is today one of the world's leading manufacturer of analytical instruments and laboratory equipment that has made an impact on the world-wide market with Italian products renowned for innovation, design and premium connectivity. VELP works according to **ISO 9001**, **ISO14001** and **ISO 45001** Quality System Certification.

Our instruments are manufactured in Italy according to the IEC 1010-1 and CE regulation.

Our product lines:

Analytical instruments

Elemental Analyzers
Digestion Units
Distillation Units
Solvent Extractors
Fiber Analyzers
Dietary Fiber Analyzers
Oxidation Stability Reactor
Consumables

Laboratory Equipment

Magnetic Stirrers
Heating Magnetic Stirrers
Heating Plates
Overhead stirrers
Vortex mixers
Dispersers
COD Thermoreactors
BOD and Respirometers
Cooled Incubators
Flocculators
Overhead Shakers
Turbidimeter
Open Circulating Baths
Pumps

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VELP Scientifica Srl