Instruction Manual
Manuale di istruzioni
Manuel d'instructions
Manual de instrucciones
Bedienungsanleitung
指导手册

FC4S - FC6S Flocculators

F105A0111, F105A0112

**General Information / Informazioni Generali / Informations Générales / Información General /
Allgemeine Hinweise / 一般信息**

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This unit has been designed and manufactured in compliance with the following standards:

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Safety requirements for electrical equipment for measurement, control and for laboratory use

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IEC/EN 61010-1

Electrical equipment for laboratory use

UL 61010-1

General requirement - Canadian electrical code

CAN/CSA-C22.2 No.61010-1

VELP reserves the right to modify the characteristics of its products with the aim to constantly improving their quality.

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During chemical-physical treatments of waters aimed to coagulate and precipitate pollutants, chemical coagulants are evaluated and their dosing optimized by jar tests. In relation to the variable composition of wastewaters the chemical dosing and the test frequency are to be adapted to local conditions. The jar test equipment manufactured by VELP is characterized by an electronic speed control with an independent speed settable for each place in order to optimize the result of settling so to reduce chemical consumption. FC4S and FC6S allow to pre-select 10 different speeds for each stirring point using the relative selectors located on the front panel (10-15-30-45-60-90-120-150-200-300). The steel structure epoxy painted is studied and tested in order to give the instrument an excellent resistance against the attack of chemical agents and corrosion in general. The stainless steel stirring rods are adjustable in height by a self-blocking chuck. The examined samples can be lighted by a back led lamp controlled by a switch located on front panel.

2. Assembly and installation

Upon receipt and after having removed the packaging, please check the integrity of the instrument. The box includes:

- Flocculator
- Plug for power supply
- Switching VELP 100-240V/12V without plug
- Instruction manual

2.1 Electrical connection

Before connecting the instrument to the power supply, make sure that the values on the rating plate correspond to those of the power supply. Connect the unit to the power supply using the switching.

Ensure that the socket and the relative cut-off device conform to current safety norms and easy to reach.

2.2 Start-up

Rotate the speed knob completely to the left. Place the beakers containing the sample and then, set the speed.

3. Operating controls

It's possible to select the desired stirrer speed by the knob "Stirrer rpm" on the frontal panel. It's possible to choose the desired values: 10-15-30-60-90-120-150-200-300 rpm by using the information written on the analogical scale around the knob.

At the end of the work cycle set the stirring speed knob on "0" rpm and turn off the light switch. If the instrument is not used for a long period it's better to disconnect the external power supply from the plug.

4. Maintenance

No routine or extraordinary maintenance is necessary apart from periodically cleaning of the unit as described in this manual. In compliance with the product guarantee law, repairs to our units must be carried out in our factory, unless previously agreed otherwise with local distributors. The instrument must be transported in its original packaging and any indications present on the original packaging must be followed (e.g. palletized).

It is the responsibility of the user, to properly decontaminate the unit in case of hazardous substances remaining on the surface or interior of the device. If in doubt about the compatibility of a cleaning or decontamination product, contact the manufacturer or distributor.

4.1 Cleaning

Disconnect the unit from the power supply and use a cloth dampened with an non-inflammable non-aggressive detergent.

5. Technical data

Power supply	100-240 V / 50-60 Hz
Power	18 W (FC4S) 23 W (FC6S)
Dimensions (WxHxD)	655x404x296 mm (FC4S) 935x404x296 mm (FC6S)
Weight (with plug)	12.5 Kg (FC4S) 18 Kg (FC6S)
Speed range	10-15-30-45-60-90-120-150-200-300 rpm
Construction material	Epoxy painted metal structure
Environmental ambient temperature	5 - 40 °C (41 - 104°F)
Storage temperature range	-10... + 60 °C (14 - 140°F)
Max humidity	80%
Pollution degree	2

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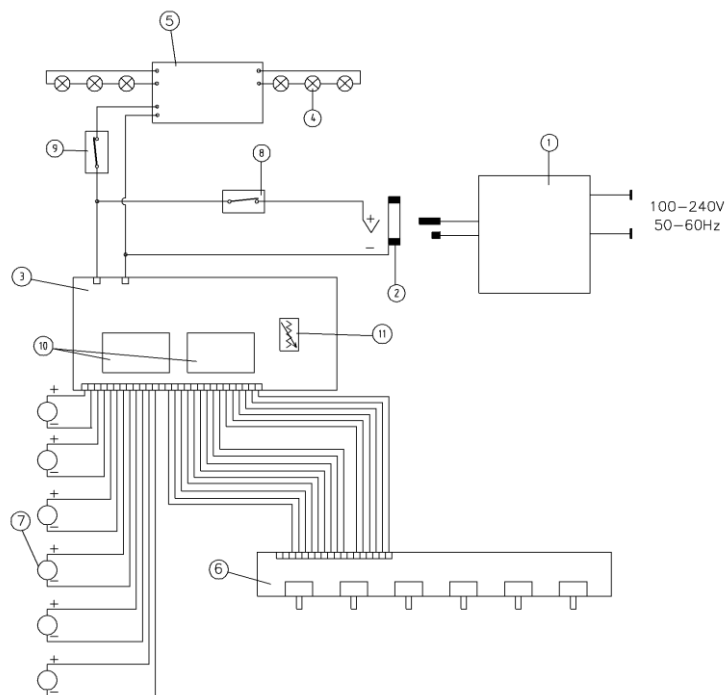
We, the manufacturer VELP Scientifica, under our responsibility declare that the product is manufactured in conformity with the following standards:

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2015/863/EU (RoHS III)

2012/19/EU (WEEE)

and satisfies the essential requirements of the following directives:

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9. Declaration of conformity ^{UK}CA

We, the manufacturer VELP Scientifica, under our responsibility declare that the product is manufactured in conformity with the following regulations:

S.I. 2016/1101 Electrical Equipment (Safety) Regulations 2016

S.I. 2016/1091 Electromagnetic Compatibility Regulations 2016

according to the relevant designated standards:

EN 61010-1 Safety requirements for electrical equipment for measurement, control, and laboratory use
Part 1: General requirements

EN 61010-2-051 Particular requirements for laboratory equipment for mixing and stirring

EN 61326-1 Electrical equipment for measurement, control and laboratory use - EMC requirements -
Part 1: General requirements

and satisfies the essential requirements of regulations:

S.I. 2008/1597 Supply of Machinery (Safety) Regulations 2008

S.I. 2012/3032 Restriction of the Use of Certain Hazardous Substances in Electrical and Electronic
Equipment Regulations 2012

S.I. 2013/3113 Waste Electrical and Electronic Equipment Regulations 2013

plus modifications.

Thank you for having chosen VELP!

Established in 1983, VELP is today one of the world's leading manufacturer of analytical instruments and laboratory equipment that has made an impact on the world-wide market with Italian products renowned for innovation, design and premium connectivity. VELP works according to **ISO 9001**, **ISO14001** and **OHSAS 18001** Quality System Certification.

Our instruments are manufactured in Italy according to the IEC 1010-1 and CE regulation.

Our product lines:

Analytical instruments

Elemental Analyzers
Digestion Units
Distillation Units
Solvent Extractors
Fiber Analyzers
Dietary Fiber Analyzers
Oxidation Stability Reactor
Consumables

Laboratory Equipment

Magnetic Stirrers
Heating Magnetic Stirrers
Heating Plates
Overhead stirrers
Vortex mixers
Homogenizers
COD Thermoreactors
BOD and Respirometers
Cooled Incubators
Flocculators
Overhead Shakers
Turbidimeter
Radiation Detector
Open Circulating Baths
Pumps

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VELP Scientifica Srl