**Instruction Manual**  
**Manuale di istruzioni**  
**Manuel d'instructions**  
**Manual de instrucciones**  
**Bedienungsanleitung**  
**指导手册**

**FP4 Portable Flocculator**

F105A0117

**General Information / Informazioni Generali / Informations Générales / Información General /  
Allgemeine Hinweise / 一般信息**

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**This unit has been designed and manufactured in compliance with the following standards:**

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Safety requirements for electrical equipment for measurement, control and for laboratory use

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**IEC/EN 61010-1**

Electrical equipment for laboratory use

**UL 61010-1**

General requirement - Canadian electrical code

**CAN/CSA-C22.2 No.61010-1**

VELP reserves the right to modify the characteristics of its products with the aim to constantly improving their quality.

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Flocculators are studied both to optimize the dosing of coagulants in order to separate the pollutions in water treatment plants, thanks to the laboratory valuation results obtained with the so-called Jar Test.

It is possible to set the stirring speed (1) in analogical way and the stirring time (2) by knobs placed on the frontal panel. The examined samples can be lighted by a back fluorescent lamp controlled by a switch located on front panel. FP4 was developed for optimum use in the field; the instrument can run both connected to a normal battery and a car's cigarette lighter.

The steel structure epoxy painted is studied and tested in order to give to the instrument an excellent resistance to the attack of chemical agents and to corrosion in general. The external power supply allows the instrument to be connected to all frequencies and voltages in the world (100÷240V / 50÷60Hz).

It's possible to select the stirring speed (20-40-50-100-200 rpm), the operating time and the lighting of central column by the frontal panel. The stirring speed is the same for each position. It is possible to set the time of running from 0 to 30 minutes and the central column start up. FP4 is characterized by a non-skid base so as to keep the beakers stable during operation.

## 2. Assembly and installation

Upon receipt and after having removed the packaging, please check the integrity of the instrument. The box includes:

- FP4 Portable Flocculator
- Switching VELP 100-240V/12V without plug
- Plug for power supply
- Instruction manual

### 2.1 Electrical connection

Before connecting the instrument to the power supply, make sure that the values on the rating plate correspond to those of the power supply. Connect the unit to the power supply using the transformer supplied.

Ensure that the socket and the relative cut-off device conform to current safety norms and easy to reach.

### 2.2 Start-up

Rotate the speed knob completely to the left. Place the beakers containing the sample and then, set the speed and time by turning the knobs.

### 2.3 Information about construction materials

Structure	Aluminum	Frontal panel label	PET	Feet	MQ/MVQ
-----------	----------	---------------------	-----	------	--------

## 3. Operating controls

It's possible to select the desired stirrer speed by the knob "Stirrer rpm" on the frontal panel. It's possible to choose the desired values: 0-20-40-50-100-200 rpm by using the information written on the analogical scale around the knob.

The start-up and the selection of the running time with automatic turn off are made by the relative timer knob:

by turning un-clockwise the timer knob on the manual modality the instrument runs in continuous way;

by turning clockwise the timer knob it's possible to select the time of running up to 30 minutes; the stirring automatically stops as the time finishes.

At the end of the work cycle set the stirring speed knob on "0" rpm, the timer knob on "OFF" and turn off the light of central column. If the instrument is not used for a long period it's better to disconnect the external power supply from the plug.

## 4. Maintenance

No routine or extraordinary maintenance is necessary apart from periodically cleaning the unit as described in this manual. In compliance with the product guarantee law, repairs to our units must be carried out in our factory, unless previously agreed otherwise with local distributors. The instrument must be transported in its original packaging and any indications present on the original packaging must be followed (e.g. palletized).

It is the responsibility of the user, to properly decontaminate the unit in case of hazardous substances remaining on the surface or interior of the device. If in doubt about the compatibility of a cleaning or decontamination product, contact the manufacturer or distributor.

### 4.1 Cleaning

Disconnect the unit from the power supply and use a cloth dampened with an non-inflammable non-aggressive detergent.

## 5. Technical data

External power supply - inlet	AC 100÷240V ; 50-60 Hz ; 0.5A
External power supply - outlet	DC 12V ; 1.25A
Power	6 W
Dimensions (WxHxD)	250x305x250mm
Weight (with plug)	3.8 Kg
Speed range	0-20-40-50-100-200 rpm
Construction material	Epoxy painted metal structure
Environmental ambient temperature	5 - 40 °C (41 - 104°F)
Storage temperature range	-10... + 60 °C (14 - 140°F)
Max humidity	80%
Pollution degree	2

## 6. Accessories / Spare parts

A00001000	Plastic Becher 1000 ml
A00001001	Glass Becher 1000 ml
A00001005	Carrying case

10000694	Black knob
10000235	Foot

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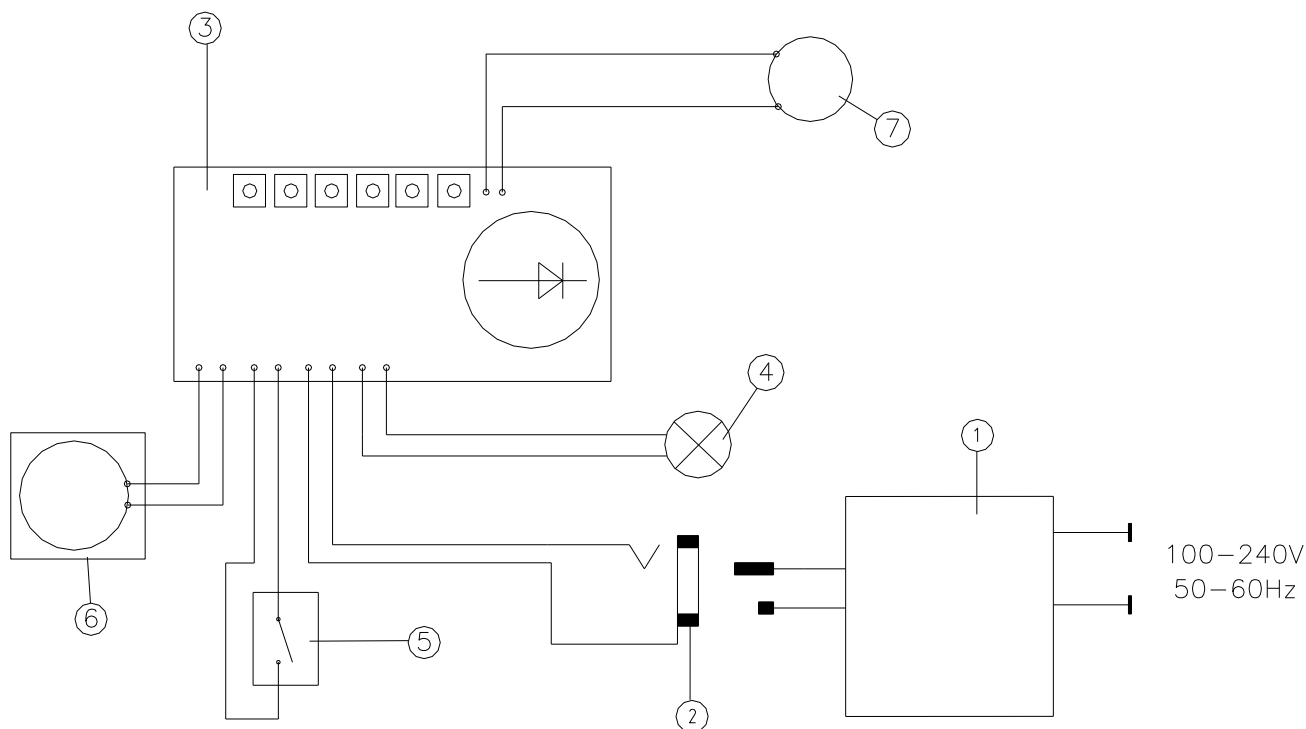
We, the manufacturer VELP Scientifica, under our responsibility declare that the product is manufactured in conformity with the following standards:

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EN 61010-2-051

EN 61326-1

and satisfies the essential requirements of the following directives:

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## 9. Declaration of conformity <sup>UK</sup>CA

We, the manufacturer VELP Scientifica, under our responsibility declare that the product is manufactured in conformity with the following regulations:

S.I. 2016/1101 Electrical Equipment (Safety) Regulations 2016

S.I. 2016/1091 Electromagnetic Compatibility Regulations 2016

according to the relevant designated standards:

EN 61010-1 Safety requirements for electrical equipment for measurement, control, and laboratory use  
Part 1: General requirements

EN 61010-2-051 Particular requirements for laboratory equipment for mixing and stirring

EN 61326-1 Electrical equipment for measurement, control and laboratory use - EMC requirements -  
Part 1: General requirements

and satisfies the essential requirements of regulations:

S.I. 2008/1597 Supply of Machinery (Safety) Regulations 2008

S.I. 2012/3032 Restriction of the Use of Certain Hazardous Substances in Electrical and Electronic  
Equipment Regulations 2012

S.I. 2013/3113 Waste Electrical and Electronic Equipment Regulations 2013

plus modifications.

## Thank you for having chosen VELP!

Established in 1983, VELP is today one of the world's leading manufacturer of analytical instruments and laboratory equipment that has made an impact on the world-wide market with Italian products renowned for innovation, design and premium connectivity. VELP works according to **ISO 9001**, **ISO14001** and **OHSAS 18001** Quality System Certification.

Our instruments are manufactured in Italy according to the IEC 1010-1 and CE regulation.

Our product lines:

### Analytical instruments

Elemental Analyzers  
Digestion Units  
Distillation Units  
Solvent Extractors  
Fiber Analyzers  
Dietary Fiber Analyzers  
Oxidation Stability Reactor  
Consumables

### Laboratory Equipment

Magnetic Stirrers  
Heating Magnetic Stirrers  
Heating Plates  
Overhead stirrers  
Vortex mixers  
Homogenizers  
COD Thermoreactors  
BOD and Respirometers  
Cooled Incubators  
Flocculators  
Overhead Shakers  
Turbidimeter  
Radiation Detector  
Open Circulating Baths  
Pumps

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VELP Scientifica Srl