Instruction Manual
Manuale di istruzioni
Manuel d'instructions
Manual de instrucciones
Bedienungsanleitung
指导手册

FTC 120 Cooled Incubator

F10300305

General Information / Informazioni Generali / Informations Générales / Información General / Allgemeine Hinweise / 一般信息

Before using the unit, please read the following instruction manual carefully.
Prima dell'utilizzo dello strumento si raccomanda di leggere attentamente il seguente manuale operativo.
Avant d'utiliser l'instrument, il est recommandé de lire attentivement le présent manuel d'instructions.
Antes de utilizar el instrumento, le recomendamos que lea con atención el siguiente manual de instrucciones.
Bitte lesen Sie vor Inbetriebnahme des Geräts diese Bedienungsanleitung sorgfältig durch.
在使用本装置之前，请仔细阅读以下使用说明书。

There is a small quantity of isobutane gas (R600a) in the refrigerating system.
Lo strumento contiene una piccola quantità di gas isobutano (R600a) nel circuito refrigerante.
Il y a une petite quantité de gaz isobutane (R600a) dans le système de réfrigération.
Hay una pequeña cantidad de gas isobutano (R600a) en el sistema de refrigeración.
In der Kälteanlage befindet sich eine geringe Menge Isobutangas (R600a).
制冷系统中存在少量的异丁烷气体 (R600a)。

Do not dispose of this equipment as urban waste, in accordance with EEC directive 2012/19/UE.
Non smaltire l'apparecchiatura come rifiuto urbano, secondo quanto previsto dalla Direttiva 2012/19/UE.
Ne pas recycler l'appareil comme déchet solide urbain, conformément à la Directive 2012/19/UE.
No tirar el equipo en los residuos urbanos, como exige la Directiva 2012/19/UE.
Dieses Gerät unterliegt der Richtlinie 2012/19/UE und darf nicht mit dem normalen Hausmüll entsorgt werden.
根据 EEC 指令 2012/19/UE，不要将本设备作为城市垃圾处理。

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It is responsibility of the user appropriately decontaminate the instrument in case of dangerous substances fall on or in it accordingly to the safety datasheet of substances used and to the current laboratories' safety standards. It is not possible to decontaminate the product under steam.

It is also responsibility of the user to use substances for cleaning or decontaminating which do not react with internal parts of the instrument or with the material contained in it. In case of doubts on the compatibility of a cleaning solution, contact the manufacturer or local distributor.

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Be aware that liquids don't get in touch with power supply or with internal electronic components.

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Verify that the power supply is connected to a power plug (easily accessible) with the requirements stated on the label of the instrument.

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Use only power cables with grounding.

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Do not use the instrument after recognition of malfunctioning. Contact the nearest service center.

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The personal protective equipment must be compatible with the risks arising from the material being processed

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Do not place explosive substances with flammable propellants inside the instrument

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Be careful and do not damage the refrigerant circuit at the back of the instrument

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The inside of the instrument is equipped with 230V - 50Hz electrical power sockets to power the Velp Scientifica devices for BOD analysis. Do not switch on other electrical appliances inside the instrument unless approved by Velp Scientifica.

Do not connect other appliances with supply voltages other than 230V and with absorption greater than 3A.

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When you dispose the instrument, do so only in an authorized waste disposal center and do not expose it to flame.

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The instrument contains a small amount of isobutylene refrigerant (R600a) in the refrigerant circuit, which is a natural gas with high environmental compatibility but flammable. Ensure that each component of the refrigerant circuit is not damaged during transport or

installation. Refrigerant R600a is flammable. In case of breakage of the refrigerant circuit, avoid open flames, sparks and ventilate the room in which the appliance is located.

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Do not place hot objects near the plastic parts of the instrument.

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The instrument is heavy. Be careful when moving it.

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Do not place objects on the internal temperature probe positioned in the middle of the left internal wall of the instrument.

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The instrument is equipped with a forced internal circulation of air, so make sure that the products positioned inside do not obstruct this circulation or are kept detached from the rear and front internal walls.

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The internal shelves are grid in order to facilitate the forced circulation of air, do not replace them with others of the closed type in order not to compromise the homogeneity of internal temperature.

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For the electrical connection, carefully follow the instructions provided in the following manual. It is advisable to wait at least four hours before connecting the appliance to allow the liquid refrigerant gas inside the refrigerant circuit to flow into the compressor.

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Adequate air circulation should be around the appliance, failure leads to abnormal overheating which can cause performance loss.

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If possible, the back of the product must be against a wall to avoid touching hot parts (compressor and condenser).

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The instrument must not be placed near radiators or heat sources

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Never cover the condenser fins on the back of the instrument.

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The unit offers excellent thermal stability and homogeneity thanks to continuous forced ventilation and electronic temperature control.

The cooled incubator is designed to run at a constant temperature of 20 °C.

The internal temperature is measured by the internal probe and it is shown on the display.

The unit is equipped with two internal sockets, controlled by a switch on the frontal panel, to power instruments placed inside the incubator.

In accordance with current regulations, VELP incubators are CFC-free (free from chloro fluoro carbons).

2. Installation

- Unpacking
 - Check the integrity of the unit after unpacking
- The box includes
 - FTC 120 Cooled Incubator
 - Instruction manual
 - Power supply cable
- Installation
 - Make sure that the technical specification of the unit and the power supply line are the same
 - Connect the instrument to an easily accessible socket (compliant with the current safety norms), using only the provided power supply cable

NOTE: do not install the equipment near sources of heat (ovens, etc.). The unit should be installed in a well-ventilated area, avoid niches or small rooms. To ensure good air circulation within the unit, do not place objects against the inside back panel.

NOTE: if the instrument has been transported or stored horizontally, leave it in the upright position for at least 24 hours before switching it on. This will allow the cooling circuit to settle. If this procedure is not carried out, the refrigerating system could be damaged. This damage is not covered by warranty.

3. Operation

Turn the unit on using the main switch on the frontal panel.

Automatically, the unit starts the thermoregulation to achieve and maintain a constant temperature of 20 °C.

The internal temperature is measured by the internal probe and it is shown on the display.

Use the dedicated switch on the frontal panel to control the internal sockets.

Error messages

When the display shows an error message, the instrument's functions are stopped automatically.

A1	Internal temperature too low
A2	Internal temperature too high
A3	Internal probe broken

To remove the error message, restart the instrument.

If the alarm persists on the display, please contact VELP Scientifica's technical service department.

4. Maintenance

No routine or extraordinary maintenance is necessary except periodically cleaning.

Cleaning

Disconnect the unit from the power supply and use a cloth dampened with a non-flammable non-aggressive detergent.

Repair

Repairs must be carried out by authorized Velp personnel only.

The instrument must be transported using the original shockproof packaging when freight forwarders, couriers or other are taking care of the unit transport. Follow the indications on the original packaging (e.g. place the instrument on a pallet).

It is the responsibility of the user, to properly decontaminate the unit in case of hazardous substances remaining on the surface or interior of the device. If in doubt about the compatibility of a cleaning or decontamination product, contact the manufacturer or distributor.

5. Technical data

Power supply	230 V – 50/60 Hz
Usable internal volume	109 l
Power	120 W
Dimensions (WxHxD)	540x912x550 mm (21.3x35.9x21.7 in)
Weight	36 Kg (79.4 lbs)
Internal temp. display	2 digits display
Temperature setting	None
Setting range temperature	Fixed at 20°C
Temperature probe	Thermistor
Fuse diam. 5x20mm	5 A
Display resolution	1 °C
Sample temperature stability	± 0.5 °C
Sample homogeneity temperature	± 0.5 °C
Max. altitude of installation	4000 m (13123 ft) above sea level
Cooling regulator type	Electronic ON-OFF
Max humidity admitted	80%
Storage temp. range	-10 ÷ +60 °C (14 ÷ 140 °F)
Environmental temp. range	5 ÷ 40 °C (41 ÷ 104 °F)
Max load (Kg) for each shelf	10 kg (homogeneously)

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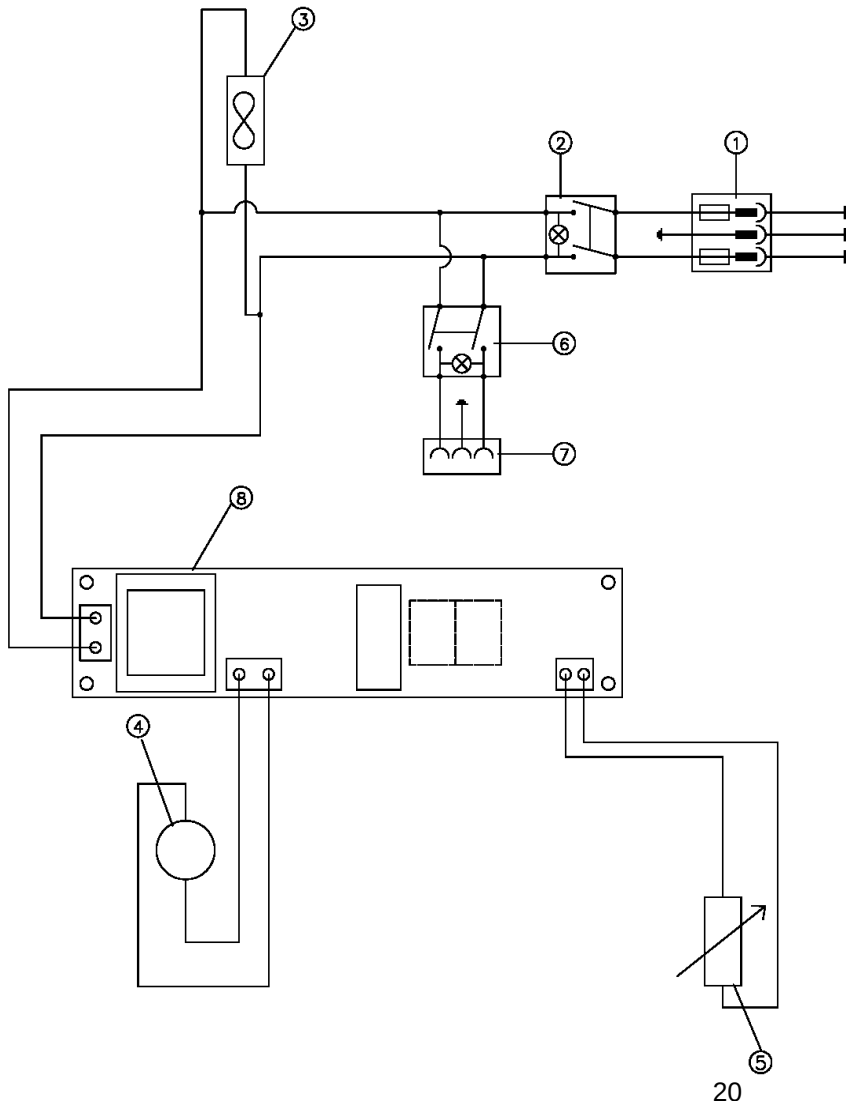
We, the manufacturer VELP Scientifica, under our responsibility declare that the product is manufactured in conformity with the following standards:

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EN 61010-1	Safety requirements for electrical equipment for measurement, control, and laboratory use - Part 1: General requirements
EN 61010-2-010	Particular requirements for laboratory equipment for the heating of materials
EN 61326-1	Electrical equipment for measurement, control and laboratory use. EMC requirements. General requirements

and satisfies the essential requirements of the following directives:

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9. Declaration of conformity ^{UK}CA

We, the manufacturer VELP Scientifica, under our responsibility declare that the product is manufactured in conformity with the following regulations:

S.I. 2016/1101 Electrical Equipment (Safety) Regulations 2016

S.I. 2016/1091 Electromagnetic Compatibility Regulations 2016

according to the relevant designated standards:

EN 61010-1 Safety requirements for electrical equipment for measurement, control, and laboratory use
Part 1: General requirements

EN 61010-2-010 Particular requirements for laboratory equipment for the heating of material

EN 61326-1 Electrical equipment for measurement, control and laboratory use - EMC requirements -
Part 1: General requirements

and satisfies the essential requirements of regulations:

S.I. 2008/1597 Supply of Machinery (Safety) Regulations 2008

S.I. 2012/3032 Restriction of the Use of Certain Hazardous Substances in Electrical and Electronic
Equipment Regulations 2012

S.I. 2013/3113 Waste Electrical and Electronic Equipment Regulations 2013

plus modifications.

Thank you for having chosen VELP!

Established in 1983, VELP is today one of the world's leading manufacturer of analytical instruments and laboratory equipment that has made an impact on the world-wide market with Italian products renowned for innovation, design and premium connectivity. VELP works according to **ISO 9001**, **ISO14001** and **ISO 45001** Quality System Certification.

Our instruments are manufactured in Italy according to the IEC 1010-1 and CE regulation.

Our product lines:

Analytical instruments

Elemental Analyzers
Digestion Units
Distillation Units
Solvent Extractors
Fiber Analyzers
Dietary Fiber Analyzers
Oxidation Stability Reactor
Consumables

Laboratory Equipment

Magnetic Stirrers
Heating Magnetic Stirrers
Heating Plates
Overhead stirrers
Vortex mixers
Dispersers
COD Thermoreactors
BOD and Respirometers
Cooled Incubators
Flocculators
Overhead Shakers
Turbidimeter
Open Circulating Baths
Pumps

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VELP Scientifica Srl