



**ermes**

CONNECT  
YOUR PRODUCT  
**NOW**

**OHS Advance**

[www.velp.com](http://www.velp.com)

# How to Connect to VELP Ermes Platform

## OHS Advance

In order to speed up the connection, be sure to get in contact with your IT specialist to get all your laboratory Wi-Fi connection details.

### Requirements:


1. **Device:** PC, Tablet, Smartphone with Wi-Fi connection
2. **Browser:** Chrome, Safari, Firefox, Edge, UC browser

## FIRST STEP | ACTIVATE VELP ERMES ACCOUNT

- Go to [www.velp.com](http://www.velp.com) through your internet browser.
- In the Smart Lab menu click on “Configure Your VELP Ermes Account”. Fill in the form and accept Terms and Conditions.
- You will receive an email containing your ID, Password and a link to proceed to VELP Ermes authentication at the email address that you entered during the registration.
- At the first login on VELP Ermes platform a new password will be requested: *remember to change it!*

## SECOND STEP | CONNECT TO YOUR WIFI NETWORK

- Turn on your OHS Advance
  - A. Press the menu key
  - B. Rotate the knob to select Set-up and press the knob
  - C. Rotate the knob to select Wi-Fi and press the knob
  - D. Rotate the knob to select AP and press the knob to confirm (see chapter 7.5)

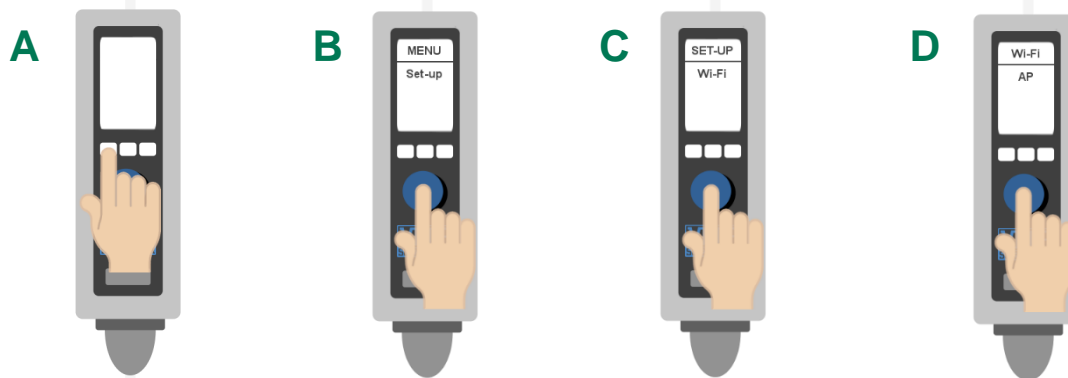

- From your device select the OHS\_(serial number) available on the Wi-Fi list, in order to connect the instrument.
- Open a browser from your device (PC, tablet, smartphone), enter 192.168.240.1 address. Use **admin** as username and password. A pop up will appear and proceed as following:
  1. Select Wireless and network configuration
  2. Press “Select an existing WiFi network name”
  3. Choose your WiFi network name\*
  4. Insert your passphrase (your password) \*
  5. Insert your passphrase into “confirm passphrase” to confirm and press next
  6. Press “Save and apply”

Wait until the “This site can’t be reached” appears on your screen.

\*If necessary, please contact your IT administrator.

### THIRD STEP | REGISTRATION ON VELP ERMES PLATFORM

**A**



Press the menu key and rotate the knob to select Service

**B**




From Service menu select ERMES.

**C**



On the display you will see OTC and "----".

**D**

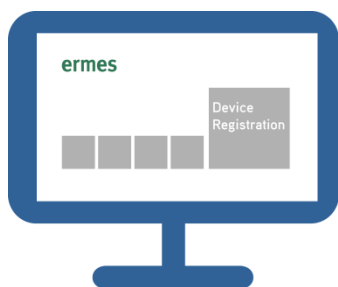

Reconnect your device to your WiFi network. Login on VELP Ermes with your ID and Password received by email, and select "Device Registration".

**E**



Enter the OHS Advance product code (you can find this information on the silver label on the back of it), serial number and purchase date. Press Next.

**F**




On the display of OHS Advance will appear a One Time Code (OTC) of four digits. Enter this number, Press Next and wait for connection.

**G**



Give a name to your OHS Advance and enter the address where your instrument is located. Press Next

**H**

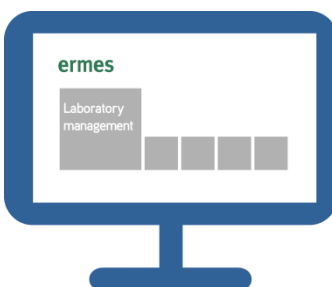

Assign your OHS Advance to a laboratory or create a new one selecting "Laboratory Management".

**I**



Now from "Instruments" you can monitor and control your connected OHS Advance.



## Always **Connected** and **Easy to Use**

Monitor and manage multiple instruments 24/7 from your device. Real time visualization of your analysis and of the instrument working conditions. Immediate notifications and alerts.



## **Safely** Manage **Your Data**

VELP Ermes collects and stores your data with the maximum level of encryption following the highest cyber-security standards. You are now able to have access to your data anytime, anywhere. VELP Ermes will allow you to easily generate and share reports of your analyses.



## Your Instrument **Always Updated**

Get software updates (OTA) immediately and benefit of all new features coming from VELP continuous product innovation.



## Enhanced **Service Support**

Faster and smarter Service Support thanks to the direct contact with VELP specialists. With VELP Ermes we can guide you to get the maximum from your instrument.