***Instruction Manual***  
***Manuale di istruzioni***  
***Manuel d'instructions***  
***Manual de instrucciones***  
***Bedienungsanleitung***  
***指导手册***

## **MSL 8 Magnetic Stirrer**

F203A0513

### **General Information / Informazioni Generali / Informations Générales / Información General / Allgemeine Hinweise / 一般信息**

Before using the unit, please read the following instruction manual carefully.

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Do not dispose of this equipment as urban waste, in accordance with EEC directive 2002/96/CE.

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**This unit must be used for laboratory applications indoor only.** The manufacturer declines all responsibility for any use of the unit that does not comply with these instructions. If the product is used in a not specified way by the manufacturer or with non specified accessories, product's safety may be compromised.

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**This unit has been designed and manufactured in compliance with the following standards:**

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Safety requirements for electrical equipment for measurement, control and for laboratory use

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**IEC/EN 61010-1  
IEC/EN 61010-2-051**

Electrical equipment for laboratory use

**UL 61010-1**

General requirement - Canadian electrical code

**CAN/CSA-C22.2 No.61010-1**

VELP reserves the right to modify the characteristics of its products with the aim to constantly improving their quality.

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The plug disconnects the instrument. Therefore, place the instrument where it can be quickly disconnected.

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The stirred solution may release toxic, dangerous or poisonous gases. Adequate safety measures must be taken, in accordance with the safety regulations in force, including the presence of hood and personal protective equipment (masks, gloves, goggles, etc.).

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Do not use with explosive and dangerous materials for which the equipment is not designed. The stirrer must not be used in explosive atmospheres, in bain-marie and to stir combustible liquids that have a low combustion temperature. The minimum fire point of flammable solution is 110 °C. Only small amounts (< 50 ml) of flammable liquid can be used with the device.

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It is responsibility of the user appropriately decontaminate the instrument in case of dangerous substances fall on or in it. It is also responsibility of the user to use safety substances for cleaning or decontaminating, which do not react with internal parts of the instrument or with the material contained in it. In case of doubts on the compatibility of a cleaning solution, contact the manufacturer or local distributor.

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In case of black-out, once the power is back the device will restart automatically.

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MSL 8 is able to stir up to 8L. Due to the new brushless motor, the stirrer stays cold even after several days of continuous operation. This feature makes it particularly appreciated in microbiology and biochemistry. Epoxy painted metal structure, with stainless steel cover, is studied and tested in order to give to the instrument an excellent resistance to the attack of chemical agents and to corrosion in general. The stirring speed can be set within a range of 40 – 1500 rpm using the knob on the front panel.

**NOTE:** it's important to choose the most suitable magnetic stirring bar related to the quantity and the liquid to be stirred as well as to the type of the beaker.

## 2. Assembly and installation

- Unpacking
  - Check the integrity of the unit after unpacking.
- The box includes
  - MSL 8 magnetic stirrer
  - Instruction manual
  - Plug for power supply
  - Switching VELP 100-240V/12V without plug
- First installation
  - Place the unit on non-flammable surface at least 30 cm from the wall.
  - Make sure that the rating value of the instrument corresponds to the one of the power supply
  - Ensure that the socket provided and the relative cut-off device are compliant with the current safety norms and easy to reach. Use only the power supply provided with the instrument.
  - Place the flasks containing the sample and a suitable magnetic stirring bar on the stirring plate.

## 3. Working

<b>Commissioning</b>	<ul style="list-style-type: none"> <li>➤ Rotate the speed control knob to zero position. Insert the transformer into the socket.</li> <li>➤ Place the flasks containing the sample and a suitable magnetic stirring bar on the stirring plate.</li> </ul>
<b>Stirring</b>	<ul style="list-style-type: none"> <li>➤ Adjust speed set point by turning the speed control knob.</li> <li>➤ Speed increases until set point is achieved.</li> <li>➤ A microprocessor ensures constant speed (counter-reaction).</li> <li>➤ Switch off the stirring by turning the knob to zero position.</li> </ul>

## 4. Maintenance

<b>Maintenance</b>	<ul style="list-style-type: none"> <li>➤ No routine or extraordinary maintenance is necessary;</li> <li>➤ Repairs must be carried out by authorized Velp personnel only;</li> <li>➤ Instrument must be transported in its original packaging any indications present on the original packaging must be followed (e.g. palletized);</li> <li>➤ It is the responsibility of the user, to properly decontaminate the unit in case of hazardous substances remaining on the surface or interior of the device. If in doubt about the compatibility of a cleaning or decontamination product, contact the manufacturer or distributor.</li> </ul>
<b>Cleaning</b>	<ul style="list-style-type: none"> <li>➤ Disconnect the unit from the power supply and use a cloth dampened with a non-flammable non-aggressive detergent.</li> </ul>

## 5. Technical data

<b>General features</b>	Model	F203A0513
	Power supply	0.5 A / 100-240 V / 50-60 Hz (+/-10%)
	Dimensions (WxHxD)	168x60x183 mm
	Weight	0,9 kg
	Power input	12 Vdc 4,5 W
	Construction material (structure)	Epoxy painted metal structure and stainless steel
	Working in continuous	Admitted
	Noisiness	<< 60 dBa
	Environmental temperature admitted	+5...+40 °C
	Storage temperature admitted	-10...+60 °C
	Max humidity	80%
	Overvoltage category	II
	Pollution degree CEI EN61010-1	2
	Max altitude	Up to 2000 m
<b>Stir</b>	Stirring capacity	8 L H <sub>2</sub> O
	Speed range	40 - 1500rpm
	Motor type	BLDC
	Speed selection	Analogic

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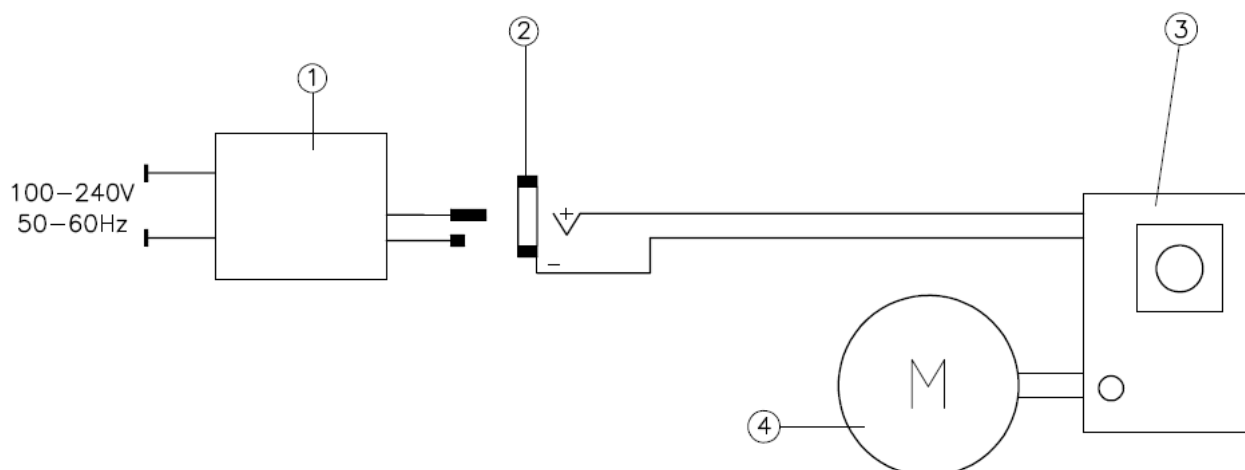
We, the manufacturer VELP Scientifica, under our responsibility declare that the product is manufactured in conformity with the following standards:

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2015/863/EU (RoHS III)

2012/19/EU (WEEE)

and satisfies the essential requirements of the following directives:

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## 9. Declaration of conformity <sup>UK</sup> <sub>CA</sub>

We, the manufacturer VELP Scientifica, under our responsibility declare that the product is manufactured in conformity with the following regulations:

S.I. 2016/1101 Electrical Equipment (Safety) Regulations 2016

S.I. 2016/1091 Electromagnetic Compatibility Regulations 2016

according to the relevant designated standards:

EN 61010-1 Safety requirements for electrical equipment for measurement, control, and laboratory use  
Part 1: General requirements

EN 61010-2-051 Particular requirements for laboratory equipment for mixing and stirring

EN 61326-1 Electrical equipment for measurement, control and laboratory use - EMC requirements -  
Part 1: General requirements

and satisfies the essential requirements of regulations:

S.I. 2008/1597 Supply of Machinery (Safety) Regulations 2008

S.I. 2012/3032 Restriction of the Use of Certain Hazardous Substances in Electrical and Electronic  
Equipment Regulations 2012

S.I. 2013/3113 Waste Electrical and Electronic Equipment Regulations 2013

plus modifications.

## Thank you for having chosen VELP!

Established in 1983, VELP is today one of the world's leading manufacturer of analytical instruments and laboratory equipment that has made an impact on the world-wide market with Italian products renowned for innovation, design and premium connectivity. VELP works according to **ISO 9001**, **ISO14001** and **ISO 45001** Quality System Certification.

Our instruments are manufactured in Italy according to the IEC 1010-1 and CE regulation.

Our product lines:

### Analytical instruments

Elemental Analyzers  
Digestion Units  
Distillation Units  
Solvent Extractors  
Fiber Analyzers  
Dietary Fiber Analyzers  
Oxidation Stability Reactor  
Consumables

### Laboratory Equipment

Magnetic Stirrers  
Heating Magnetic Stirrers  
Heating Plates  
Overhead stirrers  
Vortex mixers  
Dispersers  
COD Thermoreactors  
BOD and Respirometers  
Cooled Incubators  
Flocculators  
Overhead Shakers  
Turbidimeter  
Open Circulating Baths  
Pumps

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VELP Scientifica Srl