Instruction Manual
Manuale di istruzioni
Manuel d'instructions
Manual de instrucciones
Bedienungsanleitung
指导手册

MULTISTIRRER 6 & 15 DIGITAL Magnetic Stirrers

F203A0179, F203A0180

**General Information / Informazioni Generali / Informations Générales / Información General /
Allgemeine Hinweise / 一般信息**

Before using the unit, please read the following instruction manual carefully.

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Do not dispose of this equipment as urban waste, in accordance with EEC directive 2002/96/CE.

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This unit must be used for laboratory applications indoor only. The manufacturer declines all responsibility for any use of the unit that does not comply with these instructions. If the product is used in a not specified way by the manufacturer or with non specified accessories, product's safety may be compromised.

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This unit has been designed and manufactured in compliance with the following standards:

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Safety requirements for electrical equipment for measurement, control and for laboratory use

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IEC/EN 61010-1

IEC/EN 61010-2-051

Electrical equipment for laboratory use

UL 61010-1

General requirement - Canadian electrical code

CAN/CSA-C22.2 No.61010-1

VELP reserves the right to modify the characteristics of its products with the aim to constantly improving their quality.

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The plug disconnects the instrument. Therefore, place the instrument where it can be quickly disconnected.

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The stirred solution may release toxic, dangerous or poisonous gases. Adequate safety measures must be taken, in accordance with the safety regulations in force, including the presence of hood and personal protective equipment (masks, gloves, goggles, etc.).

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Do not use with explosive and dangerous materials for which the equipment is not designed. The stirrer must not be used in explosive atmospheres.

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It is responsibility of the user appropriately decontaminate the instrument in case of dangerous substances fall on or in it.

It is also responsibility of the user to use safety substances for cleaning or decontaminating, which do not react with internal parts of the instrument or with the material contained in it. In case of doubts on the compatibility of a cleaning solution, contact the manufacturer or local distributor.

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In case of black-out, once the power is back the device will restart automatically.

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Do not use with explosive and dangerous materials for which the equipment is not designed. The stirrer must not be used in explosive atmospheres, in bain-marie and to stir combustible liquids that have a low combustion temperature. The minimum fire point of flammable solution is 200 °C. Only small amounts (< 50 ml) of flammable liquid can be used with the device.

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MULTISTIRRER Digital is able to stir up to 6 beakers having a maximum diameter of 85 mm (MULTISTIRRER Digital 6) or up to 15 beakers having a maximum diameter of 64 mm (MULTISTIRRER Digital 15) at the same time.

Thanks to the brushless motors, the stirrer stays cold even after several days of continuous operation

Epoxy painted metal structure, with stainless steel cover, are studied and tested in order to give to the instrument an excellent resistance to the attack of chemical agents and to corrosion in general.

The stirring speed can be set using the frontal knob within a range of 80 - 1500 rpm.

The stirrer display shows the speed set point.

The instrument is equipped with advanced functions as auto-reverse of the stirring direction, intermittent mode programs and integrated timer to meet users' workflow and safety lock feature to avoid possible alteration of the working conditions.

NOTE: *it's important to choose the most suitable magnetic stirring bar related to the quantity and the liquid to be stirred as well as to the type of the beaker.*

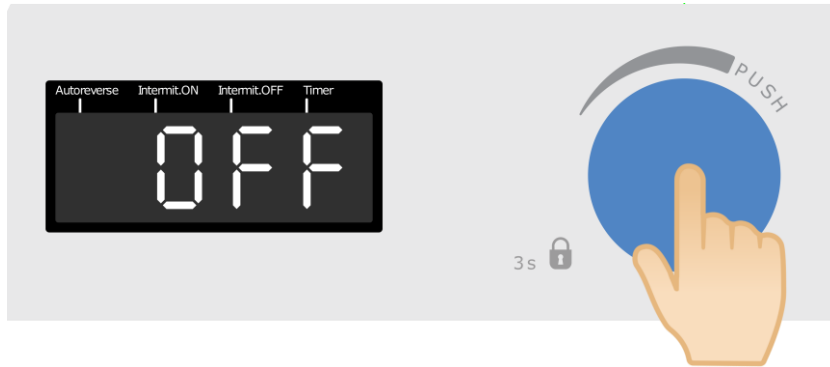
For MULTISTIRRER Digital 6 is suggested the stirring bar A00001056 (6x35 mm); for MULTISTIRRER Digital 15 the A00001057 (6x20 mm).

2. Installation


- Unpacking
 - Check the integrity of the unit after unpacking
- The box includes
 - MULTISTIRRER Digital magnetic stirrer
 - Instruction manual
 - Plug for power supply
 - Switching power supply 100-240 V / 12 V
- Installation
 - Place the unit on a non-flammable surface
 - Make sure that the technical specification of the unit and the power supply line are the same
 - Place the flasks containing the sample and a suitable magnetic stirring bar on the stirring plate
 - Connect the instrument to an easily accessible socket (compliant with the current safety norms), using only the provided power supply and dedicated plug

3. Operation

- Push the knob to start stirring.

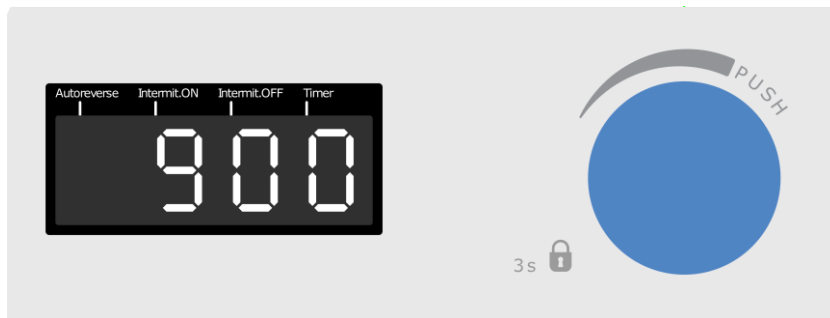
- Rotate the knob to set the stirring speed.


Note: push the knob to stop stirring.

Autoreverse, intermittent mode and timer functions

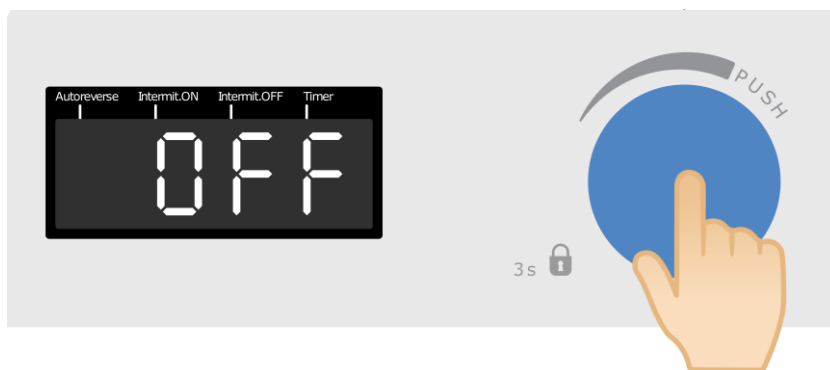- While the display shows OFF, push the knob for 3 seconds.

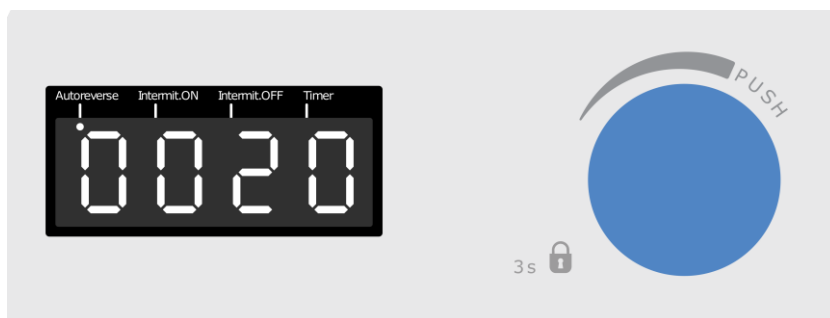
- Rotate the knob to set the autoreverse time (minutes/seconds).

The rotation direction will change after each set interval time.


Note: select NNSS to disable the function.

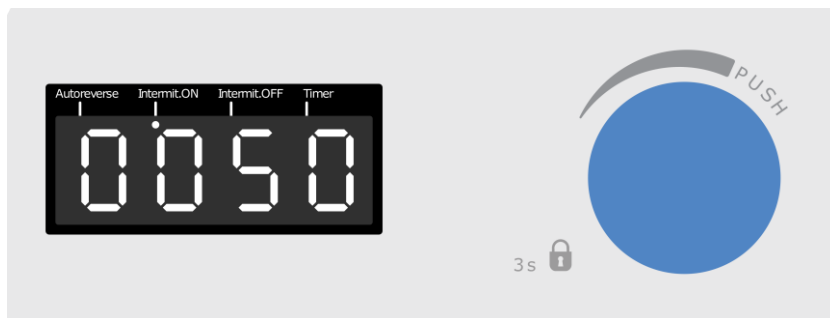- Push the knob. Then rotate it to set the run time of the intermittent mode (minutes/seconds).

Note: select *NNSS* to disable the function.

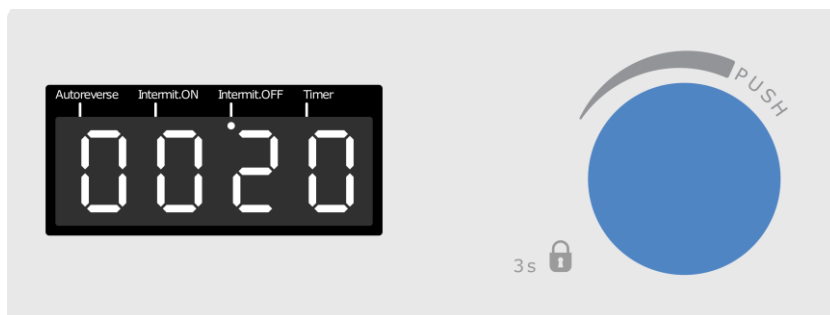
- Push the knob. Then rotate it to set the pause time of the intermittent mode (minutes/seconds).

The stirring function will be interrupted cyclically. The run time and the pause time should both be set.

Note: select *NNSS* to disable the function.

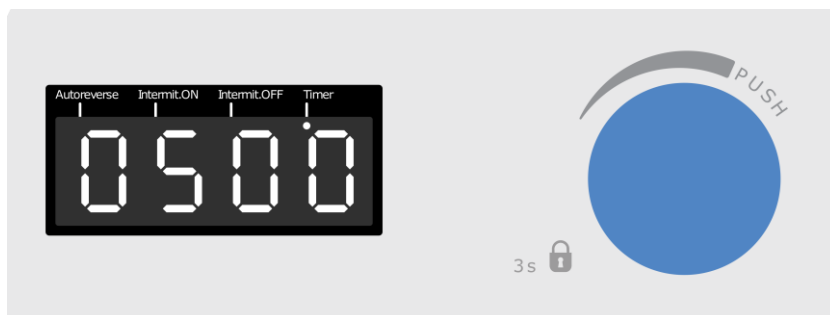- Push the knob. Then rotate it to set the timer (hours/minutes).

After the timer has elapsed, the stirring will be stopped and the display will show *End*.

Note: select *HHnn* to disable the function.

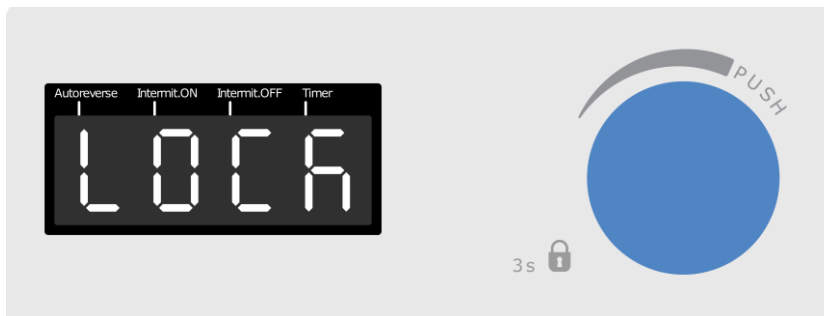
- Push the knob or wait 10 seconds to confirm.

Dedicated blinking leds show the active functions.

Note: each function can be set individually or in combination with the others.

Lock function

- While the unit is working, push the knob for 3 seconds to enable the lock function.

If the lock function is enabled, it is not possible to change the work conditions. The display shows *Lock* if the knob is pressed or rotated.

Note: push the knob for 3 seconds to disable the function.

4. Maintenance

No routine or extraordinary maintenance is necessary except periodically cleaning.

Cleaning

Disconnect the unit from the power supply and use a cloth dampened with a non-flammable non-aggressive detergent.

Repair

Repairs must be carried out by authorized Velp personnel only.

The instrument must be moved using the original shockproof packaging when freight forwarders, couriers or other are taking care of the unit transport. Follow the indications on the original packaging (e.g. place the instrument on a pallet). It is the responsibility of the user, to properly decontaminate the unit in case of hazardous substances remaining on the surface or interior of the device. If in doubt about the compatibility of a cleaning or decontamination product, contact the manufacturer or distributor.

5. Technical data

		MULTISTIRRER Digital 6	MULTISTIRRER Digital 15
General features	Models	F203A0179	F203A0180
	Voltage	100 – 240 V / 50 – 60 Hz	
	Power input	12 VDC 4.1 W	12 VDC 9 W
	Current consumption	0.35 A	0.75 A
	Dimensions (WxHxD)	230x51.5x370 mm	
	Weight	1.75 kg	2.1 kg
	Construction material	Epoxy painted metal structure and stainless steel	
	Timer	1 min ÷ 99 h 59 min	
	Autoreverse / Intermittent mode	5 s ÷ 99 min 55s	
	Working in continuous	Admitted	
	Noisiness	< 35 dBA	
	Environmental temperature admitted	+5...+40 °C	
	Storage temperature admitted	-10...+60 °C	
	Max humidity	80%	
	Overvoltage category	II	
	Pollution degree CEI EN61010-1	2	
	Max altitude	2000 m	
Stir	Stirring capacity	6x400 ml H ₂ O	15x250 ml H ₂ O
	Speed range	0 ÷ 1500 rpm	
	Motors type	BLDC	
	Speed selection	10 rpm	

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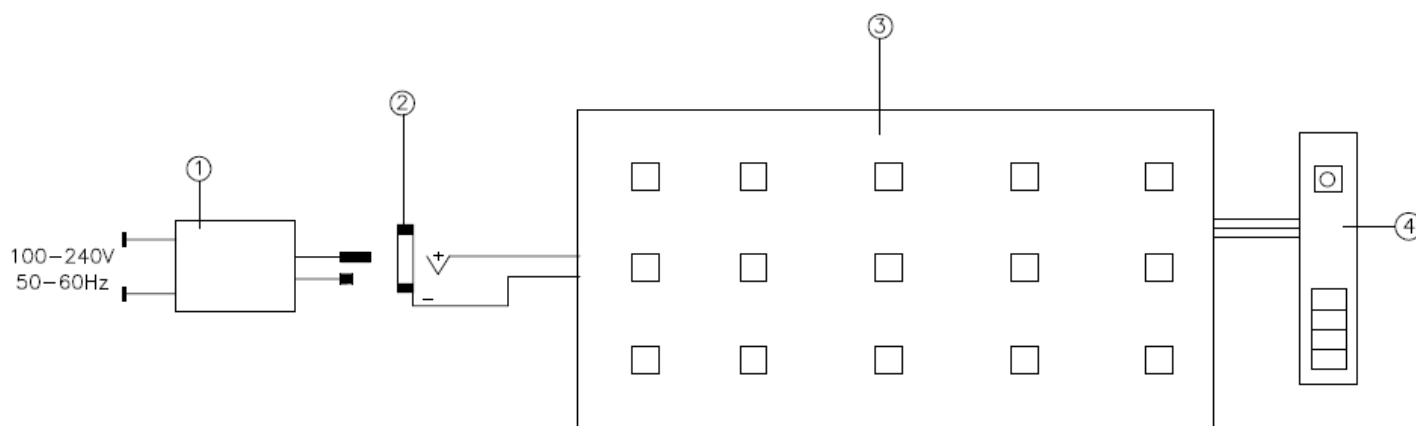
We, the manufacturer VELP Scientifica, under our responsibility declare that the product is manufactured in conformity with the following standards:

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2015/863/EU (RoHS III)

2012/19/EU (WEEE)

and satisfies the essential requirements of the following directives:

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9. Declaration of conformity

We, the manufacturer VELP Scientifica, under our responsibility declare that the product is manufactured in conformity with the following regulations:

S.I. 2016/1101 Electrical Equipment (Safety) Regulations 2016

S.I. 2016/1091 Electromagnetic Compatibility Regulations 2016

according to the relevant designated standards:

EN 61010-1 Safety requirements for electrical equipment for measurement, control, and laboratory use
Part 1: General requirements

EN 61010-2-051 Particular requirements for laboratory equipment for mixing and stirring

EN 61326-1 Electrical equipment for measurement, control and laboratory use - EMC requirements -
Part 1: General requirements

and satisfies the essential requirements of regulations:

S.I. 2008/1597 Supply of Machinery (Safety) Regulations 2008

S.I. 2012/3032 Restriction of the Use of Certain Hazardous Substances in Electrical and Electronic
Equipment Regulations 2012

S.I. 2013/3113 Waste Electrical and Electronic Equipment Regulations 2013

plus modifications.

Thank you for having chosen VELP!

Established in 1983, VELP is today one of the world's leading manufacturer of analytical instruments and laboratory equipment that has made an impact on the world-wide market with Italian products renowned for innovation, design and premium connectivity. VELP works according to **ISO 9001**, **ISO14001** and **ISO 45001** Quality System Certification.

Our instruments are manufactured in Italy according to the IEC 1010-1 and CE regulation.

Our product lines:

Analytical instruments

Elemental Analyzers
Digestion Units
Distillation Units
Solvent Extractors
Fiber Analyzers
Dietary Fiber Analyzers
Oxidation Stability Reactor
Consumables

Laboratory Equipment

Magnetic Stirrers
Heating Magnetic Stirrers
Heating Plates
Overhead stirrers
Vortex mixers
Dispersers
COD Thermoreactors
BOD and Respirometers
Cooled Incubators
Flocculators
Overhead Shakers
Turbidimeter
Open Circulating Baths
Pumps

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VELP Scientifica Srl