Instruction Manual
Manuale di istruzioni
Manuel d'instructions
Manual de instrucciones
Bedienungsanleitung
指导手册

MULTISTIRRER 6 & 15 Magnetic Stirrers

F203A0177, F203A0178

**General Information / Informazioni Generali / Informations Générales / Información General /
Allgemeine Hinweise / 一般信息**

Before using the unit, please read the following instruction manual carefully.

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Do not dispose of this equipment as urban waste, in accordance with EEC directive 2002/96/CE.

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This unit must be used for laboratory applications indoor only. The manufacturer declines all responsibility for any use of the unit that does not comply with these instructions. If the product is used in a not specified way by the manufacturer or with non specified accessories, product's safety may be compromised.

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This unit has been designed and manufactured in compliance with the following standards:

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Safety requirements for electrical equipment for measurement, control and for laboratory use

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IEC/EN 61010-1
IEC/EN 61010-2-051

Electrical equipment for laboratory use

UL 61010-1

General requirement - Canadian electrical code

CAN/CSA-C22.2 No.61010-1

VELP reserves the right to modify the characteristics of its products with the aim to constantly improving their quality.

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The plug disconnects the instrument. Therefore, place the instrument where it can be quickly disconnected.

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The stirred solution may release toxic, dangerous or poisonous gases. Adequate safety measures must be taken, in accordance with the safety regulations in force, including the presence of hood and personal protective equipment (masks, gloves, goggles, etc.).

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Do not use with explosive and dangerous materials for which the equipment is not designed. The stirrer must not be used in explosive atmospheres.

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It is responsibility of the user appropriately decontaminate the instrument in case of dangerous substances fall on or in it. It is also responsibility of the user to use safety substances for cleaning or decontaminating, which do not react with internal parts of the instrument or with the material contained in it. In case of doubts on the compatibility of a cleaning solution, contact the manufacturer or local distributor.

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Do not use with explosive and dangerous materials for which the equipment is not designed. The stirrer must not be used in explosive atmospheres, in bain-marie and to stir combustible liquids that have a low combustion temperature. The minimum fire point of flammable solution is 200 °C. Only small amounts (< 50 ml) of flammable liquid can be used with the device.

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In case of black-out, once the power is back the device will restart automatically.

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MULTISTIRRER is able to stir up to 6 beakers having a maximum diameter of 85 mm (MULTISTIRRER 6) or up to 15 beakers having a maximum diameter of 64 mm (MULTISTIRRER 15) at the same time. The stirrer stays cold even after several days of continuous operation. This feature makes it particularly appreciated in microbiology and biochemistry. Epoxy painted metal structure, with stainless steel cover, is studied and tested in order to give to the instrument an excellent resistance to the attack of chemical agents and to corrosion in general. The stirring speed can be set within a range of 80 - 1500 rpm using the knob on the front panel. The Stirrer led shows when the stirring is running.

NOTE: it's important to choose the most suitable magnetic stirring bar related to the quantity and the liquid to be stirred as well as to the type of the beaker. For MULTISTIRRER 6 is suggested the stirring bar A00001056 (6x35 mm); for MULTISTIRRER 15 the A00001057 (6x20 mm).

2. Assembly and installation

Upon receipt and after having removed the packaging, please check the integrity of the instrument. The box includes:

- Magnetic stirrer MULTISTIRRER
- Plug for power supply
- Switching VELP 100-240V/12V without plug
- Instruction manual

2.1 Electrical connection

After having unpacked the instrument, place the unit on the laboratory bench.

Before connecting the instrument to the power supply, make sure that the values on the rating plate correspond to those of the power supply. Connect the unit to the power supply using the transformer supplied.

Ensure that the socket and the relative cut-off device conform to current safety norms and easy to reach.

2.2 Start-up

Rotate the speed knob completely to the left. Place the flasks containing the sample and a suitable magnetic stirring bar on the stirring plate. Then, set the speed by turning the speed knob.

3. Operating controls

Turn the stirring speed knob on the front panel in order to start the stirring function.

It is possible to select speeds between 80 and 1500 rpm. When the led is lighted it means that the stirring is on.

4. Maintenance

No routine or extraordinary maintenance is necessary apart from periodically cleaning the unit as described in this manual. In compliance with the product guarantee law, repairs to our units must be carried out in our factory, unless previously agreed otherwise with local distributors. The instrument must be transported in its original packaging and any indications present on the original packaging must be followed (e.g. palletized).

It is the responsibility of the user, to properly decontaminate the unit in case of hazardous substances remaining on the surface or interior of the device. If in doubt about the compatibility of a cleaning or decontamination product, contact the manufacturer or distributor.

4.1 Cleaning

Disconnect the unit from the power supply and use a cloth dampened with an non-inflammable non-aggressive detergent.

5. Technical data

Power	MULTISTIRRER 6: 12 Vdc 3,6 W	MULTISTIRRER 15: 12 Vdc 9 W
Admitted power supply	0.5 A / 100-240 V / 50-60 Hz (+/-10%)	
Dimensions (WxHxD)	230x51.5x370 mm	
Weight (with plug)	MULTISTIRRER 6: 1,75 kg	MULTISTIRRER 15: 2,1 kg
Speed range	80 ÷ 1500 rpm	
Stirring capacity	MULTISTIRRER 6: 6x400 ml	MULTISTIRRER 15: 15x250 ml
Operating mode	Continuous	
Construction material	Epoxy painted metal structure and stainless steel	
Environmental ambient temperature	5 - 40 °C (41 - 104°F)	
Storage temperature range	-10... + 60 °C (14 - 140°F)	
Pollution degree	2	
Overvoltage category	II	
Sound level	35 dBa	
Max humidity	80%	
Altitude	Up to 2000 m.	

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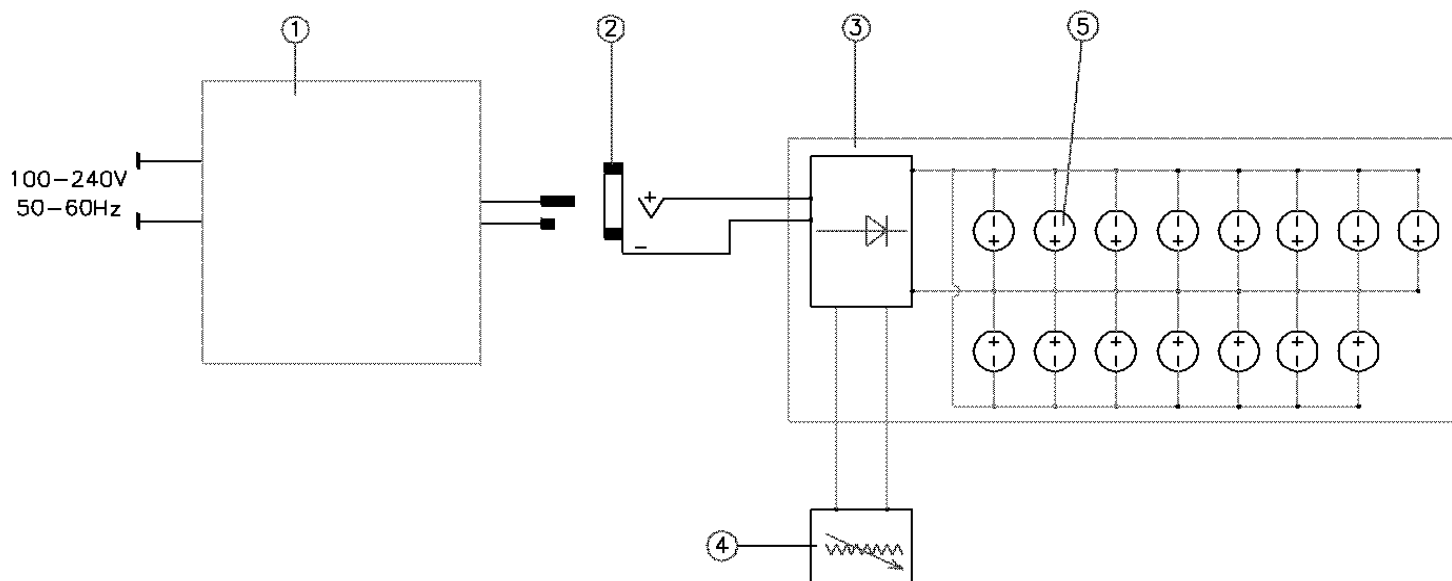
We, the manufacturer VELP Scientifica, under our responsibility declare that the product is manufactured in conformity with the following standards:

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9. Declaration of conformity ^{UK}CA

We, the manufacturer VELP Scientifica, under our responsibility declare that the product is manufactured in conformity with the following regulations:

S.I. 2016/1101 Electrical Equipment (Safety) Regulations 2016

S.I. 2016/1091 Electromagnetic Compatibility Regulations 2016

according to the relevant designated standards:

EN 61010-1 Safety requirements for electrical equipment for measurement, control, and laboratory use
Part 1: General requirements

EN 61010-2-051 Particular requirements for laboratory equipment for mixing and stirring

EN 61326-1 Electrical equipment for measurement, control and laboratory use - EMC requirements -
Part 1: General requirements

and satisfies the essential requirements of regulations:

S.I. 2008/1597 Supply of Machinery (Safety) Regulations 2008

S.I. 2012/3032 Restriction of the Use of Certain Hazardous Substances in Electrical and Electronic
Equipment Regulations 2012

S.I. 2013/3113 Waste Electrical and Electronic Equipment Regulations 2013

plus modifications.

Thank you for having chosen VELP!

Established in 1983, VELP is today one of the world's leading manufacturer of analytical instruments and laboratory equipment that has made an impact on the world-wide market with Italian products renowned for innovation, design and premium connectivity. VELP works according to **ISO 9001**, **ISO14001** and **ISO 45001** Quality System Certification.

Our instruments are manufactured in Italy according to the IEC 1010-1 and CE regulation.

Our product lines:

Analytical instruments

Elemental Analyzers
Digestion Units
Distillation Units
Solvent Extractors
Fiber Analyzers
Dietary Fiber Analyzers
Oxidation Stability Reactor
Consumables

Laboratory Equipment

Magnetic Stirrers
Heating Magnetic Stirrers
Heating Plates
Overhead stirrers
Vortex mixers
Dispersers
COD Thermoreactors
BOD and Respirometers
Cooled Incubators
Flocculators
Overhead Shakers
Turbidimeter
Open Circulating Baths
Pumps

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VELP Scientifica Srl