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This unit must be used for laboratory applications only and for internal use only.

The manufacturer declines all responsibility for any use of the unit that does not comply with these instructions.

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VELP reserves the right to modify the characteristics of its products to constantly improve their quality.

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The VELP OHS/ControllerSoft software is designed to control working parameters such as temperature, speed (rpm) and torque of VELP OHS Series overhead stirrers and VELP Controller Advance.

2. Assembling and installing

Upon receipt and after having removed the packaging, please check the integrity of the instrument. The box includes:

- USB cable
- OHS/ControllerSoft™ Software

The following VELP instruments are enabled to work with the software by connecting to the PC via USB:

OHS 100 Advance 230V/50-60Hz	F20100481
OHS 100 Advance 115V/60Hz	F20110481
OHS 100 Advance 115V/60Hz CG-2054-100	F20110481CG
OHS 200 Advance 230V/50-60Hz	F20100480
OHS 200 Advance 115V/60Hz	F20110480
OHS 60 Advance 230V/50-60Hz	F20100482
OHS 60 Advance 115V/60Hz	F20110482
Controller Advance 115-230V/50-60Hz	F201A0500
Controller Advance CG-2033-V	F201A0500CG

The OHS/ControllerSoft™ Software is able to manage up to 4 units independently.

The OHS/ControllerSoft™ Software requires an activation code in order to establish the connection with the instrument. To do so is mandatory to have the serial number of the VELP instrument and the serial number of the PC where the software is installed.

The USB stick contains the drivers that have to be installed to recognize the instrument at the first connection to the PC.

3. OHS/ControllerSoft™

3.1 Installation

General rules for installing the OHS/ControllerSoft™ Software:



In order to install the OHS/ControllerSoft™_Software correctly, administrator rights for the PC must be available. The PC on which the OHS/ControllerSoft™ Software will be installed must not be overloaded with other software. This is to prevent unstable operation of the OHS/ControllerSoft™ Software.

To install the OHS/ControllerSoft™ Software:

1. Place the USB stick in the PC. A folder should open automatically.
2. If no folder automatically open, access to the USB and run **setup.exe** on the USB stick folder.
3. The installation guide is then displayed. Click **Next** and follow the instructions. The OHS/Controllersoft™ icon will be created on the screen automatically.
4. Connect the VELP instrument (OHS advance models and/or Controller advance) to the PC, using the USB cable.
5. Install the drivers, opening the Driver folder on the USB and following the instructions that appear on the screen.
6. Re-boot the computer to complete the installation.


3.2 Activation

1. Open the software and go in Help/Software activation/Activation form and fill with the data request. .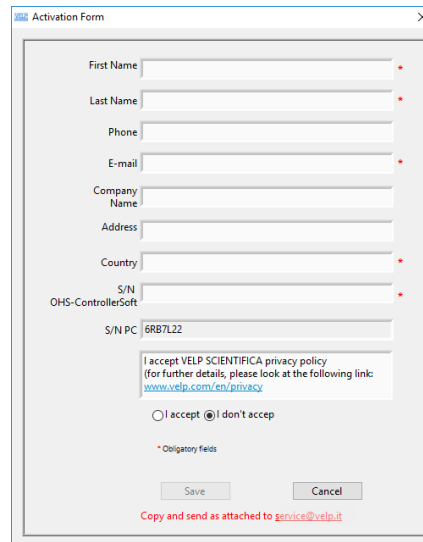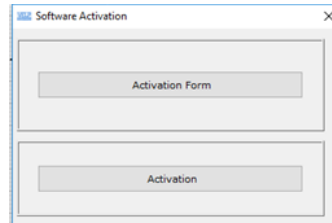


The S/N PC is filled automatically by the software

2. Save the data. The software created a request file in C:\OHS-ControllerSoft\Config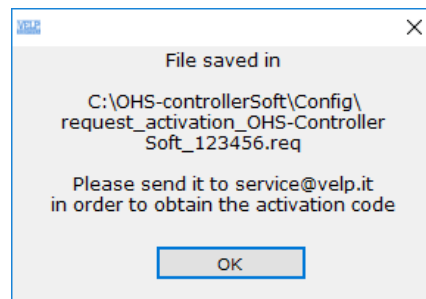

3. Send this file to service@velp.it to obtain the activation file.
4. When received it, go in the Help/Software activation/Activation and load the file received. Software will show if the activation run well: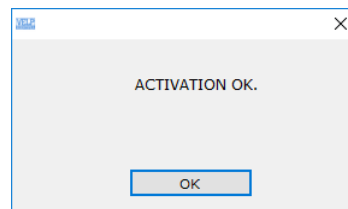

Software can be installed in more than one PC but it can be used only to review the files previously created.

4. OHS/Controllersoft™ Quick Start Guide

This page shows an abstract of the OHS/ControllerSoft™


The purpose is to help the user identify the various possibilities offered by the instrument and how to reach them easily.

Menu	Submenu	Description
File	Save data	Data of temperature, rpm and torque of all the instruments connected are saved in a dedicated file (.csv) These data saved are the ones visualized until the saving moment. (a file for each instrument connected)
	Reset data	Cleaning of the graph by the data visualized until that moment.
	Exit	Exit from the program.
Record	Record	Start to record the data of the temperature, rpm and torque in a dedicated file (.csv)
	Review	Load the file previously recorded
Service	Find device	Connect the instruments to the OHS/ControllerSoft™
	Name table	Give the name to the instruments connected
Help	Ermes	Information about ermes solution
	Software activation	Activate the software
	About	How to contact VELP SCIENTIFICA

5. Instruments and OHS/Controllersoft™ Connection

In order to connect the instrument to the PC, is necessary to use the USB cable, supplied with the software.


1. After connecting it to the PC and activating the software, select **Find device** from the **Service** menu. The connection is automatically found.
2. When the unit is connected to the OHS/ControllerSoft™ the software shows the S/N in the right box and the graph starts to visualize the temperature, speed (rpm) and torque.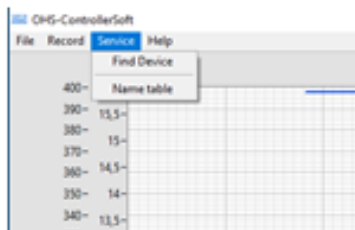

6. OHS/Controllersoft™ Workflow

6.1 Name the instrument

It is possible to give the name to the instrument connected, filling the table in **Service/Name table**: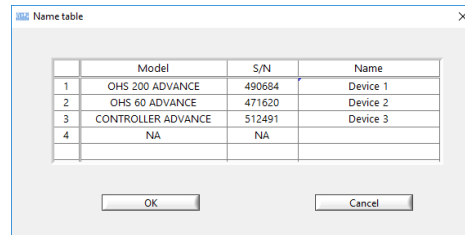

	Model	S/N	Name
1	OHS 200 ADVANCE	490684	Device 1
2	OHS 60 ADVANCE	471620	Device 2
3	CONTROLLER ADVANCE	512491	Device 3
4	NA	NA	


6.2 Visualize the data

The graph of the main screen shows as default the data related to the temperature, rpm and torque. From the right box it is possible to select or unselect the data to be visualized. The graph visualizes the data of the last 10 working hours, without saving.

In **File/Save data** it is possible to save in a dedicated file (.csv) the data visualized until the saving moment.

In **File/Reset data** it is possible to clean the graph.


In **Record/Review** it is possible to load the file previously recorded, pressing the button LOAD.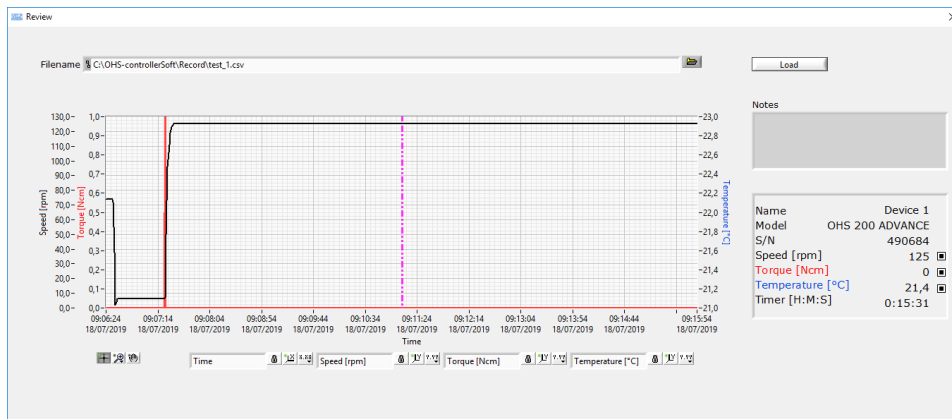

On the main screen is visualized also the working time or the countdown if it is set on the instrument.


6.3 Record the data

In order to start to record the data, select **Record/Record** menu. The following panel will be opened:

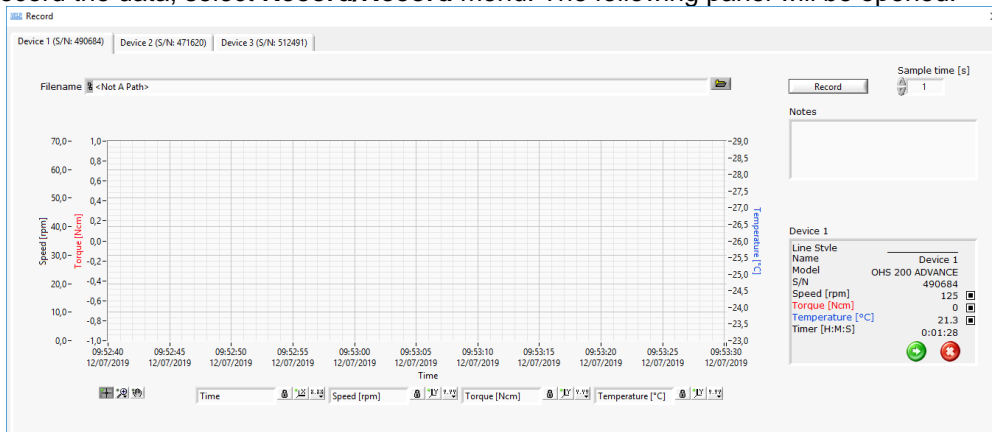Press the RECORD button to start the recording. The green indication "Recording" in the Recording window and in the main window shows that the recording step is active.

Press STOP button to stop the recording.

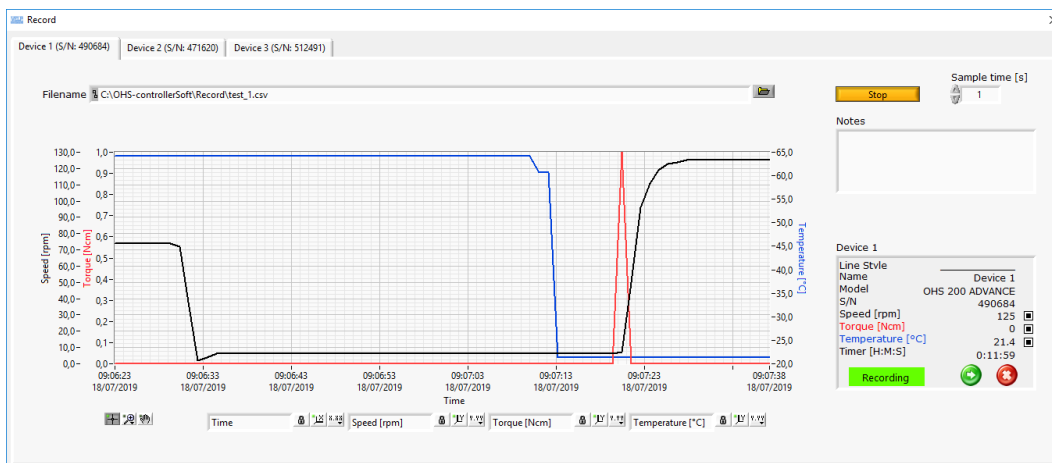In this panel is opened one “recording” window for each instrument in order to manage the instruments separately. It is possible to choose the sampling time, writing a whole number from 1 second (set as default)

6.4 Start the instrument

It is possible to start the instrument by the OHS/Controllersoft™ pressing the icon  in the main screen or on the recording window. The instrument will start to work as per the parameters set on the instrument.

6.5 Stop the instrument

It is possible to stop the instrument by the OHS/Controllersoft™ pressing the icon  in the main screen or on the recording window.

7. PC Technical requirements

Microsoft Window 7 (SP1), 8 or 10.
Pentium IV processor 2 GHz or more
RAM 4 GB
50 MB of hard disk memory free, dedicated
Minimum n° 1 USB-port (for communication between the instrument and PC)
Minimum recommended graphic resolution: 1366x768

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The image shows two screenshots of a software activation process. The top screenshot is a window titled "Software Activation" with two buttons: "Activation Form" and "Activation". The bottom screenshot is the "Activation Form" window, which contains the following fields: First Name, Last Name, Phone, E-mail, Company Name, Address, Country, S/N OHS-ControllerSoft, and S/N PC (with the value 6RB7L22). Below the fields is a privacy policy notice: "I accept VELP SCIENTIFICA privacy policy (for further details, please look at the following link: www.velp.com/en/privacy)". There are radio buttons for "I accept" and "I don't accept". A legend indicates that red asterisks mark obligatory fields. At the bottom, there are "Save" and "Cancel" buttons, and a red instruction: "Copy and send as attached to service@velp.it".

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10005112 Cavo USB A/Micro B 1.5m

Thank you for having chosen VELP!

Established in 1983, VELP is today one of the world's leading manufacturer of analytical instruments and laboratory equipment that has made an impact on the world-wide market with Italian products renowned for innovation, design and premium connectivity. VELP works according to **ISO 9001**, **ISO14001** and **ISO 45001** Quality System Certification.

Our instruments are manufactured in Italy according to the IEC 1010-1 and CE regulation.

Our product lines:

Analytical instruments

Elemental Analyzers
Digestion Units
Distillation Units
Solvent Extractors
Fiber Analyzers
Dietary Fiber Analyzers
Oxidation Stability Reactor
Consumables

Laboratory Equipment

Magnetic Stirrers
Heating Magnetic Stirrers
Heating Plates
Overhead stirrers
Vortex mixers
Dispersers
COD Thermoreactors
BOD and Respirometers
Cooled Incubators
Flocculators
Overhead Shakers
Turbidimeter
Open Circulating Baths
Pumps

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VELP Scientifica Srl