Instruction Manual
Manuale di istruzioni
Manuel d'instructions
Manual de instrucciones
Bedienungsanleitung
指导手册

OHS 200 DIGITAL Overhead stirrers

F20100490, F20110490

OHS 100 DIGITAL Overhead stirrers

F20100491, F20110491

OHS 60 DIGITAL Overhead stirrers

F20100492, F20110492

OHS 40 DIGITAL Overhead stirrers

F20100493, F20110493

OHS 20 DIGITAL Overhead stirrers

F20100494, F20110494

General Information / Informazioni Generali / Informations Générales / Información General / Allgemeine Hinweise / 一般信息

Before using the unit, please read the following instruction manual carefully.

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Do not dispose of this equipment as urban waste, in accordance with EEC directive 2002/96/CE.

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This unit must be used for laboratory applications indoor only. The manufacturer declines all responsibility for any use of the unit that does not comply with these instructions. If the product is used in a not specified way by the manufacturer or with not specified accessories, product's safety may be compromised.

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IEC/EN 61010-1 IEC/E 61010-2-051

Electrical equipment for laboratory use
General requirement - Canadian electrical code

UL 61010-1
CAN/CSA-C22.2 No.61010-1

VELP reserves the right to modify the characteristics of its products with the aim to constantly improving their quality.
Nell'impegno di migliorare costantemente la qualità dei prodotti, VELP si riserva la facoltà di variane le caratteristiche.
Dans le but d'améliorer constamment la qualité de ses produits, VELP se réserve le droit d'apporter des modifications aux caractéristiques de ceux-ci.
VELP se reserva el derecho de modificar las características de sus productos con el objetivo de mejorar constantemente su calidad.
VELP behält sich zum Zwecke der ständigen Verbesserung der Produktqualität das Recht auf Änderung der Geräteeigenschaften vor.
VELP 保留修改其产品特性的权利，目的是不断提高其质量。

Safety Regulations / Norme di Sicurezza / Consignes de Sécurité / Advertencias de Seguridad / Sicherheitshinweise / 安全条例

The plug disconnects the instrument. Therefore, place the instrument where it can be quickly disconnected. / La spina è il mezzo di disconnessione dell'apparecchio. Pertanto, non posizionare l'apparecchio in modo che sia difficile azionare il mezzo di disconnessione. / Le bouchon est le moyen de déconnexion de l'appareil. Par conséquent, placer l'appareil où il peut être rapidement débranché. / El tapón es el medio de desconexión del dispositivo. No coloque el dispositivo en una forma que es difícil de desconectar. / Der Stecker trennt das Gerät. Daher Stellen Sie das Instrument, wo es schnell getrennt werden kann. / 该插头可以断开仪器的连接。因此，要把仪器放在可以快速断开的地方。

The values indicated on the rating plate of the instrument must correspond to those of the power supply.
I valori di tensione indicato sulla targhetta del modello e quello di rete devono coincidere.
Les valeurs indiquées sur la plaque signalétique de l'appareil doivent correspondre à ceux de l'alimentation.
Los valores de tensión indicados en la placa y que de la red debe ser los mismos.
Die angegebene Spannung Wert auf dem Typenschild und das Netzwerk muss gleich sein.
仪表的铭牌上标明的数值必须与电源的数值一致。

Position the instrument on a flat surface, with a distance from the wall of 30 cm (at least).
Posizionare lo strumento su superfici piane, ad una distanza dalle pareti di almeno 30 cm.
Positionner l'appareil sur une surface plat, avec une distance de la paroi de 30 cm (au moins).
Coloque la unidad sobre una superficie plana, con una distancia de la pared de 30 cm (por lo menos).
Stellen Sie das Gerät auf einer ebenen Fläche mit einem Abstand zur Wand von 30 cm (mindestens).
将仪器放置在一个平坦的表面上，与墙壁的距离为 30 厘米（至少）。

Fasten the unit to the support rod (A00000366 or A00000369) using the double clamp (A00001301). Secure the receptacle using the ribbon clamp (A00001302).
Fissare saldamente lo strumento allo stativo (A00000366 o A00000369) mediante il morsetto doppio (A00001301) e il recipiente di agitazione con l'apposita cinghia (A00001302).
Fixez l'unité à le statif (A00000366 ou A00000369) con la noix de fixation double (A00001301). Fixer le réceptacle à l'attache souple avec ruban (A00001302).
Asegurar firmemente la herramienta a la barra de soporte (A00000366 o A00000369) con la abrazadera doble (A00001301) y el recipiente con la abrazadera de cinta (A00001302).
Befestigen Sie das Gerät an der H-Stativ (A00000366 oder A00000369) mit der Kreuzmuffe (A00001301). Sichern Sie die Aufnahme mit der Spannhalter (A00001302).
使用双夹子（A00001301）将仪器固定在支撑杆（A00000366 或 A00000369）上。用带状夹子（A00001302）固定插座。

Safe working conditions are ensured only when the accessories described in the dedicated chapter are used.
Il funzionamento sicuro è garantito soltanto con gli accessori descritti nel relativo capitolo.
Conditions de travail sûres sont assurées que lorsque les accessoires décrits dans le chapitre dédié sont utilisés.

El funcionamiento es seguro sólo con los accesorios descritos en el capítulo correspondiente.
Sichere Arbeitsbedingungen sind gewährleistet, wenn das Zubehör in dem Kapitel beschrieben verwendet werden.
只有在**使用**专用章节中描述的附件时，才能确保安全的工作条件。

The working speed set on the instrument must be such as to avoid wobbling and/or splashes.
Il numero di giri impostato deve escludere eventuali squilibri dell'agitatore e possibili spruzzi del prodotto agitato.
Le nombre de tours de l'ensemble agitateur doit exclure les déséquilibres et les éclaboussures du produit agité.
El número de revoluciones del agitador debe excluir cualquier desequilibrio y posible de salpicar de el producto agitado.
Die Arbeitsgeschwindigkeit des Gerätes muss gesetzt sein, wie Wackeln und / oder Spritzer zu vermeiden.
仪器上设定的工作速度必须能够避免摇晃和/或飞溅。

Do not use with explosive or dangerous materials for which the equipment is not designed. The stirrer must not be used in explosive atmospheres, in bain-marie or to stir harmful liquids prior using protective measures according to the safety standards of the processed products and/or in force in the laboratories including personal protective equipment and the presence of an extraction hood which ensures at least 10-fold air change in accordance with the standards EN 14175 and DIN 12924.

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It is dangerous to run the unit with the stirring blade turning in free air. Always place the stirring shaft in the receptacle before turning the unit on.

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It is responsibility of the user appropriately decontaminate the instrument in case of dangerous substances fall on or in it.

It is also responsibility of the user to use safety substances for cleaning or decontaminating, which do not react with internal parts of the instrument or with the material contained in it. In case of doubts on the compatibility of a cleaning solution, contact the manufacturer or local distributor.

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The solution may release toxic, dangerous or poisonous gases. Adequate safety measures must be taken, in accordance with the safety regulations in force, including the presence of hood and personal protective equipment (masks, gloves, goggles, etc.).

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Switch off the the stirring before opening chuck locking ring. Switch off the instrument before removing chuck.

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OHS overhead stirrers with electronic speed control, brushless motor, and advanced safety features are able to satisfy the most difficult laboratory applications in terms of viscosity and volume. The new chuck ensures higher safety for the operator and allows to use a passing rod of up to 8.5mm in diameter.

Figure 2. OHS 200 Digital

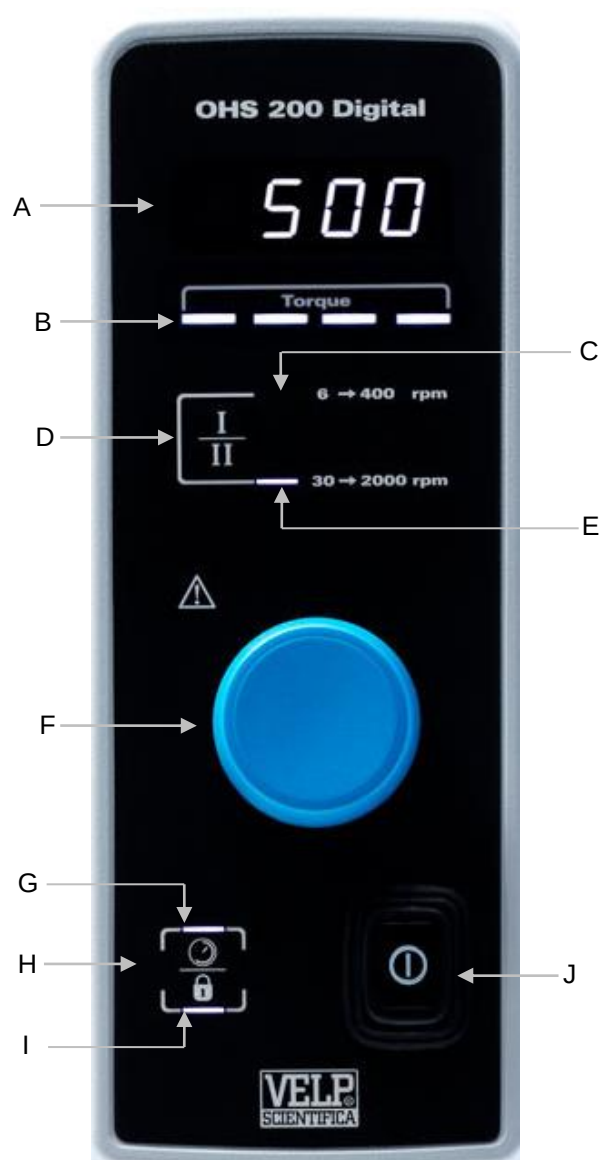
Figure 1. OHS 200 Digital - Frontal view

- A 7 Segment Display LED
- B Torque indicator bar
- C Low speed LED (only OHS 200)
- D Gear key (only OHS 200) – Timer key for other models
- E High speed LED (only OHS 200) – Timer LED for other models

- F Speed control knob
- G Timer LED (for OHS 200) - Upper lock LED for other models
- H Timer key (for OHS 200) – Lock key for all models
- I Lower Lock LED
- J Main switch
- K Smart chuck

2. Assembly and installation

- Unpacking
 - Check the integrity of the unit after unpacking.
- The box includes
 - OHS (XXX) digital overhead stirrer
 - Power supply cable
 - Instruction manual
 - Allen key and allen head screw
 - Shaft support overhead stirrer
- Installation
 - Place the unit on non-flammable surface
 - Fix the shaft support to the instrument with the allen head screw
 - Fasten the unit to the support rod (A00000366 or A00000369) using the double clamp (A00001301)
 - Secure the receptacle using the ribbon clamp (A00001302)
 - Slide the stirring shaft into the chuck and tight it
 - Make sure that the rating value of the instrument corresponds to the one of the power supply
 - Ensure that the socket provided, with grounding, is compliant with the current safety norms and easy to reach.

3. Working

Commissioning	➤ Switch on the instrument through the main switch (J) ➤ Display (A) shows software version, last set point value and OFF (if "Mode" is set on Stop, see chapter 0).
Stirring	➤ Set the operating speed range by clicking Gear key (D) (Only OHS 200). ➤ Adjust speed set point by turning the speed control knob (F). ➤ Click the knob to start stirring. ➤ Speed increases until set point achievement. ➤ A microprocessor ensures constant speed even when the viscosity changes (counter-reaction). ➤ Switch off the stirring by clicking the knob.
Timer	➤ Click Timer key (D or H). HH:MM appears on display. ➤ Rotate speed control knob (F) to set the desired value. Click to confirm. ➤ Once timer is set, timer LED lights up: - (G) for OHS 200 - (E) for other models ➤ The LED lights continuously until the instrument starts to stir. When the instrument stirs, the LED flashes. ➤ It's possible to visualize (or modify) countdown in each moment during the analysis by clicking timer key. ➤ When countdown finishes, "End" is visualized on the display.
Gear (only for OHS 200)	➤ Set the operating speed range by clicking Gear button (D). ➤ The gear set has its relative LED (C or E) lighted on. ➤ It's possible to select between two gears: - Gear I: Low speed (6 – 400 rpm) High torque (till 200 Ncm). - Gear II: High speed (30 – 2000 rpm) Low torque (till 40 Ncm).
Torque	➤ The torque indicator bar (B), made by four LEDs, shows the intensity of the torque applied
Lock	➤ Holding the Lock key (H) for 3 seconds it's possible to lock the control panel. ➤ Unlock the control panel by holding key H for 3 seconds. ➤ When the instrument is locked, - LED (I) lights on for OHS 200 - LED (G) and (I) light on for other models

4. External Connections

USB	➤ Each model has USB connection for software upgrading by PC. Software version, dedicated program to be installed on the PC and software installation guide must be requested by e-mail to service@velp.it
------------	--

5. Menu

Hold the control knob (F) for 3 seconds to enter into the setting mode when stirring is off.

Turning the control knob (F) it is possible to pass from one of the following parameters to the next one. Click the knob to modify:

Parameter shown on display	Default value		Range	Description
NOdE	StOP		StOP - run	Set starting mode: ➤ Stop: when the instrument is switched on through the main switch, OFF is on display ➤ Run: when the instrument is switched on, it restarts to work with the last set points set
LINT	OHS 200	400 (I) 2000 (II)	100 – 400 100 - 2000	Set the maximum value of speed set point speed
	OHS 100	1300	100 – 1300	
	OHS 60	2000	100 – 2000	
	OHS 40	2000	100 - 2000	
	OHS 20	2000	100 - 2000	
COun				Running time in H. If “.” Is showed before the number, the running time is in Day.
UPGr				To install a new software version
rSEt	nO		YES - nO	To restore default setting parameters
CAI	nO		YES - nO	To reset torque zero value

To leave settings mode wait 10 seconds without touching any knob or click the lock key.

6. SmartChuck™

Smart chuck (K) allows to fix stirring shaft with one hand.

- Open the chuck: From work position (Figure 3), turn Locking Ring 90° to the left (Figure 4). Pull Locking Ring down (Figure 5).
- Close the chuck: Push Locking Ring up (Figure 6) and turn it 90° to the right.

Figure 3. Working position

Figure 4. Intermediate position

Figure 5. Open position

Figure 6. Intermediate position

7. Error messages

When the display shows an error message, the stirring function stops automatically.

Error code	Cause
AL1	Motor doesn't start stirring
AL2	High internal motor temperature
AL3	Motor overload
AL4	High driver temperature
AL5	Safety relay fault

To remove the error message, disconnect the instrument from the power supply.

If alarm persists on the display, please contact VELP Scientifica's technical service department. service@velp.it

8. Maintenance

Maintenance	➤ No routine or extraordinary maintenance is necessary. ➤ Repairs must be carried out by authorized Velp personnel only. ➤ Instrument must be transported in its original packaging any indications present on the original packaging must be followed (e.g. palletized). ➤ It is the responsibility of the user, to properly decontaminate the unit in case of hazardous substances remaining on the surface or interior of the device. If in doubt about the compatibility of a cleaning or decontamination product, contact the manufacturer or distributor.
Cleaning	➤ Disconnect the unit from the power supply and use a cloth dampened with a non-flammable non-aggressive detergent.

9. Technical data

	Models		F20100490, F20100491, F20100492, F20100493	F20110490, F20110491, F20110492, F20110493
General features	Power supply		230 V / 50-60 Hz (+/-10%)	115V / 60 Hz (+/-10%)
	Dimensions (WxHxD)	OHS200	90x315x235 mm (3.54x12.40x9.25 in)	90x315x235 mm (3.54x12.40x9.25 in)
		OHS 100 – OHS60 – OHS40 – OHS20	90x285x235 mm (3.54x11.22x9.25 in)	90x285x235 mm (3.54x11.22x9.25 in)
	Weight	OHS200	4,6 kg (10,14 lb)	4,6 kg (10,14 lb)
		OHS100	4,1 kg (9,04 lb)	4,1 kg (9,04 lb)
		OHS60 – OHS40 – OHS20	4,1 kg (9,04 lb)	4,1 kg (9,04 lb)
	Power input		190 W	190 W
	Construction material (structure)		Aluminum	Aluminum
	Working in continuous		Admitted	Admitted
	Settable restart modality		Stop or work	Stop or work
	Noisiness		<< 60 dBa	<< 60 dBa
	Environmental temperature admitted		+5...+40 °C	+5...+40 °C
	Storage temperature admitted		-10...+60 °C	-10...+60 °C
	Max humidity		80%	80%
	Level of electrical protection CEI EN60529		IP 54	IP 54
	Overvoltage category		II	II
	Pollution degree CEI EN61010-1		2	2
	Max altitude		2000 m	2000 m
Stir	Stirring capacity	OHS200	100 l H ₂ O	100 l H ₂ O
		OHS100	100 l H ₂ O	100 l H ₂ O
		OHS60	40 l H ₂ O	40 l H ₂ O
		OHS40	25 l H ₂ O	25 l H ₂ O
		OHS20	25 l H ₂ O	25 l H ₂ O
	Programmable speed range	OHS200	6-400rpm (I) – 30-2000rpm (II)	6-400rpm (I) – 30-2000rpm (II)
		OHS100	30-1300rpm	30-1300rpm
		OHS60	30-2000rpm	30-2000rpm
		OHS40	30-2000rpm	30-2000rpm
		OHS20	30-2000rpm	30-2000rpm
	Motor type		BLDC	BLDC
	Speed selection		1 rpm step	1 rpm step
	Stirring alarm		Motor fault	Motor fault
	Motor rating output		150 W	150 W
Torque	Max torque admitted	OHS200	200 Ncm (I) – 40 Ncm (II)	200 Ncm (I) – 40 Ncm (II)
		OHS100	100 Ncm	100 Ncm
		OHS60	60 Ncm	60 Ncm
		OHS40	40 Ncm	40 Ncm
		OHS20	20 Ncm	20 Ncm
Counters	Motor counter		Working hours	Working hours

10. Accessories / Spare parts

A00000366	Support rod and base OHS	A00001307	Stirring shaft with propeller
A00000369	Telescopic support rod and base OHS	A00001308	Stirring shaft with paddle, 6 holes
A00000391	OHS/ControllerSoft	A00001309	Stirring shaft with turbine
A00001300	Support rod and base	A00001310	Stirring shaft with turbo propeller
A00001301	Double clamp	A00001311	Stirring shaft with anchor
A00001302	Ribbon clamp	A00001312	Chuck guard OHS
A00001304	Stirring shaft with floating blade		
A00001305	Stirring shaft with folding blade	10007175	Knob 35D blu
A00001306	Stirring shaft with fixed blade		

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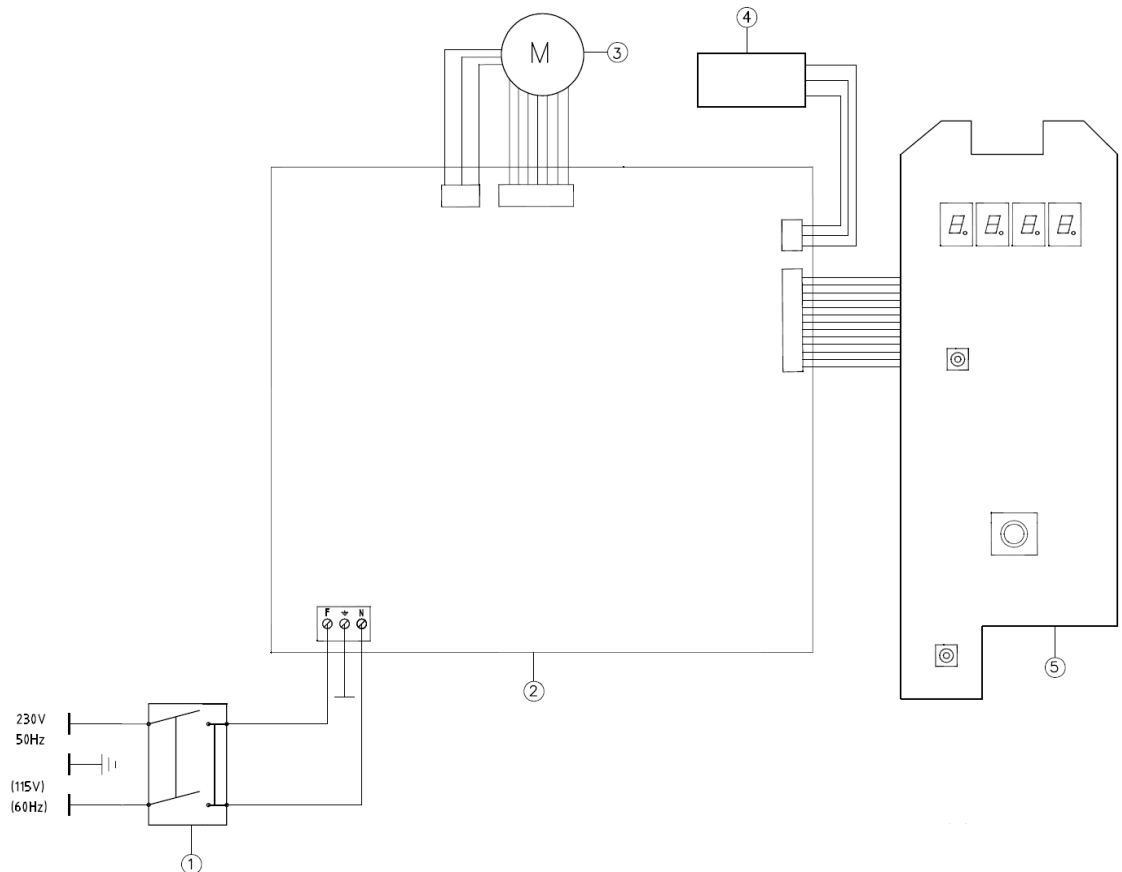
We, the manufacturer VELP Scientifica, under our responsibility declare that the product is manufactured in conformity with the following standards:

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EN 61010-1	Safety Requirements for Electrical Equipment for Measurement, Control, and Laboratory Use - Part 1: General Requirements
EN 61010-2-051	Safety requirements for electrical equipment for measurement, control and laboratory use - Part 2-051: Particular requirements for laboratory equipment for mixing and stirring
EN 61326-1	Electrical equipment for measurement, control and laboratory use - EMC requirements -
2015/863/EU (RoHS III)	Restriction of the use of certain hazardous substances
2012/19/EU (WEEE)	On waste electrical and electronic equipment

and satisfies the essential requirements of the following directives: / e soddisfa i requisiti essenziali delle direttive: / et qu'il satisfait les exigences essentielles des directives: / y cumple con los requisitos esenciales de las directivas: / und den Anforderungen folgender Richtlinien entspricht: / 并满足以下指令的基本要求。

- Machinery directive 2006/42/EC / Macchine 2006/42/CE / Machines 2006/42/CE / Máquinas 2006/42/CE / Maschinen 2006/42/EG / 机械指令 2006/42/EC
- Electromagnetic compatibility directive 2014/30/EU / Compatibilità elettromagnetica 2014/30/EU / Compatibilité électromagnétique 2014/30/EU / Compatibilidad electromagnética 2014/30/EU / Elektromagnetische Verträglichkeit 2014/30/EU / 电磁兼容性指令 2014/30/EU
- plus modifications / più modifiche / plus modifications / más sucesivas modificaciones / in der jeweils gültigen Fassung / 加上变化

1. Declaration of conformity

We, the manufacturer VELP Scientifica, under our responsibility declare that the product is manufactured in conformity with the following regulations:

- | | |
|----------------|--|
| S.I. 2016/1101 | Electrical Equipment (Safety) Regulations 2016 |
| S.I. 2016/1091 | Electromagnetic Compatibility Regulations 2016 |

according to the relevant designated standards:

- | | |
|----------------|---|
| EN 61010-1 | Safety requirements for electrical equipment for measurement, control, and laboratory use
Part 1: General requirements |
| EN 61010-2-051 | Particular requirements for laboratory equipment for mixing and stirring |
| EN 61326-1 | Electrical equipment for measurement, control and laboratory use - EMC requirements -
Part 1: General requirements |

and satisfies the essential requirements of regulations:

- | | |
|----------------|---|
| S.I. 2008/1597 | Supply of Machinery (Safety) Regulations 2008 |
| S.I. 2012/3032 | Restriction of the Use of Certain Hazardous Substances in Electrical and Electronic Equipment
Regulations 2012 |
| S.I. 2013/3113 | Waste Electrical and Electronic Equipment Regulations 2013 |

plus modifications.

Thank you for having chosen VELP!

Established in 1983, VELP is today one of the world's leading manufacturer of analytical instruments and laboratory equipment that has made an impact on the world-wide market with Italian products renowned for innovation, design and premium connectivity. VELP works according to **ISO 9001**, **ISO14001** and **OHSAS 18001** Quality System Certification.

Our instruments are manufactured in Italy according to the IEC 1010-1 and CE regulation.

Our product lines:

Analytical instruments

Elemental Analyzers
Digestion Units
Distillation Units
Solvent Extractors
Fiber Analyzers
Dietary Fiber Analyzers
Oxidation Stability Reactor
Consumables

Laboratory Equipment

Magnetic Stirrers
Heating Magnetic Stirrers
Heating Plates
Overhead stirrers
Vortex mixers
Homogenizers
COD Thermoreactors
BOD and Respirometers
Cooled Incubators
Flocculators
Overhead Shakers
Turbidimeter
Radiation Detector
Open Circulating Baths
Pumps

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VELP Scientifica Srl