Instruction Manual
Manuale di istruzioni
Manuel d'instructions
Manual de instrucciones
Bedienungsanleitung
指导手册

PW - LH - LS - ES Overhead stirrers

F201A0150, F201A0156, F201A0151, F201A0152

**General Information / Informazioni Generali / Informations Générales / Información General /
Allgemeine Hinweise / 一般信息**

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This unit has been designed and manufactured in compliance with the following standards:

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Electrical equipment for laboratory use

UL 61010-1

General requirement - Canadian electrical code

CAN/CSA-C22.2 No.61010-1

VELP reserves the right to modify the characteristics of its products with the aim to constantly improving their quality.

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The values indicated on the rating plate of the instrument must correspond to those of the power supply.

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Position the instrument on a flat surface, with a distance from the wall of 30 cm (at least).

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Safe working conditions are ensured only when the accessories described in the dedicated chapter are used.

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The working speed set on the instrument must be such as to avoid wobbling and/or splashes.

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Do not use the instrument in explosive environment or with dangerous or flammable substances. The unit must not be immersed in water.

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It is dangerous to run the unit with the stirring blade turning in free air. Always place the stirring shaft in the receptacle before turning the unit on.

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The unit is fitted with two fuses (2xT1,6 A L 250 V – 2xT800 mA x ES), found in the socket on the back. To replace one or more disconnect the mains cable and, using a screwdriver, lift up the small cover on the fuse box.

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It is responsibility of the user appropriately decontaminate the instrument in case of dangerous substances fall on or in it.

It is also responsibility of the user to use safety substances for cleaning or decontaminating, which do not react with internal parts of the instrument or with the material contained in it. In case of doubts on the compatibility of a cleaning solution, contact the manufacturer or local distributor.

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The solution may release toxic, dangerous or poisonous gases. Adequate safety measures must be taken, in accordance with the safety regulations in force, including the presence of hood and personal protective equipment (masks, gloves, goggles, etc.).

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Overhead stirrers are generally used for laboratory applications where the solubilization, mixing, suspension, emulsifying and homogenization of samples is required.

The PW, LH, LS and ES overhead stirrers with electronic speed control and advanced safety features are able to satisfy the most diverse needs of the laboratory in terms of both viscosity and volume.

The chuck is keyless and ensures a secure and easy-to-fit mount for stirring shafts of up to 10mm in diameter. The stirring speed is electronically controlled. The shaft bearings and belt drive ensure silent running.

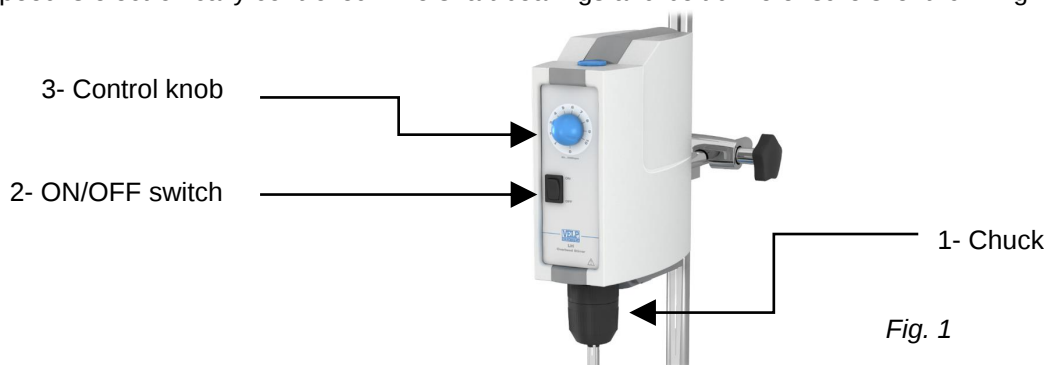
Fig. 1

2. Assembly and installation

Check the integrity of the unit after unpacking. The box includes:

- Overhead stirrer
- Instruction manual
- Allen key
- Allen head screw
- Shaft
- Power supply cable

2.1 Electrical connection

After unpacking the instrument, place the unit on the laboratory bench. Before connecting the instrument to the power supply, make sure that the values on the rating plate correspond to those of the power supply.

Ensure that the socket and the relative cut-off device conform to current safety norms and are easy to reach.

2.2 Start-up

Assemble the shaft using the Allen head screw. Fasten the instrument securely to the the H-Stand (A00000366 or A00000369) with the double clamp (A00001301). Secure the receptacle using the ribbon clamp (A00001302) to avoid damage.

2.3 Information about construction materials

Housing	Die casting Polymer	Structure	Aluminum	Control panel	PET
---------	---------------------	-----------	----------	---------------	-----

3. Operating controls

Turn the unit on using the mains switch. Turn the potentiometer knob to be select the stirring speed. Stirring speeds are indicated on the analogue scale around the knob. **NOTE:** The speed setting is saved when the instrument is turned off or if there is a power failure. When the instrument is turned on again, stirring starts at the previously set speed.

3.1 Overload protection

The current and voltage are electronically limited in order to safeguard the motor and the electronic parts against electric surges.

In the event of abnormal conditions arising the unit shuts down automatically.

Over Load Current surges

The stirrer can compensate for eventual current surges for a limited period of time.

If the situation persists, the speed of the motor is gradually reduced, when a speed of 0 is reached the unit shuts down.

Over Curr Rpm error

This error occurs if the chuck is accidentally blocked or if the number of revs exceeds the max. permitted value.

Over Temp Temperature error

This error occurs if the unit is used for heavy-duty work for prolonged periods of time and at an ambient temperature that exceeds the max. permitted temperature.

Before resetting the instrument allow it to cool down.

Over Volt Voltage surge

The software limits the power supply of the motor to the nominal value of the motor itself (60V).

A high speed setting or a heavy load on the chuck may prompt a higher voltage. In this case the voltage is limited to the nominal value of the motor (60 Volt) and the number of revs is reduced accordingly.

If for any reason the power supply of the motor exceeds the maximum threshold value of 70V, "OVERLOAD" the unit will shut down.

NOTE: To reset the unit remove the load from the chuck and turn the unit off and on again using the mains switch.

4. Maintenance

No routine or extraordinary maintenance is necessary apart from periodically cleaning the unit as described in this manual. In compliance with the product guarantee law, repairs to our units must be carried out in our factory, unless previously agreed otherwise with local distributors. The instrument must be transported in its original packaging and any indications present on the original packaging must be followed (e.g. palletized).

It is the responsibility of the user, to properly decontaminate the unit in case of hazardous substances remaining on the surface or interior of the device. If in doubt about the compatibility of a cleaning or decontamination product, contact the manufacturer or distributor.

4.1 Cleaning

Disconnect the unit from the power supply and use a cloth dampened with a mild, non-flammable detergent.

5. Technical data

Code Model	F201A0150 PW	F201A0156 LH	F201A0151 LS	F201A0152 ES
Admitted power supply	110-230V 50/60Hz(+/-10%)	110-230V 50/60Hz(+/-10%)	110-230V 50/60Hz(+/-10%)	110-230V 50/60Hz(+/-10%)
Max. input / output power	190 W / 435 W	190 W / 280 W	120 W / 180 W	30 W / 45 W
Dimensions (WxHxD)	80x230x196 mm (3.1x9.0x7.7 in)	80x230x196 mm (3.1x9.0x7.7 in)	80x215x196 mm (3.1x8.5x7.7 in)	80x160x200 mm (3.1x6.3x7.9 in)
Weight	2,9 Kg (6,4 lbs)	2,9 Kg (6,4 lbs)	2,3 Kg (5 lbs)	1,3 Kg (2,86 lbs)
Speed range at nominal load	20 – 1200rpm	50 – 2000rpm	50 – 2000rpm	50 – 1300rpm
Speed indicator	Analogue	Analogue	Analogue	Analogue
Max torque stirrer shaft	120Ncm	80Ncm	40Ncm	15Ncm
Max stirring quantity H ₂ O	70 litres	40 litres	25 litres	15 litres
Clamping chuck clamping range	From 1 to 10mm	From 1 to 10mm	From 1 to 10mm	From 1 to 10mm
Overload and overtemperature protection	Automatic cut-off	Automatic cut-off	Automatic cut-off	Automatic cut-off
Enviromental ambient temperature	5 - 40°C (41 - 104°F)	5 - 40°C (41 - 104°F)	5 - 40°C (41 - 104°F)	5 - 40°C (41 - 104°F)
Storage temperature range	-10... + 60°C (14 - 140°F)	-10... + 60°C (14 - 140°F)	-10... + 60°C (14 - 140°F)	-10... + 60°C (14 - 140°F)
Max humidity	80%	80%	80%	80%
Noise Level	<<80dBa	<<80dBa	<<80dBa	<<80dBa
Operating mode	Continuous	Continuous	Continuous	Continuous
Overvoltage category CEI-EN61010-1	II	II	II	II
Pollution degree CEI-EN61010-1	2	2	2	2
Max altitude installation	2000m (6561 ft) asl	2000m (6561 ft) asl	2000m (6561 ft) asl	2000m (6561 ft) asl
Rapid Fuse diam. 5x20mm	2xF1,6A L 250V	2xF1,6A L 250V	2xF1,6A L 250V	2XF800mA L 250V

6. Accessories / Spare parts

A00000369	Telescopic H-Stand
A00000366	Universal H-Stand
A00001301	Double clamp
A00001302	Ribbon clamp
A00001304	Stirring shaft with floating blade
A00001305	Stirring shaft with folding blade
A00001306	Stirring shaft with fixed blade
A00001307	Stirring shaft with propeller
A00001308	Stirring shaft with paddle, 6 holes
A00001309	Stirring shaft with turbine
A00001310	Stirring shaft with turbo propeller
A00001311	Stirring shaft with anchor

			
A00000369	A00000366	A00001302	A00001301
			
A00001305	A00001306	A00001307	A00001308
			
A00001304	A00001309	A00001310	A00001311

10005266	Knob 26D blue
10000132	Gasket for overhead stirrer

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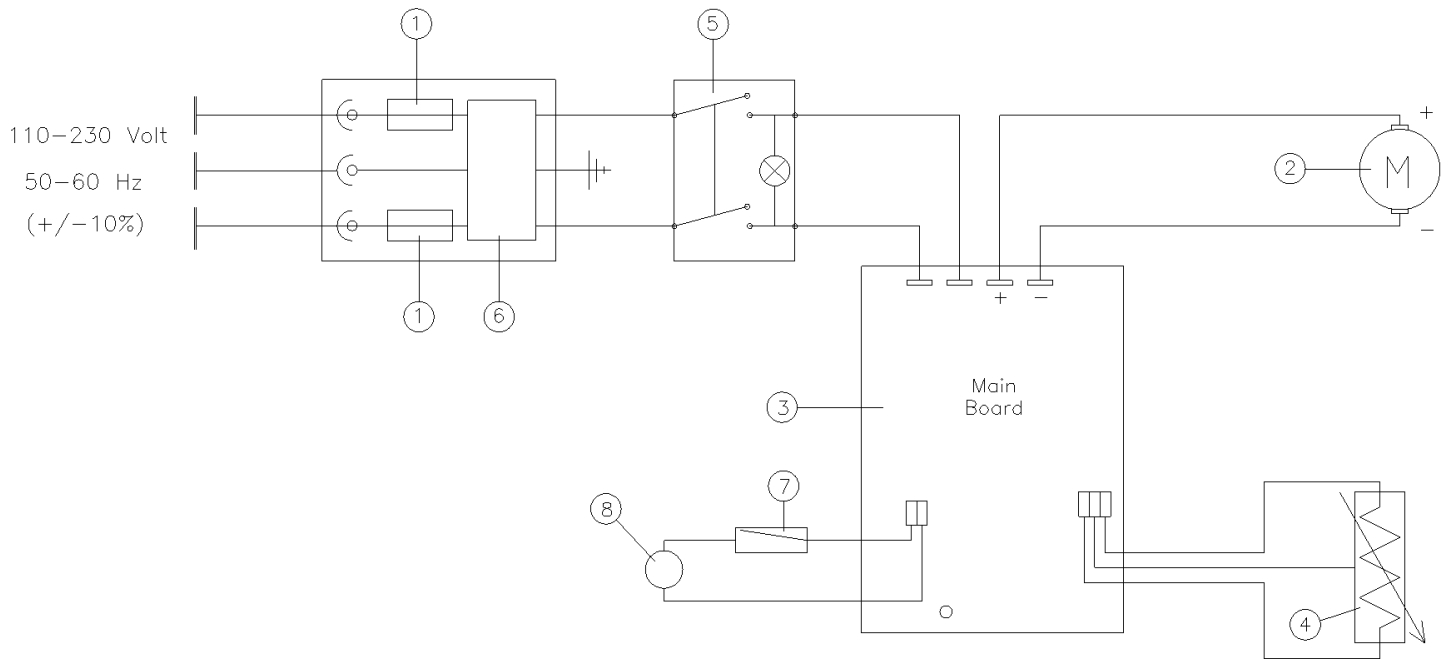
We, the manufacturer VELP Scientifica, under our responsibility declare that the product is manufactured in conformity with the following standards:

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EN 61010-2-051

EN 61326-1

and satisfies the essential requirements of the following directives:

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9. Declaration of conformity

We, the manufacturer VELP Scientifica, under our responsibility declare that the product is manufactured in conformity with the following regulations:

S.I. 2016/1101 Electrical Equipment (Safety) Regulations 2016

S.I. 2016/1091 Electromagnetic Compatibility Regulations 2016

according to the relevant designated standards:

EN 61010-1 Safety requirements for electrical equipment for measurement, control, and laboratory use
Part 1: General requirements

EN 61010-2-051 Particular requirements for laboratory equipment for mixing and stirring

EN 61326-1 Electrical equipment for measurement, control and laboratory use - EMC requirements -
Part 1: General requirements

and satisfies the essential requirements of regulations:

S.I. 2008/1597 Supply of Machinery (Safety) Regulations 2008

S.I. 2012/3032 Restriction of the Use of Certain Hazardous Substances in Electrical and Electronic
Equipment Regulations 2012

S.I. 2013/3113 Waste Electrical and Electronic Equipment Regulations 2013

plus

modifications.

Thank you for having chosen VELP!

Established in 1983, VELP is today one of the world's leading manufacturer of analytical instruments and laboratory equipment that has made an impact on the world-wide market with Italian products renowned for innovation, design and premium connectivity. VELP works according to **ISO 9001**, **ISO14001** and **OHSAS 18001** Quality System Certification.

Our instruments are manufactured in Italy according to the IEC 1010-1 and CE regulation.

Our product lines:

Analytical instruments

Elemental Analyzers
Digestion Units
Distillation Units
Solvent Extractors
Fiber Analyzers
Dietary Fiber Analyzers
Oxidation Stability Reactor
Consumables

Laboratory Equipment

Magnetic Stirrers
Heating Magnetic Stirrers
Heating Plates
Overhead stirrers
Vortex mixers
Homogenizers
COD Thermoreactors
BOD and Respirometers
Cooled Incubators
Flocculators
Overhead Shakers
Turbidimeter
Radiation Detector
Open Circulating Baths
Pumps

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