Instruction Manual
Manuale di istruzioni
Manuel d'instructions
Manual de instrucciones
Bedienungsanleitung
指导手册

ROTAX Overhead Mixer

F10600118, F10610118

General Information / Informazioni Generali / Informations Générales / Información General / Allgemeine Hinweise / 一般信息

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This unit has been designed and manufactured in compliance with the following standards:

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Safety requirements for electrical equipment for measurement, control and for laboratory use

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**IEC/EN 61010-1
IEC/EN 61010-2-051**

Electrical equipment for laboratory use

UL 61010-1

General requirement - Canadian electrical code

CAN/CSA-C22.2 No.61010-1

VELP reserves the right to modify the characteristics of its products with the aim to constantly improving their quality.

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The plug disconnects the instrument. Therefore, place the instrument where it can be quickly disconnected.

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The stirred solution may release toxic, dangerous or poisonous gases. Adequate safety measures must be taken, in accordance with the safety regulations in force, including the presence of hood and personal protective equipment (masks, gloves, goggles, etc.).

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Position the instrument on a flat surface, with a distance from the wall of 30 cm (at least).

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Do not use the instrument in explosive environment or with dangerous or flammable substances. The unit must not be immersed in water.

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It is responsibility of the user appropriately decontaminate the instrument in case of dangerous substances fall on or in it.

It is also responsibility of the user to use safety substances for cleaning or decontaminating, which do not react with internal parts of the instrument or with the material contained in it. In case of doubts on the compatibility of a cleaning solution, contact the manufacturer or local distributor.

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Safe working conditions are ensured only when the accessories described in the dedicated chapter are used.

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The German standard methods for the examination of water, waste water and sludge require the use of overhead mixing for the determination of the leachability by water of sludge, sediments and solid wastes (DIN 38414, October 1984, Group S, Sections 4). The equipment can be used in all the applications where rotating, shaking and mixing are required.

Summary of the DIN 38414 - Equipment and products

Overhead mixer, glass or plastic bottles 2000 ml, set for membrane filtration, conductimeter, pH-meter and distilled water. 100 g of solid material (dry weight) with dimensions less than 10 mm are introduced with 1 l of distilled water into 2 l glass or plastic vessels. The vessels are overturned at low speed 5-6 rpm. at room temperature during 24 hours. The eluate after membrane filtration (0.45 mm) is used for the determination of soluble pollutants.

2. Assembly and installation

Check the integrity of the unit after unpacking. The box includes:

- ROTAX
- Power supply cable
- Instruction manual

2.1 Electrical connection

Before connecting the instrument to the power supply, make sure that the values on the rating plate correspond to those of the power supply. Connect the unit to the power supply using the transformer supplied.

Ensure that the socket and the relative cut-off device conform to current safety norms and easy to reach.

2.2 Start-up

The bottles are placed first on one side platform and fixed by pushing the spring knob and fixing it by the screw clamp.

The instrument is then rotated 180° by turning on the main switch and the second group of bottles can be placed into position and fixed. To increase the life span of the motor always distribute the weight of the bottles evenly.

2.3 Arm replacement

The use with 1 l bottles is possible by the substitution of the arm holding bottles 220-280 mm high (2 l) with the complementary A00001023. The substitution is performed unscrewing the four screws fixing the normally mounted arm and replacing it with the new arm fixed by the same four screws but at the lower row. Repeat the substitution on the opposite side in order to maintain equilibrated the rotating masses.

2.4 Safety devices and warning signals

The instrument is equipped with a safety device that cuts-off the power supply to the motor when the rotating part encounters an obstacle; a warning message "AL" appears on the display. To return to normal running mode turn the instrument off and remove the obstacle before turning it on again. Should a warning message appear on the display when no obstacles are present, check that the weight of the bottles is evenly distributed between the two brackets. Always make sure that the mains switch is turned to "0" when loading and unloading bottles.

3. Operating controls

The equipment can hold up to six 2 l glass or plastic vessels, 220-280 mm high or with an optional clamp up to 8 vessels 200-270 mm high. The motor is directly connected to the rotating shaft by a gear speed reducer. The speed of rotation

controlled by the electronic card and shown on the display can be varied continuously from 2 to 30 rpm using  and .

4. Maintenance

No routine or extraordinary maintenance is necessary apart from periodically cleaning the unit as described in this manual. In compliance with the product guarantee law, repairs to our units must be carried out in our factory, unless previously agreed otherwise with local distributors. The instrument must be transported in its original packaging and any indications present on the original packaging must be followed (e.g. palletized).

It is the responsibility of the user, to properly decontaminate the unit in case of hazardous substances remaining on the surface or interior of the device. If in doubt about the compatibility of a cleaning or decontamination product, contact the manufacturer or distributor.

4.1 Cleaning

Disconnect the unit from the power supply and use a cloth dampened with a non-inflammable non-aggressive detergent.

5. Technical data

Admitted power supply	F10600118 230V-50/60Hz +/-10% - F10610118 115V-50/60Hz +/-10%
Power	100W
Dimensions (WxHxD)	665x520x470 mm
Weight	30 Kg
Speed range / precision / selection	0-30 rpm / ±1 rpm / 1 rpm a display
Operating mode	Continuous
Construction material	Epoxy painted metal structure
Environmental ambient temperature	5 - 40 °C (41 - 104°F)
Storage temperature range	-10... + 60 °C (14 - 140°F)
Pollution degree	2
Overvoltage category	II
Sound level	35 dBa
Max humidity	80%
Altitude	Up to 2000 m.
Fuses	F10600118 2xT1 A L 250 V - F10610118 2xT1 A L 250 V

6. Accessories / Spare parts

A00001023	Arm holding 8 1 l bottles	A00001022	Polyethylene bottles, 1 l
A00001024	Glass bottles with grinded glass cap, 2 l	A00001021	Polyethylene bottles, 2 l

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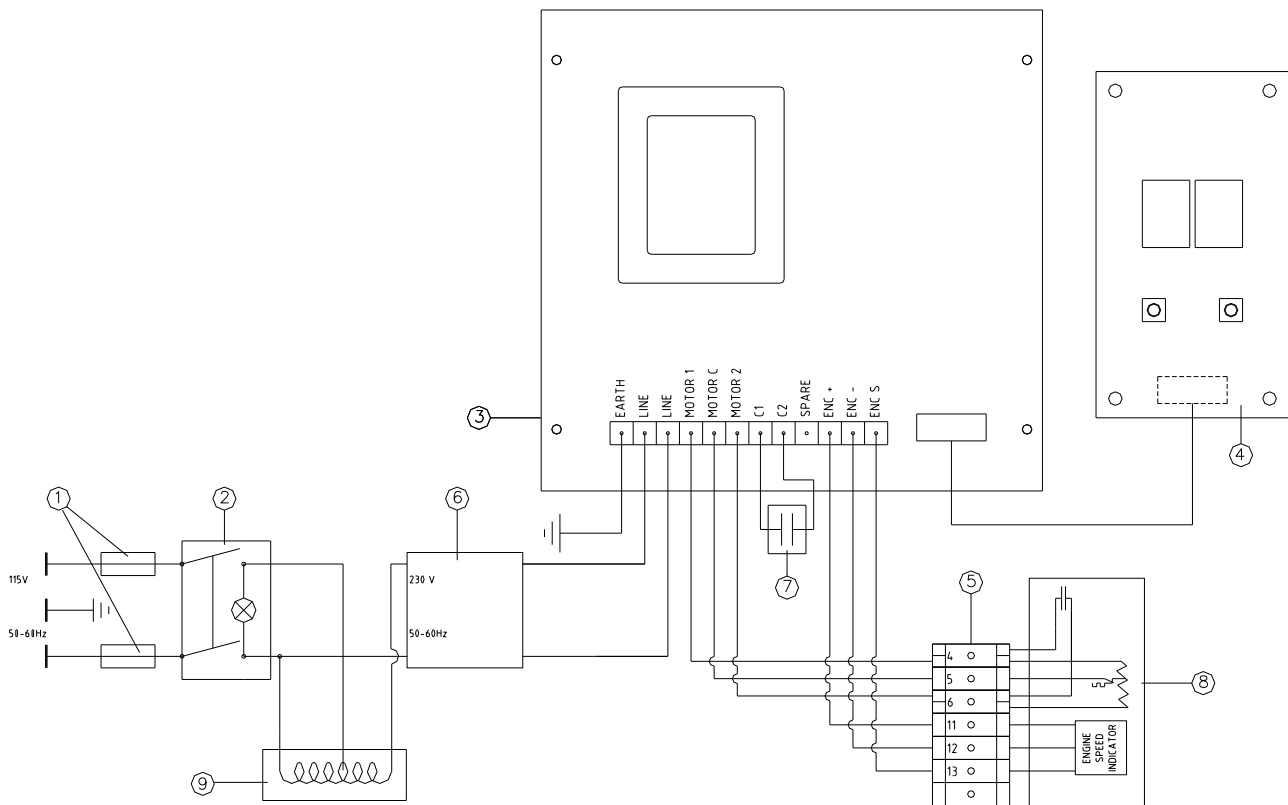
We, the manufacturer VELP Scientifica, under our responsibility declare that the product is manufactured in conformity with the following standards:

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EN 61010-2-051

EN 61326-1

and satisfies the essential requirements of the following directives:

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9. Declaration of conformity ^{UK} _{CA}

We, the manufacturer VELP Scientifica, under our responsibility declare that the product is manufactured in conformity with the following regulations:

S.I. 2016/1101 Electrical Equipment (Safety) Regulations 2016

S.I. 2016/1091 Electromagnetic Compatibility Regulations 2016

according to the relevant designated standards:

EN 61010-1 Safety requirements for electrical equipment for measurement, control, and laboratory use
Part 1: General requirements

EN 61010-2-051 Particular requirements for laboratory equipment for mixing and stirring

EN 61326-1 Electrical equipment for measurement, control and laboratory use - EMC requirements -
Part 1: General requirements

and satisfies the essential requirements of regulations:

S.I. 2008/1597 Supply of Machinery (Safety) Regulations 2008

S.I. 2012/3032 Restriction of the Use of Certain Hazardous Substances in Electrical and Electronic Equipment Regulations 2012

S.I. 2013/3113 Waste Electrical and Electronic Equipment Regulations 2013

plus modifications.

Thank you for having chosen VELP!

Established in 1983, VELP is today one of the world's leading manufacturer of analytical instruments and laboratory equipment that has made an impact on the world-wide market with Italian products renowned for innovation, design and premium connectivity. VELP works according to **ISO 9001**, **ISO14001** and **ISO 45001** Quality System Certification.

Our instruments are manufactured in Italy according to the IEC 1010-1 and CE regulation.

Our product lines:

Analytical instruments

Elemental Analyzers
Digestion Units
Distillation Units
Solvent Extractors
Fiber Analyzers
Dietary Fiber Analyzers
Oxidation Stability Reactor
Consumables

Laboratory Equipment

Magnetic Stirrers
Heating Magnetic Stirrers
Heating Plates
Overhead stirrers
Vortex mixers
Dispersers
COD Thermoreactors
BOD and Respirometers
Cooled Incubators
Flocculators
Overhead Shakers
Turbidimeter
Open Circulating Baths
Pumps

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VELP Scientifica Srl