**Instruction Manual**  
**Manuale di istruzioni**  
**Manuel d'instructions**  
**Manual de instrucciones**  
**Bedienungsanleitung**  
**指导手册**

## **SP311 Peristaltic Pump**

F40200002, F40200006, F40200012, F40200060

### **General Information / Informazioni Generali / Informations Générales / Información General / Allgemeine Hinweise / 一般信息**

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**This unit has been designed and manufactured in compliance with the following standards:**

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Safety requirements for electrical equipment for measurement, control and for laboratory use **IEC/EN 61010-1**

Electrical equipment for laboratory use **UL 61010-1**

General requirement - Canadian electrical code **CAN/CSA-C22.2 No.61010-1**

VELP reserves the right to modify the characteristics of its products with the aim to constantly improving their quality.

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## 1. Instruction

EN

Connect the pump to mains 230V/50Hz plus ground.

Stirring is turned on by the illuminated switch (A) located on the front panel.

Stirring speed is selected turning the electronic control knob (B).

The equipment is equipped by 2 fuses connected in the socket located on rear panel (C).

A blown fuse can be changed after removing rear connection socket and opening the small cover, using the tip of a screwdriver.

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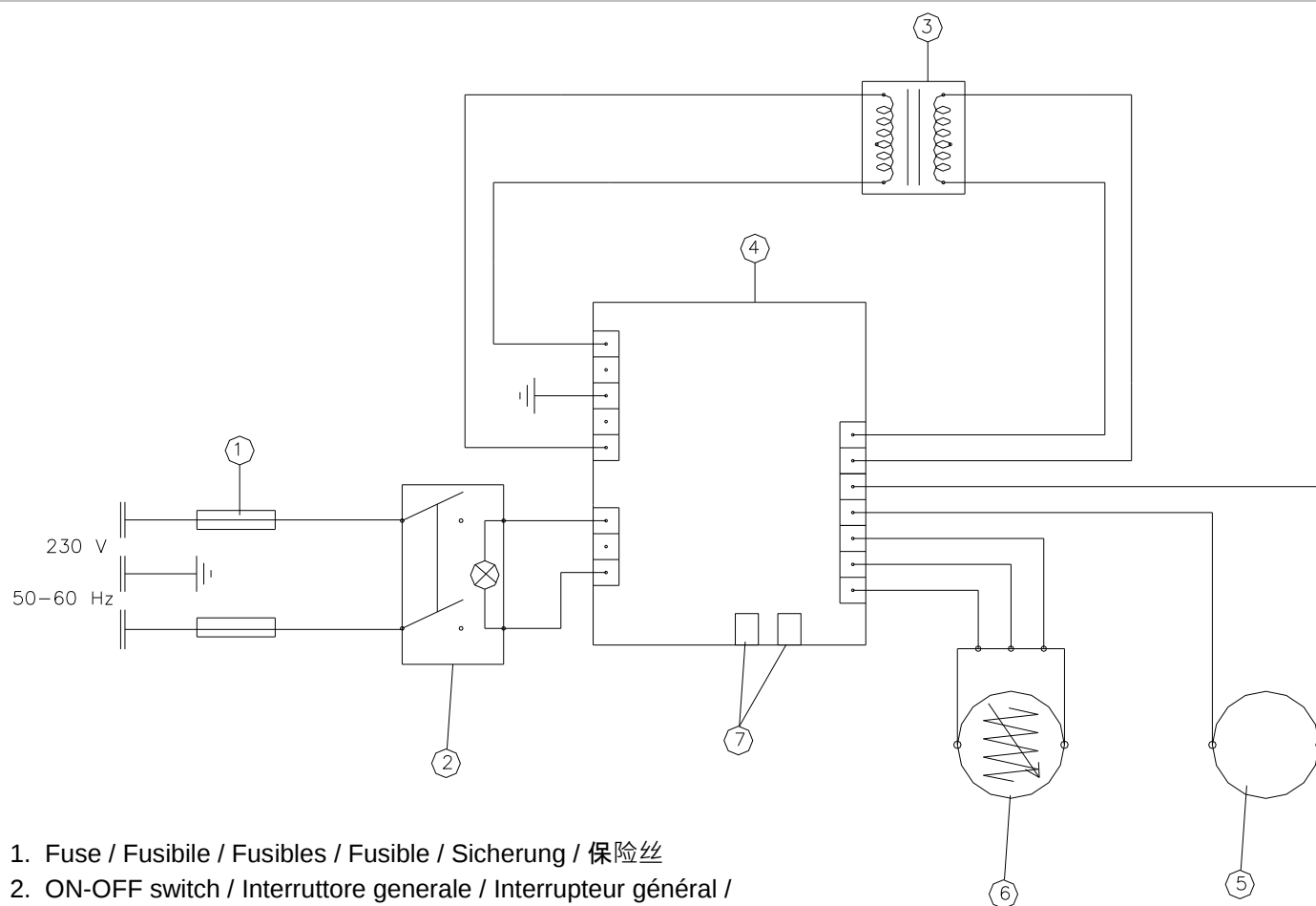
We, the manufacturer VELP Scientifica, under our responsibility declare that the product is manufactured in conformity with the following standards:

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EN 61010-1

EN 61326-1

and satisfies the essential requirements of the following directives:

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## 5. Declaration of conformity

We, the manufacturer VELP Scientifica, under our responsibility declare that the product is manufactured in conformity with the following regulations:

S.I. 2016/1101      Electrical Equipment (Safety) Regulations 2016

S.I. 2016/1091      Electromagnetic Compatibility Regulations 2016

according to the relevant designated standards:

EN 61010-1      Safety requirements for electrical equipment for measurement, control, and laboratory use  
Part 1: General requirements

EN 61326-1      Electrical equipment for measurement, control and laboratory use - EMC requirements -  
Part 1: General requirements

and satisfies the essential requirements of regulations:

S.I. 2008/1597      Supply of Machinery (Safety) Regulations 2008

S.I. 2012/3032      Restriction of the Use of Certain Hazardous Substances in Electrical and Electronic  
Equipment Regulations 2012

S.I. 2013/3113      Waste Electrical and Electronic Equipment Regulations 2013

plus modifications.

## Thank you for having chosen VELP!

Established in 1983, VELP is today one of the world's leading manufacturer of analytical instruments and laboratory equipment that has made an impact on the world-wide market with Italian products renowned for innovation, design and premium connectivity. VELP works according to **ISO 9001**, **ISO14001** and **ISO 45001** Quality System Certification.

Our instruments are manufactured in Italy according to the IEC 1010-1 and CE regulation.

Our product lines:

### Analytical instruments

Elemental Analyzers  
Digestion Units  
Distillation Units  
Solvent Extractors  
Fiber Analyzers  
Dietary Fiber Analyzers  
Oxidation Stability Reactor  
Consumables

### Laboratory Equipment

Magnetic Stirrers  
Heating Magnetic Stirrers  
Heating Plates  
Overhead stirrers  
Vortex mixers  
Dispersers  
COD Thermoreactors  
BOD and Respirometers  
Cooled Incubators  
Flocculators  
Overhead Shakers  
Turbidimeter  
Open Circulating Baths  
Pumps

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