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Before using the unit, please read the following instruction manual carefully. An improper use of the device and not as written in the manual can be dangerous for safety.

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The instrument is classified as electrical/electronic apparatus and must be disposed of accordingly. The unit is subject to waste separation and cannot be disposed of as urban waste under EEC directive 2002/96/CE. For more information please contact the relative division of your local town council.

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This unit must be used for laboratory applications only and for internal use only.

The manufacturer declines all responsibility for any use of the unit that does not comply with these instructions.

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This unit has been designed and manufactured in compliance with the following standards:

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UL 61010-1

CAN/CSA-C22.2 No.61010-1

VELP reserves the right to modify the characteristics of its products to constantly improve their quality.

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The exclusive wireless data transmission from the RESPIROMETRIC Sensor to the Wireless DataBox™ can be set at selected intervals ranging from 2 hours to 48 hours in the case of BOD/BMP analysis or from 10 minutes to 48 hours for pressure analysis. The sampling time value, depends on the total duration of analysis, till 180 days. RESPIROMETRIC Sensor automatically stores the most recent values. Data transmission is ensured even when the Wireless DataBox™ is not connected to the PC. The Wireless DataBox™ can store unlimited results for up to 50 samples.

2. Assembling and installing

Upon receipt and after having removed the packaging, please check the integrity of the instrument. The box includes:

- Wireless DataBox™
- Power USB cable
- RESPIROSoft™
- Power supply USB

The Wireless DataBox™ comes already prepared and doesn't need any action for the assembling it. Using the USB cable is possible connect it to the PC or to the power supplier when analyses are in progress. In this way is possible record the dates inside the unit without use the computer.

On the USB-pen drive is possible to find the driver if required by Windows after the first connection to the PC.

3. RESPIROSoft™

3.1 Installation

General rules for installing the RESPIROSoft™ Software:

In order to install the RESPIROSoft™ Software correctly, administrator rights for the PC must be available. The PC on which the RESPIROSoft™ Software will be installed must not be overloaded with other software. This is to prevent unstable operation of the RESPIROSoft™ Software.

To install the RESPIROSoft™ Software:

1. Place the USB-pen Drive in the USB port.
2. Start the installation running the function **setup.exe** from the USB-pen drive.
3. The installation guide is displayed. Click **Next** and follow the instructions. The RESPIROSoft™ icon will be created on the screen automatically.
4. Connect the Wireless DataBox™ to the PC.
5. After a few seconds the PC acknowledges the presence of a new hardware and will automatically start the installation procedure of the drivers. When requested, specify the path of the **Driver** folder on the USB-pen drive and proceed with installation following the instructions that appear on the screen.
6. Re-boot the computer to put the settings provided by the installation procedure into effect.

3.2 Turning on

Before turning on the instrument, all the steps described in the previous paragraphs must be carried out:

1. Turn on the PC and the screen.
2. Check that the RESPIROSoft™ has been installed.
3. Connect the Wireless DataBox™ to the PC using the USB cable.
4. Launch the RESPIROSoft™ (shortest way > double-click on the appropriate icon on the screen) (Fig. 1).
5. The "User" level password is "velp". Enter the password and press OK (Fig. 2).
6. The main window will appear (Fig. 3).

Fig. 1

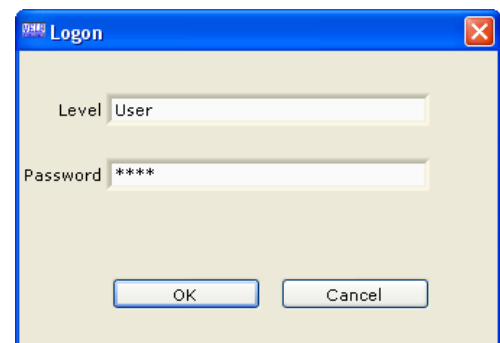
Fig. 2

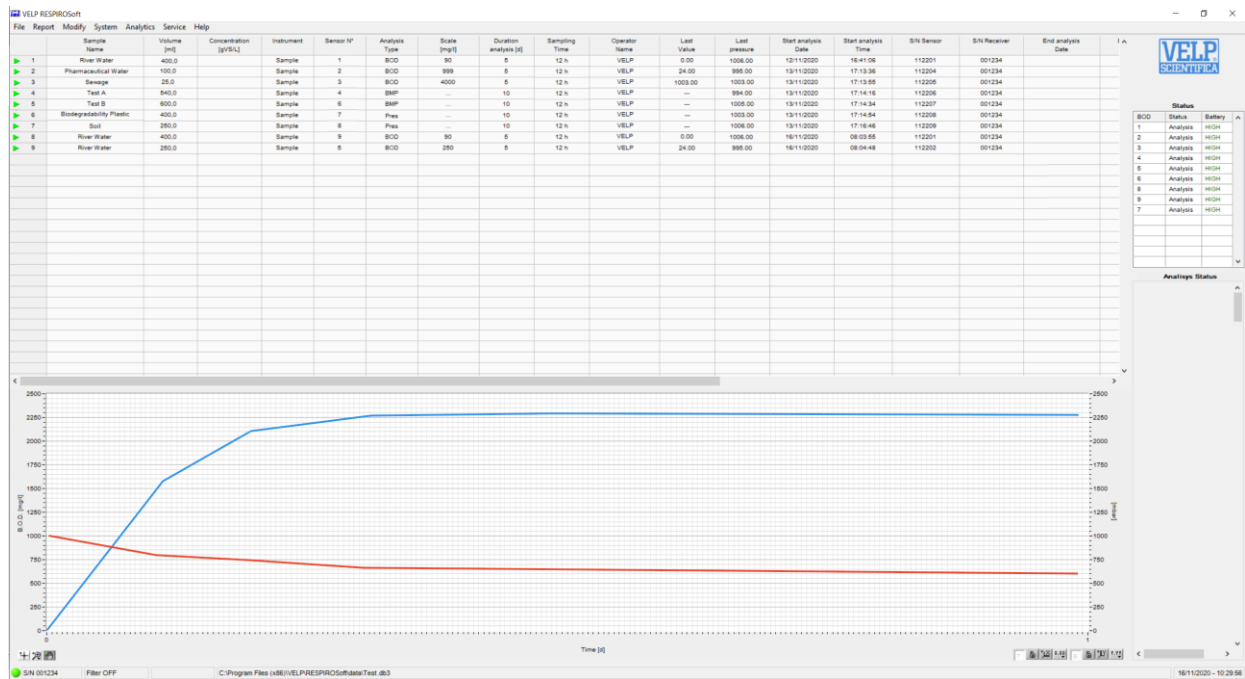
Fig. 3

4. RESPIROsoft™ Rapid Guide

This page shows an abstract of the RESPIROsoft™.

The purpose is to help the user identify the various possibilities offered by the instrument and how to reach them easily.

Menu	Submenu	Description	
File	New	Create a new analysis data base and "Laboratory information" file.	
	Open	Load an already existing database.	
	Save as	Save an analysis data base (including "Laboratory information" file) assigning a file name. If the file name already exists, it can be overwritten.	
	Laboratory information	Enter the details of the laboratory concerned. This information will be included in the test report.	
	Exit	Exit from the programme and return to the desktop.	
Report	Create Report	Create a single or multiple test report with calculation of average and RSD values.	
	Export Data	Export data of the rows selected to an .xls or .txt or csv file.	
Modify	Filter	Create database filters.	
	Standard column	Reset the order of the database columns to the default setting.	
	Configure	Configure the sequence of the database columns, show/hide columns.	
	Remove	Remove one or more rows from the database.	
	Select all	Select all the rows of the database containing a value.	
System	Languages	Change the language of the program.	
Analytics	Methods	Display, modify, create or delete analytical methods.	
	Standards	Display, modify, create or delete standards.	
	BMP Calculator	Calculator for BMP Analysis	
Service	Connect Receiver	Manual search for Wireless DataBox™.	
	Disconnect Receiver	Disconnect the Wireless DataBox™ from the PC.	
	Ermes	Enable/Disable ERMES connection	
	Start AP	Start the AP mode on Wireless DataBox	
	Acoustic signals	Select the operations that activate an acoustic signal.	
	Logon	Modify the software password and entry level.	
	Diagnostic	Perform a diagnosis of the status of the instrument.	
	Control panel	For VELP or authorized technical staff only.	
	Help	Software activation	Under construction.
		About	Contact VELP Scientifica directly by e-mail.
Manual		PDF operating manual.	

5. DataBox™ Connection

In order to connect the Wireless DataBox™ to the PC, it is necessary to use the USB cable, supplied with the unit.

1. After having connected it to the PC, select **Connect Receiver** from the **Service** menu (Fig. 4).
2. Using the “Connect Receiver” menu with the Automatic Search option, it is possible to find and connect automatically the Wireless DataBox™ (Fig. 5).
3. When the unit is connected to the RESPIROSoft™, the software shows the S/N and a green led on the left bottom (Fig. 6).

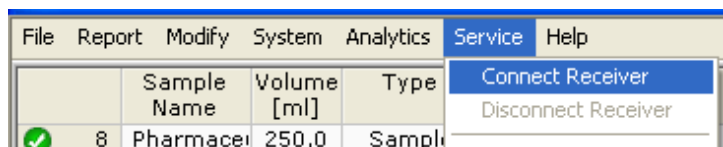
Fig. 4

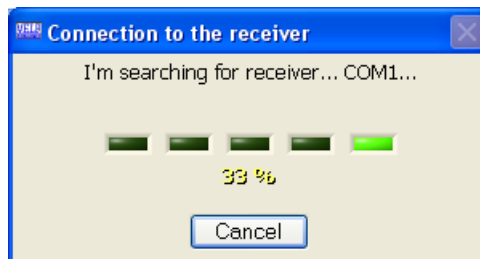
Fig. 5

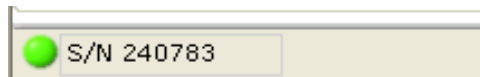
Fig. 6

6. RESPIROSoft™ Workflow

6.1 Preparing for analysis

About the sample preparation, please see the analytical methods section of RESPIROMETRIC System operating manual. In the Analytics/Methods menu, it is possible to see a summary of methods reported in the analytical methods section of RESPIROMETRIC System operating manual.

6.1.1 Filling the database

Several databases can be created and stored in the PC. Each database can store max 1000 analyses. Select **File/New** to create a new database or select **File/open** to load an already existing database. Each database is connected to a “Laboratory information” file (Fig. 7).

The information in this file will be included in the test reports. The company logo can also be uploaded in order to customize the test reports. The “Laboratory Information” file can be modified as many times as necessary.

Before starting the analysis, some of the columns must be filled in. The following database information is mandatory and **must be entered** before starting the analysis:

Sample Name	Name of sample (alphanumeric field)
Test Type	Type of analysis to be performed
Sensor N°	Number used for identify the Sensor
Analysis Type	BOD, BMP or PRESSURE
Scale	Scale used for the analysis if necessary
Duration Analysis	Total days of analysis
Sampling Time	Sampling time for the analysis

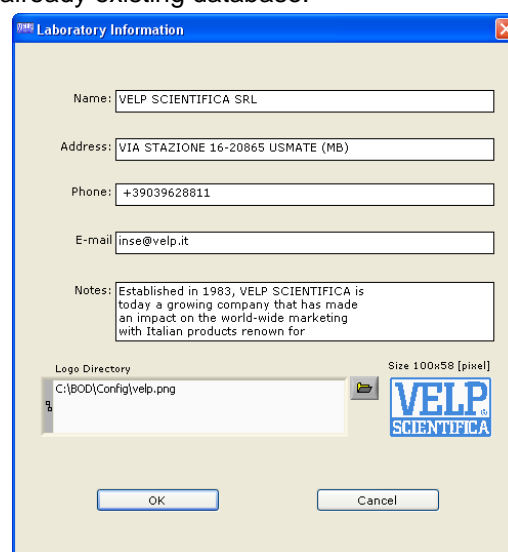
Fig. 7

Once the analysis has been carried out the content of the database columns cannot be modified. All the other database columns are filled automatically by the instrument at the end of analysis. The information can be entered once and copied using **Copy/Add** (mouse right button) to add new row. The new row will be added in sequence. In case of BMP, also Volume, Vbottle, MS and Incubator T must be compiled if the BMP value is required.

6.1.2 Choosing the sample “Test Type”

The “Type” column offers the following options (Fig. 8):

Sample	Analysis on a sample.
Standard	When is necessary to compare the last value with a range of result, e.g. with a standard in order to verify the RESPIROMETRIC Sensor

In **Analytics/Standards**, it is possible to find some standard that are suggested for the test.

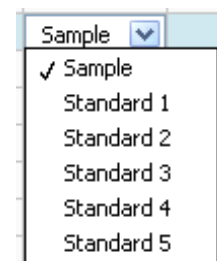
Fig. 8

6.1.3 Choosing the “Sensor N°”

To use a Sensor with RESPIROSoft™, is necessary to register it with a number that is used to identify the unit. In order to register it, it is necessary to keep pressed “SET” on the RESPIROMETRIC Sensor until appears the following screen on the PC (Fig. 9).

Give a number to the Sensor and press OK.

In the column **Sensor N°** it is possible to select the units that have been already registered.


Fig. 9

The registration of sensors is necessary only when use for the first time a RESPIROMETRIC Sensor. with RESPIROSoft™. Otherwise, if the unit was already used, the number is already registered.

6.1.4 Choosing the “Scale”

Choose the **Scale** in case of BOD analysis, with the following options (Fig. 10):

Scale	Volume of sample
A: 0 ÷ 4000 mg O ₂ /l	25 ml
B: 0 ÷ 999 mg O ₂ /l	100 ml
C: 0 ÷ 600 mg O ₂ /l	150 ml
D: 0 ÷ 250 mg O ₂ /l	250 ml
E: 0 ÷ 90 mg O ₂ /l	400 ml

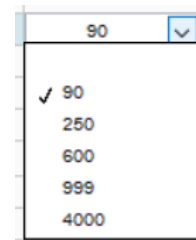6.1.5 Choosing the “Duration Analysis”

Using the **Duration Analysis**, it is possible to select different value for the duration of analysis. This value depends on the type of analysis. BOD, from 5 to 30 days; BMP, from 10 to 60 days; PRESSURE, from 1 to 180 days.

6.1.6 Choosing the “Sampling Time”

Using the **Sampling Time**, it is possible to select different period of sampling. This value depends to the Duration Analysis selected before.

6.2 Performing analysis

In order to start the analysis, please refer to the RESPIROMETRIC Sensor operating instructions.

Before performing an analysis, all the steps described above must be carried out.

The graph at the bottom of the display always refers to the analysis selected and not necessarily to the analysis in progress.

6.2.1 Stopping analysis

In order to stop the analysis, please refer to the RESPIROMETRIC Sensor operating instructions.

Once the analysis has been stopped, the icon  will be showed on the left part of analysis.

6.2.2 Alarms

In order to see all the possible alarms, please refer to the RESPIROMETRIC Sensor operating instructions.

RESPIROSoft™ gives on screen a short explanation of the problem.

6.3 Data evaluation

In order to evaluate the data obtained, the software offers the following possibilities:

- Configuring the data base
- Filtering the data base
- Exporting data

The indicator in the first column of the data base gives the status of each analysis in the series:

Indicator	Meaning
	Analysis ended correctly
	Analysis terminated manually with a reset by user.
	Analysis ended correctly but there are some notes (e.g. last value outside from range)

6.3.1 Configuring the data base

The database consists of a set of data entered by the operator as well as data calculated by the software after analysis. Select **Modify** to modify the configuration of the database.

Filter	Filter data.
Standard column	View columns in standard configuration.
Configure	Change the order of columns (up and down arrows) or conceal columns (left and right arrows). Alternatively, press the right key of the mouse and select Configure (<i>Fig. 11</i>).
Remove	Delete one or more rows from the data base. Alternatively, press the right key of the mouse and select Remove.
Select all	Selecting all the rows in the database

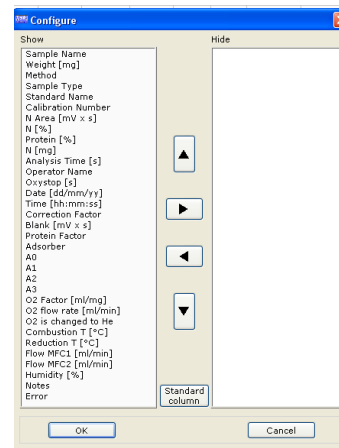
Fig. 11

6.3.2 Filtering the database

The data shown can be filtered using the filtering function. Select **Modify/filters**. The following window appears (*Fig. 12*).

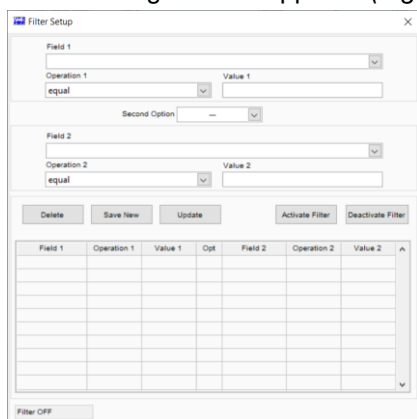
Fig. 12

Select filters as follows:

- 1- choose the fields to be filtered and the value required
- 2- press Save New in order to save the filter
- 3- Apply with Activate Filter or Deactivate Filter.

The filter status ("Filter ON" or "Filter OFF") is shown at the bottom of the main window.

6.3.3 Exporting data

Information can be exported from the data base in Excel, Text or Lims format by selecting **Report/Export Data**.

Report Type EXCEL	an .xls file is created; graphs not exported.
Report Type TXT	a .txt file is created; graphs not exported.
Report Type LIMS (<i>Fig. 13</i>)	a .csv file format is created which can be easily read by the majority of laboratory management systems (LIMS). Select the value/unit to be exported (see below). Select the configuration of the exported data from the options available. It is possible to attribute the comma, semicolon or tab as delimiting character and the comma or decimal point as the decimal indicator depending on the type of the Lims. Press OK to confirm.

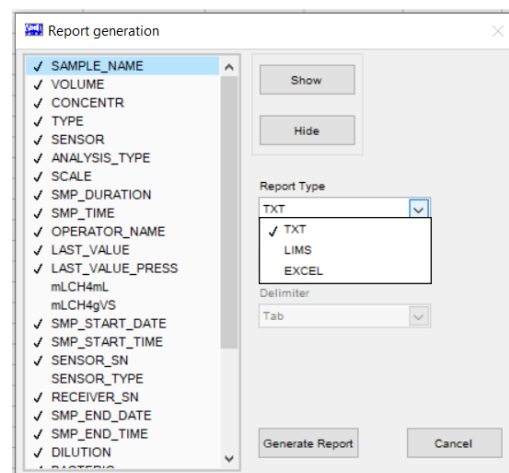
Fig. 13

6.4 Test reports

Information can also be exported from the data base to a predefined Excel file called Test report.



It is possible to create:

Test report for single analysis

Highlight a row and select **Report/Create**.

Test report for multiple analyses

Highlight the rows and select **Report/Create**.

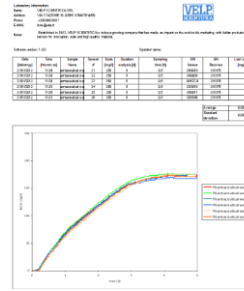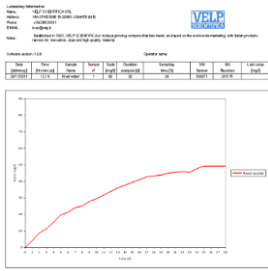NOTE: the test report is an open file and can be modified.

NOTE: if the analyses selected have not been performed under the same conditions (different scale, duration and sampling time), the test report does not display the results in Average and SD.

7. Software installation requirements

Operating system	Windows 7 (SP 1), Windows 8 or Windows 10
Hardware	Core 2 Duo, RAM 4 GB
Available space on the hard disk	100 MB
Recommended graphic resolution	1024 x 768

8. Technical Data

Dimensions (WxHxD)	125x25x70 mm (4.9x1.0x2.7 in)	
Weight	85 g (0.04 lb)	
RESPIROMETRIC Sensor management	Up to 48 per Wireless DataBox™	
Power supply	5 V – USB Cable	
External power supply	100-240V±10% 50-60Hz 200mA	
Power	0,225 W	
Overvoltage category	II	
Pollution degree	2	
Max. altitude	2000 m	
Max. humidity	85%	
Admitted room temperature	Storing -25/+65° C	Working +5/+40°C
Frequency WiFi	2.4GHz	
Protocols WiFi	802.11 b/g/n	

9. Spare Parts

10005112	USB cable A/Micro B 1.5m
40001862	Power supply 100-240V/5V no cord/plug
10003082	UE plug for power supply

10. Warranty

The unit is guaranteed against production defects for **25 months** from our invoice date. In accordance with this guarantee VELD Scientifica undertakes to repair any instruments resulting as faulty due to the quality of the materials used or poor workmanship. Faults arising due to inexpert handling/use or carelessness will not be replaced or repaired under. For more details please contact your Distributor.

Exclusions:

The warranty will be considered null and void for faults resulting from:

- inexperience and carelessness of the customer.
- repairs, maintenance or replacements carried out by unauthorized third parties.
- use of non-original spare parts.

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11. Declaration of conformity

We, the manufacturer VELP Scientifica, under our responsibility declare that the product is manufactured in conformity with the following standards:

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EN 62311:2008	Assessment of electronic and electrical equipment related to human exposure restrictions for electromagnetic fields (0 Hz - 300 GHz)
EN 301 489-1 V2.2.0	Electromagnetic compatibility and Radio spectrum Matters (ERM); ElectroMagnetic Compatibility (EMC) standard for radio equipment and services; Part 1: Common technical requirements
EN 301 489-17 V3.2.0	Electromagnetic compatibility and Radio spectrum Matters (ERM); ElectroMagnetic Compatibility (EMC) standard for radio equipment; Part 17: Specific conditions for Broadband Data Transmission Systems
EN 300 328 V1.9.1	Electromagnetic compatibility and Radio spectrum Matters (ERM); Wideband transmission systems; Data transmission equipment operating in the 2,4 GHz ISM band and using wide band modulation techniques; Harmonized EN covering the essential requirements of article 3.2 of the R&TTE Directive
2011/65/EU (RoHS)	Restriction of the use of certain hazardous substances
2012/19/EU (RAEE)	On waste electrical and electronic equipment
EN 61010-1 (2010)	Safety Requirements for Electrical Equipment for Measurement, Control, and Laboratory Use - Part 1: General Requirements.

and satisfies the essential requirements of the following directives:

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Thank you for having chosen VELP!

Established in 1983, VELP is today one of the world's leading manufacturer of analytical instruments and laboratory equipment that has made an impact on the world-wide market with Italian products renowned for innovation, design and premium connectivity. VELP works according to **ISO 9001**, **ISO14001** and **OHSAS 18001** Quality System Certification.

Our instruments are manufactured in Italy according to the IEC 1010-1 and CE regulation.

Our product lines:

Analytical instruments

Elemental Analyzers
Digestion Units
Distillation Units
Solvent Extractors
Fiber Analyzers
Dietary Fiber Analyzers
Oxidation Stability Reactor
Consumables

Laboratory Equipment

Magnetic Stirrers
Heating Magnetic Stirrers
Heating Plates
Overhead stirrers
Vortex mixers
Homogenizers
COD Thermoreactors
BOD and Respirometers
Cooled Incubators
Flocculators
Overhead Shakers
Turbidimeter
Radiation Detector
Open Circulating Baths
Pumps

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